Level

Pressure

Flow

Temperature

Liquid
Analysis

Registration

Systems
Components


Services

Solutions

Kurzanleitung / Brief Operating Instructions

Memocheck Plus CYP01D, Memocheck CYP02D

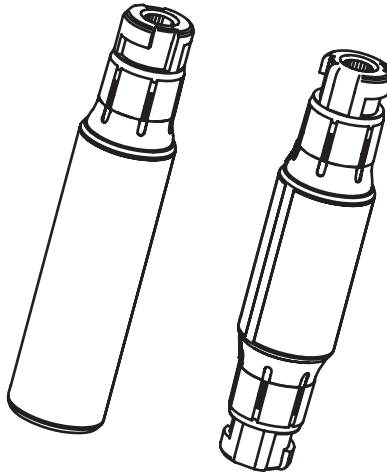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en

Tool for the qualification of measuring systems and service tool for quick checks of measuring systems

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English

| | | |
|-----------|---|-----------|
| 6 | Safety instructions | 14 |
| 6.1 | Designated use | 14 |
| 6.2 | Operational safety | 14 |
| 6.3 | Safety instructions for hazardous areas | 15 |
| 7 | Identification | 15 |
| 7.1 | Nameplate | 15 |
| 7.2 | Product structure | 16 |
| 7.3 | Identification of Ex versions | 17 |
| 7.4 | Scope of delivery | 17 |
| 7.5 | Certificates and approvals | 18 |
| 8 | Installation | 20 |
| 8.1 | Dimensions | 20 |
| 8.2 | Environment | 20 |
| 8.3 | Temperature classes of Ex versions | 20 |
| 9 | Operation | 21 |
| 9.1 | Description | 21 |
| 9.2 | Application | 22 |
| 10 | Accessories | 23 |
| 10.1 | pH glass electrodes | 23 |
| 10.2 | Transmitters | 24 |
| 10.3 | Software | 24 |

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Memocheck Plus CYP01D-*G**
Memocheck CYP02D-*G**

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Qualification tool

Memocheck Plus CYP01D

This tool serves for the qualification of measurement chains with Memosens technology and thus to verify the data communication of sensor coupling over transducers up to the process control system. Each qualification tool consists of five plug-in heads. The acquired qualification tool has been manufactured with the greatest care and observation of all technical rules. The materials used are subject to continuous quality control. Before leaving our factory, each individual plug-in head has passed a stringent test on our computerized inspection unit.

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6 Safety instructions

6.1 Designated use

Memocheck Plus CYP01D is a comprehensive tool for the qualification of measurement chains with Memosens technology. It enables you to verify data transmission from sensor coupling via transmitter to process control system. Memocheck Plus consists of five plug-in heads. Each plug-in head simulates a specific predefined sensor status.

Memocheck Plus CYP01D is **not** suitable for application directly in the process.

Memocheck CYP02D is a service tool for quick, on-site checks of measuring systems with Memosens technology.

Memocheck CYP02D is **not** suitable for application directly in the process.

Any other use than the one described here compromises the safety of persons and the entire measuring system and is, therefore, not permitted.

The manufacturer is not liable for damage caused by improper or non-designated use.

6.2 Operational safety

Memocheck has been designed and tested according to the state of the art and left the factory in perfect functioning order.

Relevant regulations and European standards have been met.

As the user, you are responsible for complying with the following safety conditions:

- Installation instructions
- Local prevailing standards and regulations.

6.3 Safety instructions for hazardous areas

Caution!

- The ATEX-certified sensor simulator types Memocheck Plus CYP01D-....G and Memocheck CYP02D-...G may only be connected via the CYK10-G..1 measuring cable to
 - the certified intrinsically safe digital sensor output circuits of the field transmitter Mycom S, type CPM153-G5..... and Mycom S, type CPM153-G6..... in accordance with the EC type examination certificate DMT 01 ATEX E 174
 - the certified intrinsically safe digital sensor output circuits of the transmitter Liquiline M, type CM42-MG..... in accordance with the EC type examination certificate EX 5 05 03 30266 012.
 The related EC declaration of conformity forms an integral part of this document.
- The FM/CSA-certified sensor simulator types Memocheck Plus CYP01D-....G and Memocheck CYP02D-...G may only be connected via the CYK10-G..1 measuring cable to
 - the certified intrinsically safe digital sensor output circuits of the field transmitter Mycom S, type CPM153-O/P/S5..... and Mycom S, type CPM153-O/P/S6..... in accordance with the control drawings of the transmitter (see transmitter documentation)
 - the certified intrinsically safe digital sensor output circuits of the transmitter Liquiline M, type CM42-MP/MS..... in accordance with the control drawings of the transmitter (see transmitter documentation).
- The electrical connection of the CYK10 measuring cable must correspond to the wiring diagram of the transmitter.
- The sensor simulators may not be operated under critical electrostatic process conditions. Strong vapour and dust streams acting directly on the plug-in system must be avoided.
- The maximum permitted cable length is 100 m.
- Ex versions are indicated by an orange-red ring in the plug-in head.


7 Identification

7.1 Nameplate

Fig. 9: Memocheck Plus CYP01D nameplate example (pH glass electrode, value: pH 14)

a0005646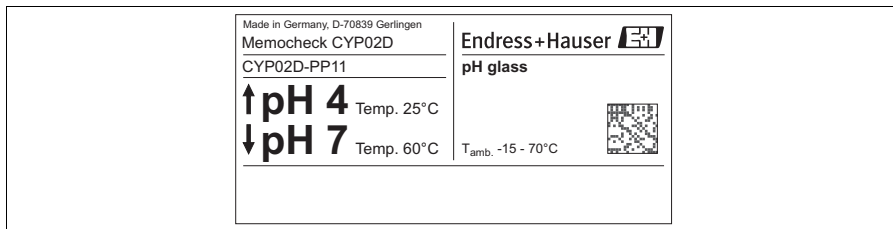

a0005647

Fig. 10: Memocheck CYP02D nameplate example (pH glass + pH glass, values: pH 4 + pH 7)

7.2 Product structure

7.2.1 Memocheck Plus CYP01D

| Parameter | | | |
|-------------|--|--|---------------------|
| PG | pH; for glass electrodes | | |
| Version | | | |
| 1 | Standard | | |
| Certificate | | | |
| A | Quality certificate | | |
| Approval | | | |
| G | ATEX II 2G EEx ia IIC T6 FM IS NI Cl. 1 Div. 1&2 Group A-D T6 CSA IS NI Cl. 1 Div. 1&2 Groups A-D T6 | | |
| 1 | Non-hazardous area | | |
| CYP01D- | | | complete order code |

7.2.2 Memocheck CYP02D

| Parameter | | | |
|-----------|---|--|---------------------|
| PP | pH glass + pH glass | | |
| Version | | | |
| 1 | Standard | | |
| Approval | | | |
| G | ATEX II 2G EEx ia IIC T6 FM IS NI Cl. 1 Div. 1&2 Groups A-D T6 CSA IS NI Cl. 1 Div. 1&2 Groups A-D T6 | | |
| 1 | Non-hazardous area | | |
| CYP02D- | | | complete order code |

7.3 Identification of Ex versions

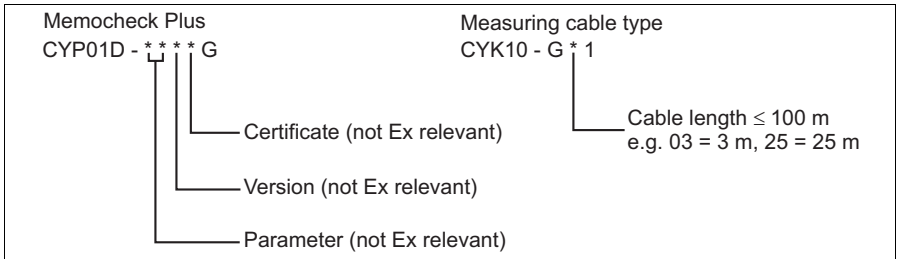
Fig. 11: Identification of CYP01D

a0007191-en

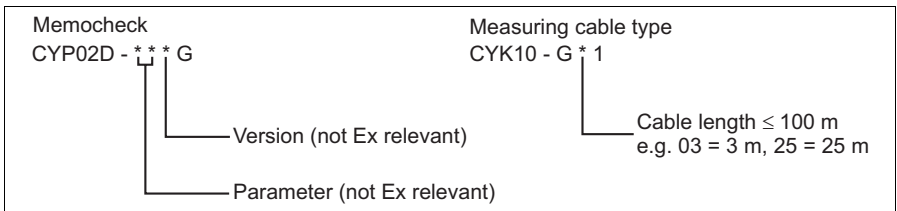
Fig. 12: Identification of CYP02D

a0007192-en

7.4 Scope of delivery

7.4.1 Memocheck Plus CYP01D

The scope of delivery includes:

- 5 Plug-in heads in a case
- 1 Quality certificate
- 1 Brief Operating Instructions


7.4.2 Memocheck CYP02D

The scope of delivery includes:

- 1 Memocheck CYP02D with 2 plug-in heads
- 1 Brief Operating Instructions

7.5 Certificates and approvals

7.5.1 Declaration of conformity for Ex versions

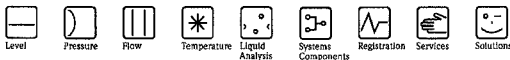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Endress+Hauser 
People for Process Automation

a0007227

7.5.2 Quality certificate of Memocheck Plus CYP01D

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Qualification tool

Memocheck Plus CYP01D

This tool serves for the qualification of measurement chains with Memosens technology and thus to verify the data communication of sensor coupling over transducers up to the process control system. Each qualification tool consists of five plug-in heads.

The acquired qualification tool has been manufactured with the greatest care and observation of all technical rules. The materials used are subject to continuous quality control.

Before leaving our factory, each individual plug-in head has passed a stringent test on our computerized inspection unit.

Serial numbers and fixed value pairs of the reference means delivered:

| | |
|---|----------------------------|
| Bestellcode / order code | CYP01D-PG1A1 |
| Seriennummer Set / serial number of set | 7C005A05E00 |
| simulierter Fehler / simulated error | Glasbruch / glass breakage |
| Prüfdatum / test date | 13.12.2005 |
| Prüfer / inspector | |
| Unterschrift / signature | |

| | Seriennummer / serial number | pH-Wert / pH value | Temperatur / temperature | Qualitätskontrolle / Quality Control |
|----|------------------------------|--------------------|--------------------------|--------------------------------------|
| 1 | 7A0D2505E00 | 0 | -10 °C | bestanden / passed |
| 2 | 7A0D2605E00 | 4 | 25 °C | bestanden / passed |
| 3 | 7A0D2705E00 | 7 | 60 °C | bestanden / passed |
| 4 | 7B002D05E00 | 10 | 90 °C | bestanden / passed |
| 5* | 7B002E05E00 | 14 | 135 °C | bestanden / passed |

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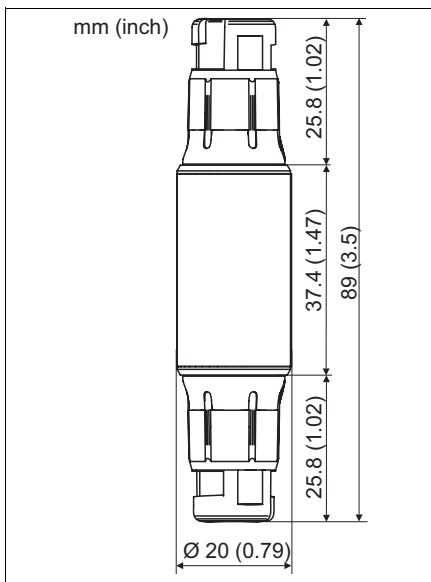
Fig. 14: Memocheck CYP02D

8.2 Environment

The permissible ambient temperature ranges from: -15 to 70 °C (5 to 158 °F)

8.3 Temperature classes of Ex versions

| Name | Type | | | | | | | Ambient temp. T_a at temperature class (T_n) | Cat. |
|----------------|------|-----|---|----|---|---|---|--|-------|
| Memocheck Plus | CYP | 01D | - | ** | * | * | G | -15 °C $\leq T_a \leq +70$ °C (T6) | II 2G |
| Memocheck | CYP | 02D | - | ** | * | | G | -15 °C $\leq T_a \leq +70$ °C (T6) | II 2G |

9 Operation

9.1 Description

9.1.1 Memocheck CYP01D

Memocheck Plus CYP01D is a set of 5 various sensor plug-in heads that provide data for the simulation of pH values. Each plug-in head simulates the state of a corresponding sensor. The following data are indicated and transmitted, they **cannot** be overwritten:

- Main measured value, e.g. pH
- Temperature
- Error message, e.g. glass breakage: One plug-in head transmits an error message.
The error message can be displayed or suppressed by switching the corresponding transmitter function, e.g. glass SCS, on or off.

Transmitted measured values for pH glass electrodes (CYP01D-PG***)

| Plug-in head | pH value | Temperature | Glass SCS |
|--------------|----------|-----------------|-----------|
| 1 | 0 | -10 °C (14 °F) | o.k. |
| 2 | 4 | 25 °C (77 °F) | o.k. |
| 3 | 7 | 60 °C (140°F) | o.k. |
| 4 | 10 | 90 °C (194 °F) | o.k. |
| 5 | 14 | 135 °C (275 °F) | Error |

Standard data set of process data

All Memocheck Plus CYP01D plug-in heads contain the complete data set of a digital sensor. This means they provide the same amount of information for display as digital sensors. The stored data set **cannot** be overwritten which guarantees a safe use as qualification tool at any time.

Among others, the following standard process data are stored in the plug-in heads:

- Total operating hours: 8760 h
- Number of sterilisations: 2
- Max. temperature: 135 °C (275 °F)
- Operating hours above 80 °C (176 °F): 14 h
- Operating hours above 100 °C (212 °F): 9 h
- Operating hours > 300 mV: 15 h
- Operating hours < -300 mV: 0 h
- First commissioning: 01.06.2005 12:00

9.1.2 Memocheck CYP02D

Memocheck CYP02D consists of two agglutinated independent plug-in heads that provide value pairs for the simulation of pH values. Each plug-in head simulates the state of a corresponding sensor. The values **cannot** be overwritten.

Available combination:

pH = 4, T = 25 °C (77 °F) for glass and pH = 7, T = 60 °C (140 °F) for glass

9.2 Application

To qualify your measuring point using Memocheck Plus, proceed as follows:

1. Unplug the Memosens coupling from your sensor (see Fig. 15).
2. If you are not already using automatic temperature compensation, set your transmitter to automatic temperature compensation.
3. Plug one of your Memocheck Plus plug-in heads into the Memosens coupling (see Fig. 16).

Caution!

Make sure that the Memocheck Plus does **not** come into contact with the process.

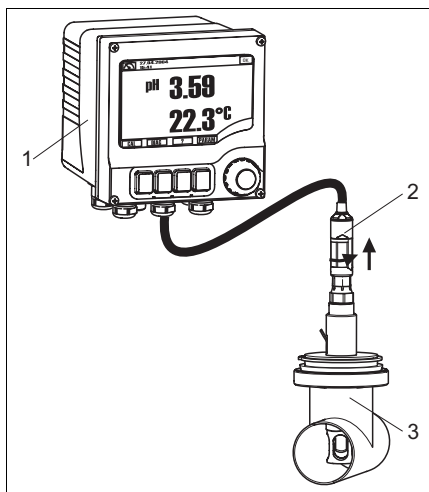
Fig. 15: Measuring point with Memosens sensor

- 1 *Liquiline M CM42 transmitter*
- 2 *Memosens coupling*
- 3 *CPA442 assembly with Memosens sensor*

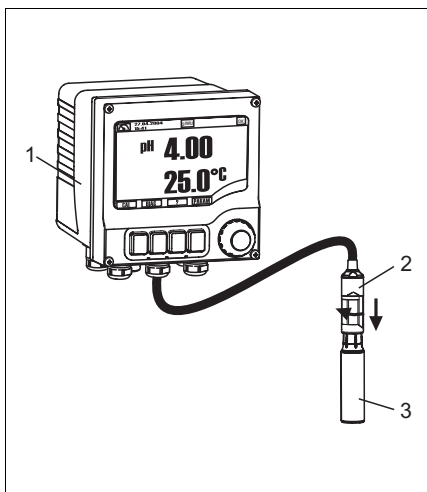
Fig. 16: Connecting Memocheck Plus CYP01D

- 1 *Liquiline M CM42 transmitter*
- 2 *Memosens coupling*
- 3 *Memocheck Plus CYP01D*

4. The fixed values of the plug-in head are automatically transmitted to the transmitter and displayed.

5. If you are using Liquiline M CM42, the transmitter switches to simulation mode ("SIMU" symbol is displayed in the status line).
6. Check if the measured value and temperature displayed by the transmitter or the process control system correspond to the given fixed values of the Memocheck Plus.
7. If you are using the fifth plug-in head of Memocheck Plus, you can check whether the error message, e.g. "glass breakage" is correctly transmitted to the transmitter or the process control system. For pH glass electrodes, you can suppress or display this error message by switching the corresponding glass SCS function on or off.
8. If you are using Liquiline M CM42, you can read out and print the Memocheck Plus data via FieldCare.
9. If the displayed values do not correspond to the given fixed values, check the following:
 - If you are using a transmitter with current outputs, make sure that the outputs are correctly wired.
 - If you are using a transmitter with Hart[®], Profibus or FOUNDATION Fieldbus communication, make sure that the communication works properly.

10 Accessories

10.1 pH glass electrodes

- Orbisint CPS11D
Digital pH sensor for process applications, with PTFE diaphragm;
Ordering acc. to product structure, see Technical Information (TI028/C07/en)
- Ceragel CPS71D
Digital pH sensor with double chamber reference system and integrated bridge electrolyte;
Ordering acc. to product structure, see Technical Information (TI245/C07/en)
- Orbipore CPS91D
Digital pH sensor with open aperture for media with high dirt load;
Ordering acc. to product structure, see Technical Information (TI375C/07/en)

10.2 Transmitters

- Liquiline M CM42-M*****
Modular two-wire transmitter, stainless steel or plastic, field or panel instrument, various Ex approvals (ATEX, FM, CSA),
Hart[®], Profibus or FOUNDATION Fieldbus available,
Ordering acc. to product structure, see Technical Information (TI381C/07/en)
- Liquisys M CPM223/253-MR/MS***
Transmitter for pH and ORP measurement, field or panel instrument,
Hart[®] or Profibus available,
Ordering acc. to product structure, see Technical Information (TI194C/07/en)
- Mycom S CPM153-*5*****
Transmitter for pH and ORP measurement, Ex or non-Ex,
Hart[®] or Profibus available
Ordering acc. to product structure, see Technical Information (TI233C/07/en)

10.3 Software

- FieldCare
Tool for Plant Asset Management
Supports Ethernet, HART, PROFIBUS, FOUNDATION Fieldbus
FieldCare Lite, order no. 56004080
FieldCare Standard, order no. SFE551-xxxx
FieldCare Professional, order no. SFE552-xxxx

www.endress.com/worldwide

Endress+Hauser 

People for Process Automation