

Technical Information

Proline Promass P 100

Coriolis flowmeter



The specialist for life sciences with an ultra-compact transmitter

Application

- Measuring principle operates independently of physical fluid properties such as viscosity or density
- Dedicated for applications under sterile conditions in the life sciences industry

Device properties

- ASME BPE, 3A and EHEDG conform & low delta ferrite
- Electropolished measuring tube in 1.4435 (316L)
- Fast recovery from CIP/SIP
- Robust, ultra-compact transmitter housing
- High ingress protection: IP69K
- Local display available

Your benefits

- Ultimate process quality – fully meets industrial requirements
- Fewer process measuring points – multivariable measurement (flow, density, temperature)
- Space-saving installation – no in/outlet run needs
- Space-saving transmitter – full functionality on smallest footprint
- Time-saving local operation without additional software and hardware – integrated web server
- Integrated verification – Heartbeat Technology™

Table of contents







Document information	4	Shock resistance	43
Symbols used	4	Vibration resistance	43
Function and system design	5	Interior cleaning	43
Measuring principle	5	Electromagnetic compatibility (EMC)	44
Measuring system	6	Process	44
Device architecture	7	Medium temperature range	44
Safety	7	Density	44
Input	7	Pressure-temperature ratings	44
Measured variable	7	Secondary containment pressure rating	48
Measuring range	8	Flow limit	48
Operable flow range	8	Pressure loss	48
Output	9	System pressure	48
Output signal	9	Thermal insulation	49
Signal on alarm	10	Heating	50
Ex connection data	12	Vibrations	50
Low flow cut off	12	Mechanical construction	51
Galvanic isolation	13	Design, dimensions	51
Protocol-specific data	13	Weight	80
Power supply	18	Materials	80
Terminal assignment	18	Process connections	82
Pin assignment, device plug	24	Operability	82
Supply voltage	26	Operating concept	82
Power consumption	26	Local display	83
Current consumption	27	Remote operation	83
Power supply failure	27	Service interface	84
Electrical connection	27	Certificates and approvals	86
Potential equalization	31	CE mark	86
Terminals	31	C-Tick symbol	86
Cable entries	32	Ex approval	87
Cable specification	32	Hygienic compatibility	87
Performance characteristics	33	Certification PROFIBUS	87
Reference operating conditions	33	Modbus RS485 certification	87
Maximum measured error	33	EtherNet/IP certification	87
Repeatability	34	Pressure Equipment Directive	87
Response time	35	Other standards and guidelines	88
Influence of ambient temperature	35	Ordering information	88
Influence of medium temperature	35	Application packages	89
Influence of medium pressure	36	Heartbeat Technology	89
Design fundamentals	36	Concentration	89
Installation	37	Accessories	89
Mounting location	37	Device-specific accessories	89
Orientation	38	Communication-specific accessories	90
Inlet and outlet runs	39	Service-specific accessories	90
Special mounting instructions	39	System components	91
Mounting Safety Barrier Promass 100	41	Documentation	91
Environment	41	Standard documentation	91
Ambient temperature range	41	Supplementary device-dependent documentation	91
Storage temperature	43		
Climate class	43		
Degree of protection	43		

Registered trademarks 92









Document information

Symbols used


Electrical symbols



Symbol	Meaning
	Direct current A terminal to which DC voltage is applied or through which direct current flows.
	Alternating current A terminal to which alternating voltage is applied or through which alternating current flows.
	Direct current and alternating current <ul style="list-style-type: none"> ▪ A terminal to which alternating voltage or DC voltage is applied. ▪ A terminal through which alternating current or direct current flows.
	Ground connection A grounded terminal which, as far as the operator is concerned, is grounded via a grounding system.
	Protective ground connection A terminal which must be connected to ground prior to establishing any other connections.
	Equipotential connection A connection that has to be connected to the plant grounding system: This may be a potential equalization line or a star grounding system depending on national or company codes of practice.

Symbols for certain types of information

Symbol	Meaning
	Permitted Indicates procedures, processes or actions that are permitted.
	Preferred Indicates procedures, processes or actions that are preferred.
	Forbidden Indicates procedures, processes or actions that are forbidden.
	Tip Indicates additional information.
	Reference to documentation Refers to the corresponding device documentation.
	Reference to page Refers to the corresponding page number.
	Reference to graphic Refers to the corresponding graphic number and page number.
	Visual inspection

Symbols in graphics

Symbol	Meaning
1, 2, 3, ...	Item numbers
1., 2., 3. ...	Series of steps
A, B, C, ...	Views
A-A, B-B, C-C, ...	Sections
	Flow direction

Symbol	Meaning
	Hazardous area Indicates a hazardous area.
	Safe area (non-hazardous area) Indicates the non-hazardous area.

Function and system design

Measuring principle

The measuring principle is based on the controlled generation of Coriolis forces. These forces are always present in a system when both translational and rotational movements are superimposed.

$$F_c = 2 \cdot \Delta m (v \cdot \omega)$$

F_c = Coriolis force

Δm = moving mass


ω = rotational velocity

v = radial velocity in rotating or oscillating system

The amplitude of the Coriolis force depends on the moving mass Δm , its velocity v in the system and thus on the mass flow. Instead of a constant rotational velocity ω , the sensor uses oscillation.

In the sensor, an oscillation is produced in the measuring tube. The Coriolis forces produced at the measuring tube cause a phase shift in the tube oscillations (see illustration):

- If there is zero flow (i.e. when the fluid stands still), the oscillation measured at points A and B has the same phase (no phase difference).
- Mass flow causes deceleration of the oscillation at the inlet of the tubes (2) and acceleration at the outlet (3).

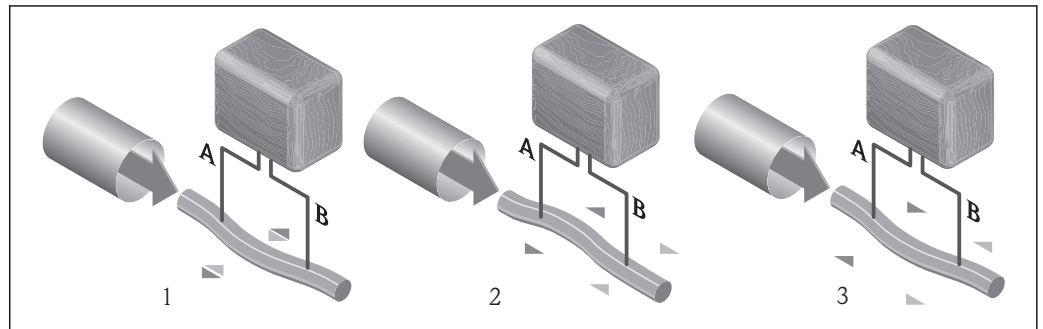

A0016772

The phase difference (A-B) increases with increasing mass flow. Electrodynamical sensors register the tube oscillations at the inlet and outlet. System balance is created by exciting an eccentrically arranged swinging mass to antiphase oscillation. The measuring principle operates independently of temperature, pressure, viscosity, conductivity and flow profile.

Density measurement

The measuring tube is continuously excited at its resonance frequency. A change in the mass and thus the density of the oscillating system (comprising measuring tube and fluid) results in a corresponding, automatic adjustment in the oscillation frequency. Resonance frequency is thus a function of medium density. The microprocessor utilizes this relationship to obtain a density signal.

Volume measurement

Together with the measured mass flow, this is used to calculate the volume flow.

Temperature measurement


The temperature of the measuring tube is determined in order to calculate the compensation factor due to temperature effects. This signal corresponds to the process temperature and is also available as an output signal.

Measuring system


The device consists of a transmitter and a sensor. If a device with Modbus RS485 intrinsically safe is ordered, the Safety Barrier Promass 100 is part of the scope of supply and must be implemented to operate the device.

One device version is available: compact version, transmitter and sensor form a mechanical unit.


Transmitter
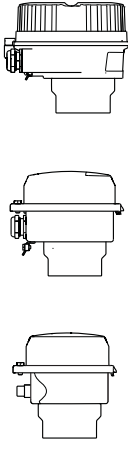
<p>Promass 100</p>  <p>A0016693</p> <p>A0016694</p> <p>A0016695</p>	<p>Device versions and materials:</p> <ul style="list-style-type: none"> ▪ Compact, aluminum coated: Aluminum, AlSi10Mg, coated ▪ Compact, hygienic, stainless: Hygienic version, stainless steel 1.4301 (304) ▪ Ultra compact, hygienic, stainless: Hygienic version, stainless steel 1.4301 (304) <p>Configuration:</p> <ul style="list-style-type: none"> ▪ Via operating tools (e.g. FieldCare) ▪ Additionally for device version with onsite display: Via Web browser (e.g. Microsoft Internet Explorer) ▪ Also for device version with 4-20 mA HART, pulse/frequency/switch output: Via Web browser (e.g. Microsoft Internet Explorer) ▪ Also for device version with EtherNet/IP output: <ul style="list-style-type: none"> - Via Web browser (e.g. Microsoft Internet Explorer) - Via Add-on Profile Level 3 for automation system from Rockwell Automation - Via Electronic Data Sheet (EDS)
---	--


Sensor
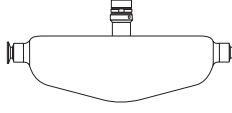
<p>Promass P</p>  <p>A0016509</p>	<ul style="list-style-type: none"> ▪ For use in sterile applications in heavily regulated industries ▪ Nominal diameter range: DN 8 to 50 (3/8 to 2") ▪ Materials: <ul style="list-style-type: none"> - Sensor: stainless steel, 1.4301 (304) - Measuring tubes: stainless steel, 1.4435 (316L) - Process connections: stainless steel, 1.4435 (316L); 1.4404 (316/316L) - Surface quality: Ra_{max}0.76 µm (30 µin) (mechanically polished) Ra_{max}0.38 µm (15 µin) (electropolished) - Delta ferrite <1%
---	---

Safety Barrier Promass 100

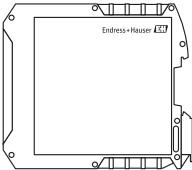 <p>A0016763</p>	<ul style="list-style-type: none"> ▪ Dual-channel safety barrier for installation in non-hazardous locations or Zone 2/Div. 2: <ul style="list-style-type: none"> - Channel 1: DC 24 V power supply - Channel 2: Modbus RS485 ▪ In addition to current, voltage and power limitation, it offers galvanic isolation of circuits for explosion protection. ▪ Easy DIN rail mounting (DIN 35mm) for installation in control cabinets
---	---

Device architecture

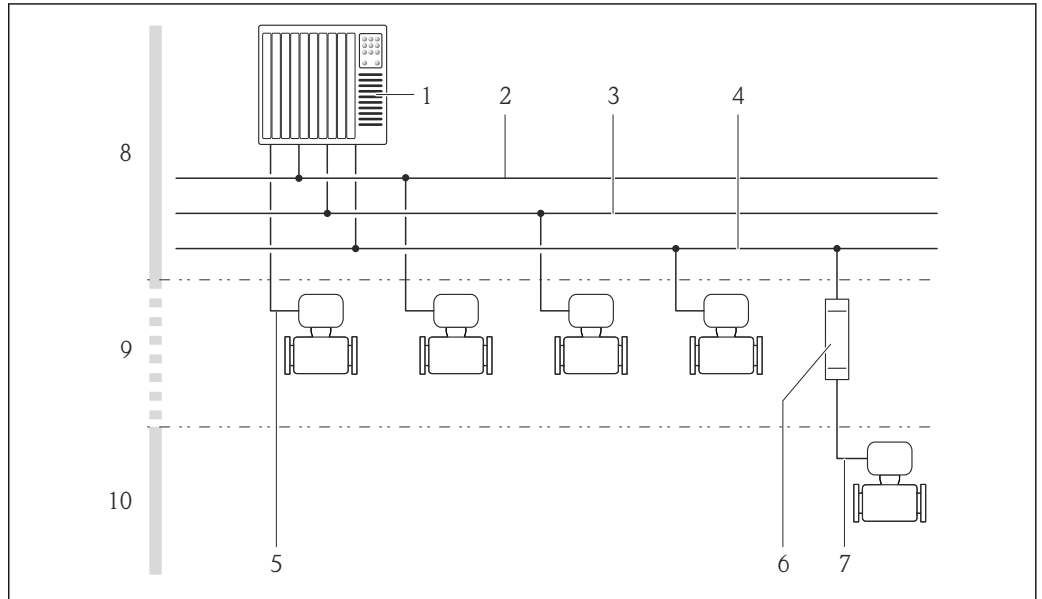

A0016779

1 Possibilities for integrating measuring devices into a system

- 1 Control system (e.g. PLC)
- 2 EtherNet/IP
- 3 PROFIBUS DP
- 4 Modbus RS485
- 5 4-20 mA HART, pulse/frequency/switch output
- 6 Safety Barrier Promass 100
- 7 Modbus RS485 intrinsically safe
- 8 Non-hazardous area
- 9 Non-hazardous area and Zone 2/Div. 2
- 10 Intrinsically safe area and Zone 1/Div. 1

Safety

IT security

We only provide a warranty if the device is installed and used as described in the Operating Instructions. The device is equipped with security mechanisms to protect it against any inadvertent changes to the device settings.

IT security measures in line with operators' security standards and designed to provide additional protection for the device and device data transfer must be implemented by the operators themselves.

Input

Measured variable

Direct measured variables

- Mass flow
- Density
- Temperature

Calculated measured variables

- Volume flow
- Corrected volume flow
- Reference density

Measuring range

Measuring ranges for liquids

DN		Measuring range full scale values $\dot{m}_{\min(F)}$ to $\dot{m}_{\max(F)}$	
[mm]	[in]	[kg/h]	[lb/min]
8	$\frac{3}{8}$	0 to 2 000	0 to 73.50
15	$\frac{1}{2}$	0 to 6 500	0 to 238.9
25	1	0 to 18 000	0 to 661.5
40	$1\frac{1}{2}$	0 to 45 000	0 to 1 654
50	2	0 to 70 000	0 to 2 573


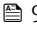
Measuring ranges for gases

The full scale values depend on the density of the gas and can be calculated with the formula below:

$$\dot{m}_{\max(G)} = \dot{m}_{\max(F)} \cdot \rho_G \cdot X$$

$\dot{m}_{\max(G)}$	Maximum full scale value for gas [kg/h]
$\dot{m}_{\max(F)}$	Maximum full scale value for liquid [kg/h]
$\dot{m}_{\max(G)} < \dot{m}_{\max(F)}$	$\dot{m}_{\max(G)}$ can never be greater than $\dot{m}_{\max(F)}$
ρ_G	Gas density in [kg/m ³] at operating conditions

DN		x
[mm]	[in]	[kg/m ³]
8	$\frac{3}{8}$	60
15	$\frac{1}{2}$	80
25	1	90
40	$1\frac{1}{2}$	90
50	2	90

 To calculate the measuring range, use the *Applicator* sizing tool (→  90)

Calculation example for gas

- Sensor: Promass P, DN 50
- Gas: Air with a density of 60.3 kg/m³ (at 20 °C and 50 bar)
- Measuring range (liquid): 70 000 kg/h
- x = 90 kg/m³ (for Promass P, DN 50)

Maximum possible full scale value:

$$\dot{m}_{\max(G)} = \dot{m}_{\max(F)} \cdot \rho_G \cdot X = 70\,000 \text{ kg/h} \cdot 60.3 \text{ kg/m}^3 : 90 \text{ kg/m}^3 = 46\,900 \text{ kg/h}$$

Recommended measuring range

"Flow limit" section (→  48)

Operable flow range


Over 1000 : 1.

Flow rates above the preset full scale value are not overridden by the electronics unit, with the result that the totalizer values are registered correctly.


Output


Output signal

Current output

Current output	4-20 mA HART (active)
Maximum output values	<ul style="list-style-type: none"> ▪ DC 24 V (no flow) ▪ 22.5 mA
Load	0 to 700 Ω
Resolution	0.38 μ A
Damping	Adjustable: 0.07 to 999 s
Assignable measured variables	<ul style="list-style-type: none"> ▪ Mass flow ▪ Volume flow ▪ Corrected volume flow ▪ Density ▪ Reference density ▪ Temperature <p> The range of options increases if the measuring device has one or more application packages.</p>

Pulse/frequency/switch output

Function	Can be set to pulse, frequency or switch output
Version	Passive, open collector
Maximum input values	<ul style="list-style-type: none"> ▪ DC 30 V ▪ 25 mA
Voltage drop	For 25 mA: \leq DC 2 V
Pulse output	
Pulse width	Adjustable: 0.05 to 2 000 ms
Maximum pulse rate	10 000 Impulse/s
Pulse value	Adjustable
Assignable measured variables	<ul style="list-style-type: none"> ▪ Mass flow ▪ Volume flow ▪ Corrected volume flow
Frequency output	
Output frequency	Adjustable: 0 to 10 000 Hz
Damping	Adjustable: 0 to 999 s
Pulse/pause ratio	1:1
Assignable measured variables	<ul style="list-style-type: none"> ▪ Mass flow ▪ Volume flow ▪ Corrected volume flow ▪ Density ▪ Reference density ▪ Temperature <p> The range of options increases if the measuring device has one or more application packages.</p>
Switch output	
Switching behavior	Binary, conductive or non-conductive
Switching delay	Adjustable: 0 to 100 s

Number of switching cycles	Unlimited
Assignable functions	<ul style="list-style-type: none"> ▪ Off ▪ On ▪ Diagnostic behavior ▪ Limit value <ul style="list-style-type: none"> - Mass flow - Volume flow - Corrected volume flow - Density - Reference density - Temperature - Totalizer 1-3 ▪ Flow direction monitoring ▪ Status <ul style="list-style-type: none"> - Partially filled pipe detection - Low flow cut off <p> The range of options increases if the measuring device has one or more application packages.</p>

PROFIBUS DP

Signal encoding	NRZ code
Data transfer	9.6 kBaud...12 MBaud

Modbus RS485

Physical interface	In accordance with EIA/TIA-485-A standard
Terminating resistor	<ul style="list-style-type: none"> ▪ For device version used in non-hazardous areas or Zone 2/Div. 2: integrated and can be activated via DIP switches on the transmitter electronics module ▪ For device version used in intrinsically safe areas: integrated and can be activated via DIP switches on the Safety Barrier Promass 100

EtherNet/IP

Standards	In accordance with IEEE 802.3
------------------	-------------------------------

Signal on alarm

Depending on the interface, failure information is displayed as follows:

Current output

4-20 mA

Failure mode	Selectable (as per NAMUR recommendation NE 43): <ul style="list-style-type: none"> ▪ Minimum value: 3.6 mA ▪ Maximum value: 22 mA ▪ Defined value: 3.59 to 22.5 mA ▪ Actual value ▪ Last valid value
---------------------	---

HART

Device diagnostics	Device condition can be read out via HART Command 48
---------------------------	--

Pulse/frequency/switch output

Pulse output	
Failure mode	Choose from: <ul style="list-style-type: none"> ▪ Actual value ▪ No pulses
Frequency output	
Failure mode	Choose from: <ul style="list-style-type: none"> ▪ Actual value ▪ Defined value: 0 to 12 500 Hz ▪ 0 Hz
Switch output	
Failure mode	Choose from: <ul style="list-style-type: none"> ▪ Current status ▪ Open ▪ Closed

PROFIBUS DP

Status and alarm messages	Diagnostics in accordance with PROFIBUS PA Profile 3.02
---------------------------	---

Modbus RS485

Failure mode	Choose from: <ul style="list-style-type: none"> ▪ NaN value instead of current value ▪ Last valid value
--------------	---

EtherNet/IP

Device diagnostics	Device condition can be read out in Input Assembly
--------------------	--

Local display



Plain text display	With information on cause and remedial measures
Backlight	Red backlighting indicates a device error.

 Status signal as per NAMUR recommendation NE 107

Operating tool

- Via digital communication:
 - HART protocol
 - PROFIBUS DP
 - Modbus RS485
 - EtherNet/IP
- Via service interface

Plain text display	With information on cause and remedial measures
--------------------	---

 Additional information on remote operation (→  83)

Web browser

Plain text display	With information on cause and remedial measures
--------------------	---

Light emitting diodes (LED)

Status information	Status indicated by various light emitting diodes The following information is displayed depending on the device version: <ul style="list-style-type: none"> ■ Supply voltage active ■ Data transmission active ■ Device alarm/error has occurred ■ EtherNet/IP network available ■ EtherNet/IP connection established
---------------------------	--


Ex connection data

These values only apply for the following device version:
Order code for "Output", option **M**: Modbus RS485, for use in intrinsically safe areas


Safety Barrier Promass 100*Safety-related values*

Terminal numbers			
Supply voltage		Signal transmission	
2 (L-)	1 (L+)	26 (A)	27 (B)
$U_{nom} = DC\ 24\ V$ $U_{max} = AC\ 260\ V$		$U_{nom} = DC\ 5\ V$ $U_{max} = AC\ 260\ V$	

Intrinsically safe values

Terminal numbers			
Supply voltage		Signal transmission	
20 (L-)	10 (L+)	62 (A)	72 (B)
$U_o = 16.24\ V$ $I_o = 623\ mA$ $P_o = 2.45\ W$ For IIC*: $L_o = 92.8\ \mu H$, $C_o = 0.433\ \mu F$, $L_o/R_o = 14.6\ \mu H/\Omega$ For IIB*: $L_o = 372\ \mu H$, $C_o = 2.57\ \mu F$, $L_o/R_o = 58.3\ \mu H/\Omega$			
* The gas group depends on the sensor and nominal diameter.			
 For an overview and for information on the interdependencies between the gas group - sensor - nominal diameter, see the "Safety Instructions" (XA) document for the measuring device			

Transmitter*Intrinsically safe values*

Order code for "Approvals"	Terminal numbers			
	Supply voltage		Signal transmission	
	20 (L-)	10 (L+)	62 (A)	72 (B)
<ul style="list-style-type: none"> ■ Option BM: ATEX II2G + IECEx Z1 Ex ia, II2D Ex tb ■ Option BO: ATEX II1/2G + IECEx Z0/Z1 Ex ia, II2D ■ Option BQ: ATEX II1/2G + IECEx Z0/Z1 Ex ia ■ Option BU: ATEX II2G + IECEx Z1 Ex ia ■ Option C2: CSA C/US IS Cl. I, II, III Div. 1 ■ Option 85: ATEX II2G + IECEx Z1 Ex ia + CSA C/US IS Cl. I, II, III Div. 1 	$U_i = 16.24\ V$ $I_i = 623\ mA$ $P_i = 2.45\ W$ $L_i = 0\ \mu H$ $C_i = 6\ nF$			
 For an overview and for information on the interdependencies between the gas group - sensor - nominal diameter, see the "Safety Instructions" (XA) document for the measuring device				

Low flow cut off

The switch points for low flow cut off are user-selectable.


Galvanic isolation

The following connections are galvanically isolated from each other:

- Outputs
- Power supply

Protocol-specific data

HART



Manufacturer ID	0x11
Device type ID	0x4A
HART protocol revision	7.0
Device description files (DTM, DD)	Information and files under: www.endress.com
HART load	Min. 250 Ω
Dynamic variables	<p>Read out the dynamic variables: HART command 3 The measured variables can be freely assigned to the dynamic variables.</p> <p>Measured variables for PV (primary dynamic variable)</p> <ul style="list-style-type: none"> ▪ Mass flow ▪ Volume flow ▪ Corrected volume flow ▪ Density ▪ Reference density ▪ Temperature <p>Measured variables for SV, TV, QV (secondary, tertiary and quaternary dynamic variable)</p> <ul style="list-style-type: none"> ▪ Mass flow ▪ Volume flow ▪ Corrected volume flow ▪ Density ▪ Reference density ▪ Temperature ▪ Totalizer 1 ▪ Totalizer 2 ▪ Totalizer 3 <p> The range of options increases if the measuring device has one or more application packages.</p> <p>Heartbeat Technology Application Package Additional measured variables are available with the Heartbeat Technology application package:</p> <ul style="list-style-type: none"> ▪ Carrier pipe temperature ▪ Oscillation amplitude 0
Device variables	<p>Read out the device variables: HART command 9 The device variables are permanently assigned.</p> <p>A maximum of 8 device variables can be transmitted:</p> <ul style="list-style-type: none"> ▪ 0 = mass flow ▪ 1 = volume flow ▪ 2 = corrected volume flow ▪ 3 = density ▪ 4 = reference density ▪ 5 = temperature ▪ 6 = totalizer 1 ▪ 7 = totalizer 2 ▪ 8 = totalizer 3 ▪ 9 = dynamic viscosity ▪ 10 = kinematic viscosity ▪ 11 = temp. compensated dynamic viscosity ▪ 12 = temp. compensated kinematic viscosity ▪ 13 = target mass flow ▪ 14 = carrier mass flow ▪ 15 = concentration

PROFIBUS DP

Manufacturer ID	0x11
Ident number	0x1561
Profile version	3.02
Device description files (GSD, DTM, DD)	Information and files under: <ul style="list-style-type: none"> ▪ www.endress.com ▪ www.profibus.org
Output values (from measuring device to automation system)	<p>Analog input 1 to 8</p> <ul style="list-style-type: none"> ▪ Mass flow ▪ Volume flow ▪ Corrected volume flow ▪ Target mass flow ▪ Carrier mass flow ▪ Density ▪ Reference density ▪ Concentration ▪ Dynamic viscosity ▪ Kinematic viscosity ▪ Temp. compensated dynamic viscosity ▪ Temp. compensated kinematic viscosity ▪ Temperature ▪ Carrier pipe temperature ▪ Electronic temperature ▪ Oscillation frequency ▪ Oscillation amplitude ▪ Frequency fluctuation ▪ Oscillation damping ▪ Tube damping fluctuation ▪ Signal asymmetry ▪ Exciter current <p>Digital input 1 to 2</p> <ul style="list-style-type: none"> ▪ Partially filled pipe detection ▪ Low flow cut off <p>Totalizer 1 to 3</p> <ul style="list-style-type: none"> ▪ Mass flow ▪ Volume flow ▪ Corrected volume flow
Input values (from automation system to measuring device)	<p>Analog output 1 to 3 (fixed assignment)</p> <ul style="list-style-type: none"> ▪ Pressure ▪ Temperature ▪ Reference density <p>Digital output 1 to 3 (fixed assignment)</p> <ul style="list-style-type: none"> ▪ Digital output 1: switch positive zero return on/off ▪ Digital output 2: perform zero point adjustment ▪ Digital output 3: switch switch output on/off <p>Totalizer 1 to 3</p> <ul style="list-style-type: none"> ▪ Totalize ▪ Reset and hold ▪ Preset and hold ▪ Stop ▪ Operating mode configuration: <ul style="list-style-type: none"> – Net flow total – Forward flow total – Reverse flow total

Supported functions	<ul style="list-style-type: none"> ▪ Identification & Maintenance Simplest device identification on the part of the control system and nameplate ▪ PROFIBUS upload/download Reading and writing parameters is up to ten times faster with PROFIBUS upload/download ▪ Condensed status Simplest and self-explanatory diagnostic information by categorizing diagnostic messages that occur
Configuration of the device address	<ul style="list-style-type: none"> ▪ DIP switches on the I/O electronics module ▪ Via operating tools (e.g. FieldCare)


Modbus RS485

Protocol	Modbus Applications Protocol Specification V1.1
Device type	Slave
Slave address range	1 to 247
Broadcast address range	0
Function codes	<ul style="list-style-type: none"> ▪ 03: Read holding register ▪ 04: Read input register ▪ 06: Write single registers ▪ 08: Diagnostics ▪ 16: Write multiple registers ▪ 23: Read/write multiple registers
Broadcast messages	Supported by the following function codes: <ul style="list-style-type: none"> ▪ 06: Write single registers ▪ 16: Write multiple registers ▪ 23: Read/write multiple registers
Supported baud rate	<ul style="list-style-type: none"> ▪ 1 200 BAUD ▪ 2 400 BAUD ▪ 4 800 BAUD ▪ 9 600 BAUD ▪ 19 200 BAUD ▪ 38 400 BAUD ▪ 57 600 BAUD ▪ 115 200 BAUD
Data transfer mode	<ul style="list-style-type: none"> ▪ ASCII ▪ RTU
Data access	Each device parameter can be accessed via Modbus RS485.  For Modbus register information (→  91)

EtherNet/IP

Protocol	<ul style="list-style-type: none"> ▪ The CIP Networks Library Volume 1: Common Industrial Protocol ▪ The CIP Networks Library Volume 2: EtherNet/IP Adaptation of CIP
Communication type	<ul style="list-style-type: none"> ▪ 10Base-T ▪ 100Base-TX
Device profile	Generic device (product type: 0x2B)
Manufacturer ID	0x49E
Device type ID	0x104A
Baud rates	Automatic ¹⁰ / ₁₀₀ Mbit with half-duplex and full-duplex detection
Polarity	Auto-polarity for automatic correction of crossed TxD and RxD pairs
Supported CIP connections	Max. 3 connections
Explicit connections	Max. 6 connections


I/O connections	Max. 6 connections (scanner)		
Configuration options for measuring device	<ul style="list-style-type: none"> ▪ DIP switches on the electronics module for IP addressing ▪ Manufacturer-specific software (FieldCare) ▪ Add-on Profile Level 3 for Rockwell Automation control systems ▪ Web browser ▪ Electronic Data Sheet (EDS) integrated in the measuring device 		
Configuration of the EtherNet interface	<ul style="list-style-type: none"> ▪ Speed: 10 MBit, 100 MBit, auto (factory setting) ▪ Duplex: half-duplex, full-duplex, auto (factory setting) 		
Configuration of the device address	<ul style="list-style-type: none"> ▪ DIP switches on the electronics module for IP addressing (last octet) ▪ DHCP ▪ Manufacturer-specific software (FieldCare) ▪ Add-on Profile Level 3 for Rockwell Automation control systems ▪ Web browser ▪ EtherNet/IP tools, e.g. RSLinx (Rockwell Automation) 		
Device Level Ring (DLR)	No		
Fix Input			
RPI	5 ms to 10 s (factory setting: 20 ms)		
Exclusive Owner Multicast		Instance	Size [byte]
	Instance configuration:	0x68	398
	O → T configuration:	0x66	64
	T → O configuration:	0x64	44
Exclusive Owner Multicast		Instance	Size [byte]
	Instance configuration:	0x69	-
	O → T configuration:	0x66	64
	T → O configuration:	0x64	44
Input only Multicast		Instance	Size [byte]
	Instance configuration:	0x68	398
	O → T configuration:	0xC7	-
	T → O configuration:	0x64	44
Input only Multicast		Instance	Size [byte]
	Instance configuration:	0x69	-
	O → T configuration:	0xC7	-
	T → O configuration:	0x64	44
Input Assembly	<ul style="list-style-type: none"> ▪ Current device diagnostics ▪ Mass flow ▪ Volume flow ▪ Corrected volume flow ▪ Density ▪ Reference density ▪ Temperature ▪ Totalizer 1 ▪ Totalizer 2 ▪ Totalizer 3 		
Configurable Input			
RPI	5 ms to 10 s (factory setting: 20 ms)		
Exclusive Owner Multicast		Instance	Size [byte]
	Instance configuration:	0x68	398
	O → T configuration:	0x66	64
	T → O configuration:	0x65	88
Exclusive Owner Multicast		Instance	Size [byte]
	Instance configuration:	0x69	-

	O → T configuration:	0x66	64
	T → O configuration:	0x65	88
Input only Multicast		Instance	Size [byte]
	Instance configuration:	0x68	398
	O → T configuration:	0xC7	-
	T → O configuration:	0x65	88
Input only Multicast		Instance	Size [byte]
	Instance configuration:	0x69	-
	O → T configuration:	0xC7	-
	T → O configuration:	0x65	88
Configurable Input Assembly	<ul style="list-style-type: none"> ■ Current device diagnostics ■ Mass flow ■ Volume flow ■ Corrected volume flow ■ Density ■ Reference density ■ Temperature ■ Totalizer 1 ■ Totalizer 2 ■ Totalizer 3 <p> The range of options increases if the measuring device has one or more application packages.</p>		
Fix Output			
Output Assembly	<ul style="list-style-type: none"> ■ Activation of reset totalizers 1-3 ■ Activation of pressure compensation ■ Activation of reference density compensation ■ Activation of temperature compensation ■ Reset totalizers 1-3 ■ External pressure value ■ Pressure unit ■ External reference density ■ Reference density unit ■ External temperature ■ Temperature unit 		
Configuration			
Configuration Assembly	<p>Only the most common configurations are listed below.</p> <ul style="list-style-type: none"> ■ Software write protection ■ Mass flow unit ■ Mass unit ■ Volume flow unit ■ Volume unit ■ Corrected volume flow unit ■ Corrected volume unit ■ Density unit ■ Reference density unit ■ Temperature unit ■ Pressure unit ■ Length ■ Totalizer 1-3: <ul style="list-style-type: none"> - Assignment - Unit - Measuring mode - Failure mode ■ Alarm delay 		

Power supply

Terminal assignment

Overview: housing version

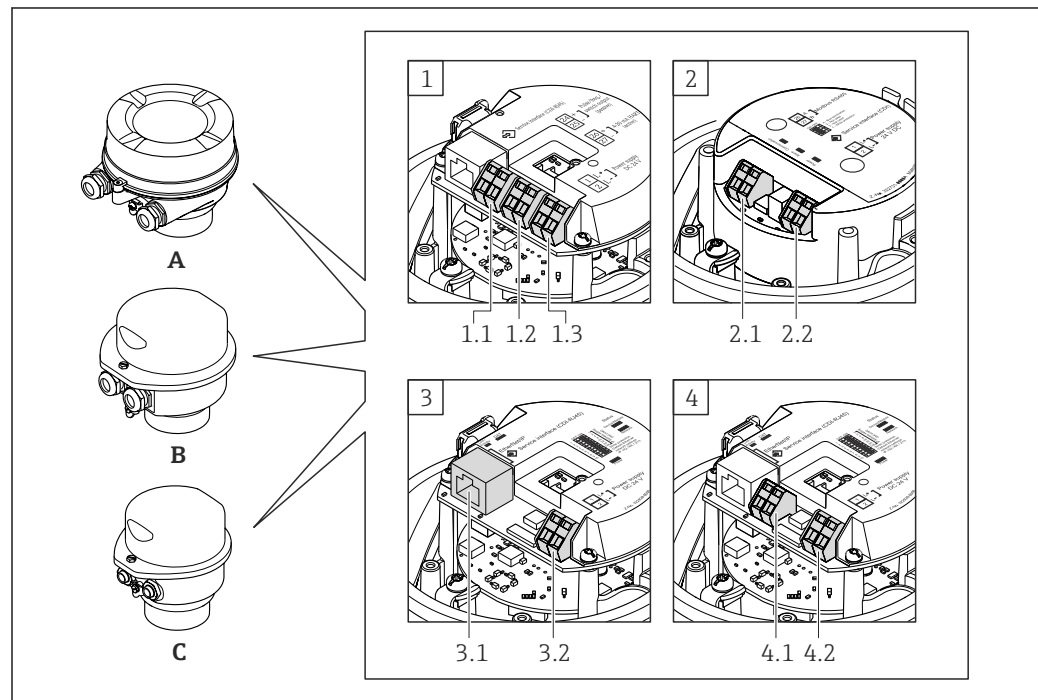

A0016770

- A Housing version: compact, aluminum coated
- B Housing version: compact, hygienic, stainless
- C Housing version: ultra compact, hygienic, stainless, M12 device plug
- 1 Connection version: 4-20 mA HART, pulse/frequency/switch output
 - 1.1 Signal transmission: pulse/frequency/switch output
 - 1.2 Signal transmission: 4-20 mA HART
 - 1.3 Supply voltage
- 2 Connection version: Modbus RS485
 - 2.1 Signal transmission
 - 2.2 Supply voltage
- 3 Connection version: EtherNet/IP
 - 3.1 Signal transmission
 - 3.2 Supply voltage
- 4 Connection version: PROFIBUS DP
 - 4.1 PROFIBUS DP
 - 4.2 Supply voltage

Transmitter

Connection version 4-20 mA HART with pulse/frequency/switch output


Order code for "Output", option **B**

Depending on the housing version, the transmitters can be ordered with terminals or device plugs.

Order code for "Housing"	Connection methods available		Possible options for order code "Electrical connection"
	Outputs	Power supply	
Options A, B	Terminals	Terminals	<ul style="list-style-type: none"> ▪ Option A: coupling M20x1 ▪ Option B: thread M20x1 ▪ Option C: thread G ½" ▪ Option D: thread NPT ½"
Options A, B	Device plug	Terminals	<ul style="list-style-type: none"> ▪ Option L: plug M12x1 + thread NPT ½" ▪ Option N: plug M12x1 + coupling M20 ▪ Option P: plug M12x1 + thread G ½" ▪ Option U: plug M12x1 + thread M20
Options A, B, C	Device plug	Device plug	Option Q: 2 x plug M12x1

Order code for "Housing":

- Option A: compact, coated aluminum
- Option B: compact, hygienic, stainless
- Option C: ultra compact, hygienic, stainless, M12 device plug

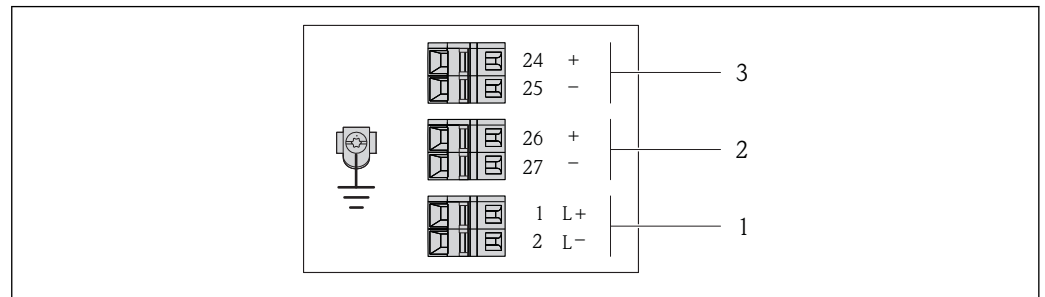

A0016888

2 Terminal assignment 4-20 mA HART with pulse/frequency/switch output

- 1 Power supply: DC 24 V
- 2 Output 1: 4-20 mA HART (active)
- 3 Output 2: pulse/frequency/switch output (passive)

Order code for "Output"	Terminal number					
	Power supply		Output 1		Output 2	
	2 (L-)	1 (L+)	27 (-)	26 (+)	25 (-)	24 (+)
Option B	DC 24 V		4-20 mA HART (active)		Pulse/frequency/switch output (passive)	

Order code for "Output":

Option B: 4-20 mA HART with pulse/frequency/switch output

PROFIBUS DP connection version

 For use in the non-hazardous area and Zone 2/Div. 2.


Order code for "Output", option **L**

Depending on the housing version, the transmitters can be ordered with terminals or device plugs.

Order code for "Housing"	Connection methods available		Possible options for order code "Electrical connection"
	Output	Power supply	
Options A, B	Terminals	Terminals	<ul style="list-style-type: none"> ▪ Option A: coupling M20x1 ▪ Option B: thread M20x1 ▪ Option C: thread G ½" ▪ Option D: thread NPT ½"
Options A, B	Device plug	Terminals	<ul style="list-style-type: none"> ▪ Option L: plug M12x1 + thread NPT ½" ▪ Option N: plug M12x1 + coupling M20 ▪ Option P: plug M12x1 + thread G ½" ▪ Option U: plug M12x1 + thread M20
Options A, B, C	Device plug	Device plug	Option Q: 2 x plug M12x1

Order code for "Housing":

- Option A: compact, coated aluminum
- Option A: compact, hygienic, stainless
- Option C: ultra compact, hygienic, stainless, M12 device plug

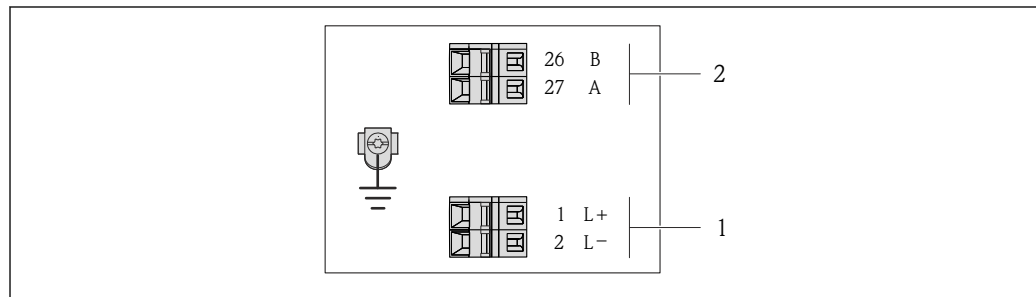

A0022716

 3 *PROFIBUS DP terminal assignment*

- 1 Power supply: DC 24 V
- 2 PROFIBUS DP

Order code for "Output"	Terminal number			
	Power supply		Output	
	2 (L-)	1 (L+)	26 (Rx/D/TxD-P)	27 (Rx/D/TxD-N)
Option L	DC 24 V		B	A

Order code for "Output":
Option L: PROFIBUS DP, for use in non-hazardous areas and Zone 2/div. 2

Modbus RS485 connection version

 For use in the non-hazardous area and Zone 2/Div. 2.


Order code for "Output", option **M**

Depending on the housing version, the transmitters can be ordered with terminals or device plugs.


Order code for "Housing"	Connection methods available		Possible options for order code "Electrical connection"
	Output	Power supply	
Options A, B	Terminals	Terminals	<ul style="list-style-type: none"> ▪ Option A: coupling M20x1 ▪ Option B: thread M20x1 ▪ Option C: thread G ½" ▪ Option D: thread NPT ½"
Options A, B	Device plug	Terminals	<ul style="list-style-type: none"> ▪ Option L: plug M12x1 + thread NPT ½" ▪ Option N: plug M12x1 + coupling M20 ▪ Option P: plug M12x1 + thread G ½" ▪ Option U: plug M12x1 + thread M20
Options A, B, C	Device plug	Device plug	Option Q: 2 x plug M12x1

Order code for "Housing":

- Option A: compact, coated aluminum
- Option B: compact, hygienic, stainless
- Option C: ultra compact, hygienic, stainless, M12 device plug
- Option C: ultra compact, stainless, M12 device plug

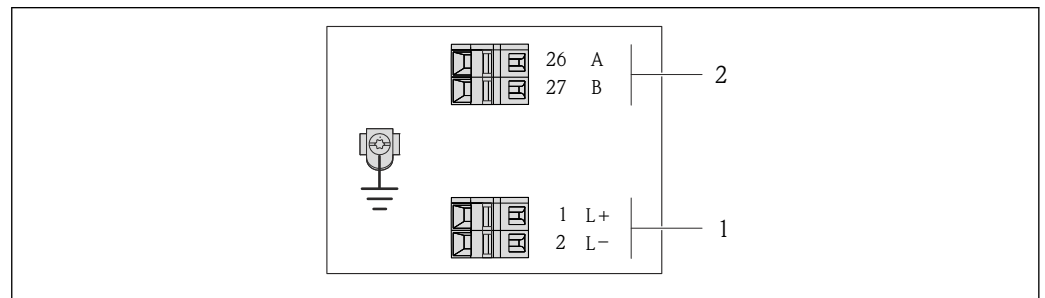

A0019528

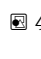
 4 Modbus RS485 terminal assignment, connection version for use in non-hazardous areas and Zone 2/Div. 2

- 1 Power supply: DC 24 V
- 2 Modbus RS485

Order code for "Output"	Terminal number			
	Power supply		Output	
	2 (L-)	1 (L+)	27 (B)	26 (A)
Option M	DC 24 V		Modbus RS485	

Order code for "Output":
Option **M**: Modbus RS485, for use in non-hazardous areas and Zone 2/Div. 2


Modbus RS485 connection version

 For use in the intrinsically safe area. Connection via Safety Barrier Promass 100.


Order code for "Output", option **M**

Depending on the housing version, the transmitters can be ordered with terminals or device plugs.

Order code for "Housing"	Connection methods available		Possible options for order code "Electrical connection"
	Output	Power supply	
Options A, B	Terminals	Terminals	<ul style="list-style-type: none"> ■ Option A: coupling M20x1 ■ Option B: thread M20x1 ■ Option C: thread G ½" ■ Option D: thread NPT ½"
A, B, C	Device plug		Option I: plug M12x1
Order code for "Housing": <ul style="list-style-type: none"> ■ Option A: compact, coated aluminum ■ Option B: compact, hygienic, stainless ■ Option C: ultra compact, hygienic, stainless, M12 device plug 			

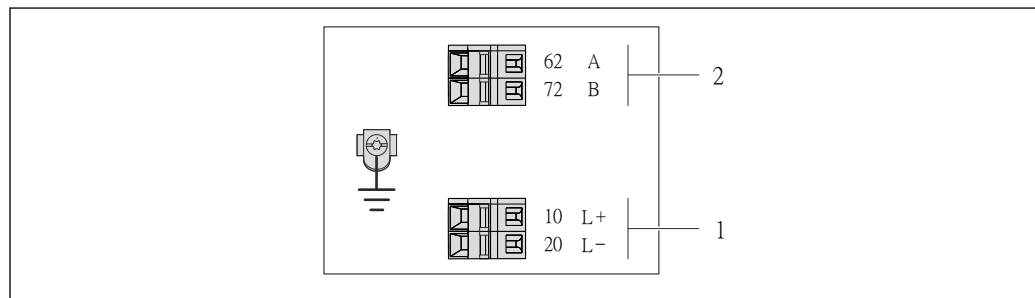

A0017053

 5 *Modbus RS485 terminal assignment, connection version for use in intrinsically safe areas (connection via Safety Barrier Promass 100)*

- 1 *Intrinsically safe power supply*
- 2 *Modbus RS485*

Order code for "Output"	20 (L-)	10 (L+)	72 (B)	62 (A)
Option M	Intrinsically safe supply voltage		Modbus RS485 intrinsically safe	
Order code for "Output": Option M : Modbus RS485, for use in intrinsically safe areas (connection via Safety Barrier Promass 100)				

EtherNet/IP connection version


Order code for "Output", option **N**

Depending on the housing version, the transmitters can be ordered with terminals or device plugs.

Order code for "Housing"	Connection methods available		Possible options for order code "Electrical connection"
	Output	Power supply	
Options A, B	Device plug	Terminals	<ul style="list-style-type: none"> ▪ Option L: plug M12x1 + thread NPT 1/2" ▪ Option N: plug M12x1 + coupling M20 ▪ Option P: plug M12x1 + thread G 1/2" ▪ Option U: plug M12x1 + thread M20
Options A, B, C	Device plug	Device plug	Option Q : 2 x plug M12x1

Order code for "Housing":

- Option **A**: compact, coated aluminum
- Option **B**: compact, hygienic, stainless
- Option **C**: ultra compact, hygienic, stainless, M12 device plug

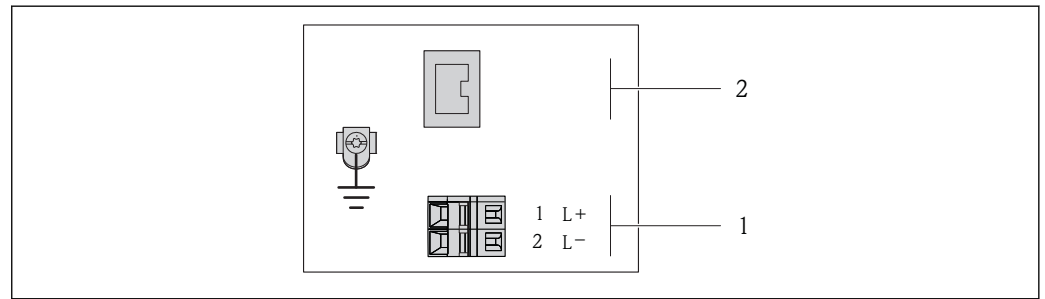


6 EtherNet/IP terminal assignment

- 1 Power supply: DC 24 V
- 2 EtherNet/IP

Order code for "Output"	Terminal number		Output Device plug M12x1
	Power supply 2 (L-)	1 (L+)	
Option N	DC 24 V		EtherNet/IP

Order code for "Output":
Option **N**: EtherNet/IP

Safety Barrier Promass 100

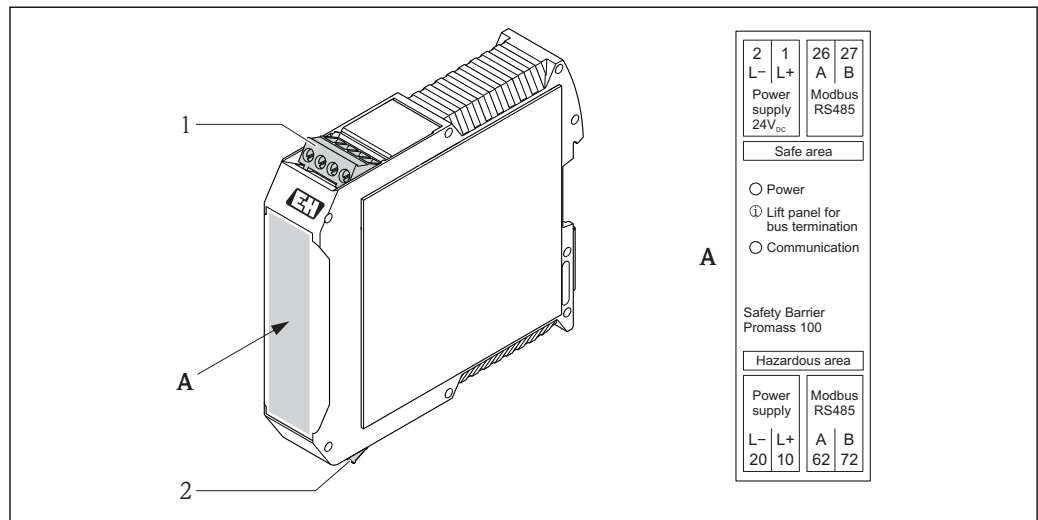

7 Safety Barrier Promass 100 with terminals

- 1 Non-hazardous area and Zone 2/Div. 2
- 2 Intrinsically safe area

Pin assignment, device plug

- i** Order codes for the M12x1 connectors, see the "Order code for **electrical connection**" column:
 - 4-20 mA HART, pulse/frequency/switch output (→ 18)
 - PROFIBUS-DP (→ 20)
 - Modbus RS485 (→ 21)
 - EtherNet/IP (→ 23)

Supply voltage

For all connection versions except MODBUS RS485 intrinsically safe (device side)


- i** Device plug MODBUS RS485 intrinsically safe with supply voltage (→ 25)

	Pin	Assignment	Coding	Plug/socket
	1	L+	DC 24 V	A
2				
3				
4	L-	DC 24 V		
5		Grounding/shielding		

- i** The following is recommended as a socket:
 - Binder, series 763, part no. 79 3440 35 05
 - Alternatively: Phoenix part no. 1669767 SAC-5P-M12MS
 - With the order code for "Output", option **B**: 4-20 mA HART, pulse/frequency/switch output
 - With the order code for "Output", option **N**: EtherNet/IP
 - When using the device in a hazardous location: Use a suitably certified socket.

4-20 mA HART with pulse/frequency/switch output

Device plug for signal transmission (device side)

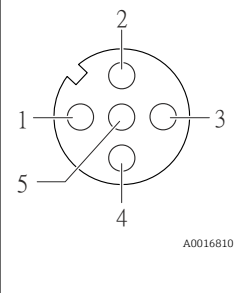
	Pin	Assignment	Coding	Plug/socket	
	1	+	4-20 mA HART (active)	A	Socket
	2	-	4-20 mA HART (active)		
	3	+	Pulse/frequency/switch output (passive)		
	4	-	Pulse/frequency/switch output (passive)		
	5		Grounding/shielding		



-  Recommended plug: Binder, series 763, part no. 79 3439 12 05
-  When using the device in a hazardous location: Use a suitably certified plug.

PROFIBUS DP

-  For use in the non-hazardous area and Zone 2/Div. 2.

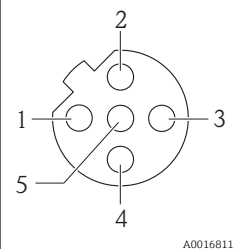
Device plug for signal transmission (device side)



	Pin	Assignment	Coding	Plug/socket	
	1			B	Socket
	2	A	PROFIBUS DP		
	3				
	4	B	PROFIBUS DP		
	5		Grounding/shielding		

-  Recommended plug: Binder, series 763, part no. 79 4449 20 05
-  When using the device in a hazardous location: Use a suitably certified plug.

MODBUS RS485

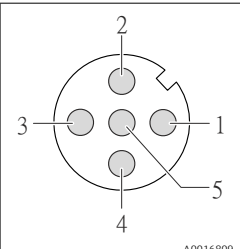
Device plug for signal transmission with supply voltage (device side), MODBUS RS485 (intrinsically safe)

	Pin	Assignment	Coding	Plug/socket	
	1	L+	Supply voltage, intrinsically safe	A	Plug
	2	A	Modbus RS485 intrinsically safe		
	3	B			
	4	L-	Supply voltage, intrinsically safe		
	5		Grounding/shielding		

-  Recommended socket: Binder, series 763, part no. 79 3439 12 05
-  When using the device in a hazardous location: Use a suitably certified socket.

Device plug for signal transmission (device side), MODBUS RS485 (not intrinsically safe)


i For use in the non-hazardous area and Zone 2/Div. 2.

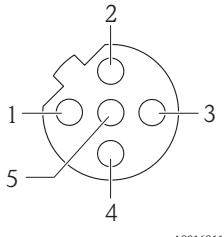 <p>A0016811</p>	Pin	Assignment		Coding	Plug/socket
	1			B	Socket
	2	A	Modbus RS485		
	3				
	4	B	Modbus RS485		
5		Grounding/shielding			

- i**
- Recommended plug: Binder, series 763, part no. 79 4449 20 05
 - When using the device in a hazardous location: Use a suitably certified plug.

EtherNet/IP

Device plug for signal transmission (device side)

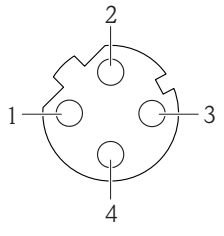 <p>A0016812</p>	Pin	Assignment		Coding	Plug/socket
	1	+	Tx	D	Socket
	2	+	Rx		
	3	-	Tx		
4	-	Rx			

- i** Recommended plug:
- Binder, series 763, part no. 99 3729 810 04
 - Phoenix, part no. 1543223 SACC-M12MSD-4Q
 - When using the device in a hazardous location: Use a suitably certified plug.

Supply voltage

Transmitter

- For device version with all communication types except Modbus RS485 intrinsically safe: DC 20 to 30 V
- For device version with Modbus RS485 intrinsically safe: power supply via Safety Barrier Promass 100

The power unit must be tested to ensure it meets safety requirements (e.g. PELV, SELV).

Safety Barrier Promass 100

DC 20 to 30 V

Power consumption

Transmitter

Order code for "Output"	Maximum Power consumption
Option B : 4-20mA HART, pulse/frequency/switch output	3.5 W
Option L : PROFIBUS DP	3.5 W
Option M : Modbus RS485, for use in non-hazardous areas and Zone 2/Div. 2	3.5 W
Option M : Modbus RS485, for use in intrinsically safe areas	2.45 W
Option N : EtherNet/IP	3.5 W

Safety Barrier Promass 100

Order code for "Output"	Maximum Power consumption
Option M : Modbus RS485, for use in intrinsically safe areas	4.8 W

Current consumption

Transmitter

Order code for "Output"	Maximum Current consumption	Maximum switch-on current
Option B : 4-20mA HART, pul./freq./switch output	145 mA	18 A (<0.125 ms)
Option L : PROFIBUS DP	145 mA	18 A (<0.125 ms)
Option M : Modbus RS485, for use in non-hazardous areas and Zone 2/Div. 2	90 mA	10 A (<0.8 ms)
Option M : Modbus RS485, for use in intrinsically safe areas	145 mA	16 A (<0.4 ms)
Option N : EtherNet/IP	145 mA	18 A (<0.125 ms)

Safety Barrier Promass 100


Order code for "Output"	Maximum Current consumption	Maximum switch-on current
Option M : Modbus RS485, for use in intrinsically safe areas	230 mA	10 A (<0.8 ms)

Power supply failure

- Totalizers stop at the last value measured.
- Depending on the device version, the configuration is retained in the device memory or in the plug-in memory (HistoROM DAT).
- Error messages (incl. total operated hours) are stored.

Electrical connection

Connecting the transmitter

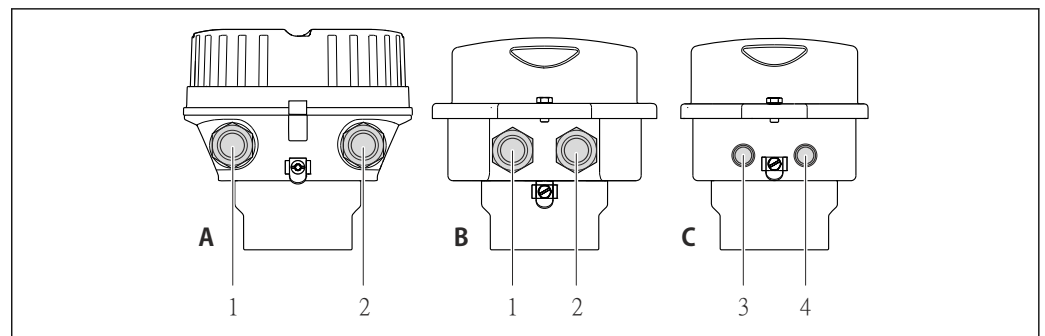


- A Housing version: compact, aluminum coated
- B Housing version: compact hygienic, stainless
- 1 Cable entry or device plug for signal transmission
- 2 Cable entry or device plug for supply voltage
- C Housing version: ultra-compact hygienic, stainless, M12 device plug
- 3 Device plug for signal transmission
- 4 Device plug for supply voltage

- Terminal assignment (→ 18)
- Pin assignment, device plug (→ 24)

In the case of device versions with a connector, the transmitter housing does not need to be opened to connect the signal cable or power supply cable.

Connection examples

Current output 4-20 mA HART

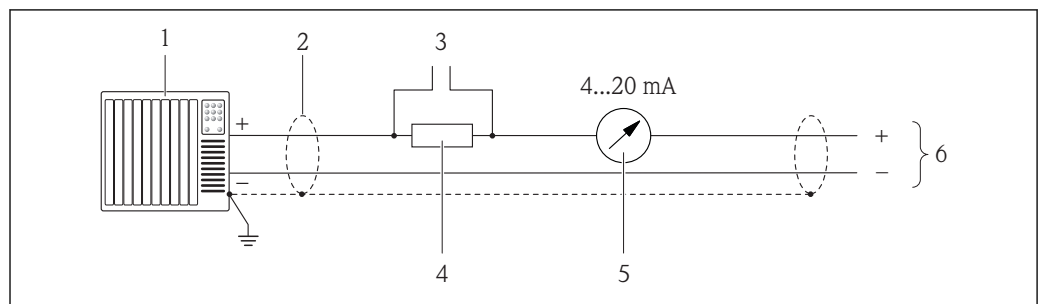


A0016800

8 Connection example for 4-20 mA HART current output (active)

- 1 Automation system with current input (e.g. PLC)
- 2 Cable shield, observe cable specifications (→ 32)
- 3 Connection for HART operating devices (→ 83)
- 4 Resistor for HART communication ($\geq 250 \Omega$): observe maximum load (→ 9)
- 5 Analog display unit: observe maximum load (→ 9)
- 6 Transmitter

Pulse/frequency output

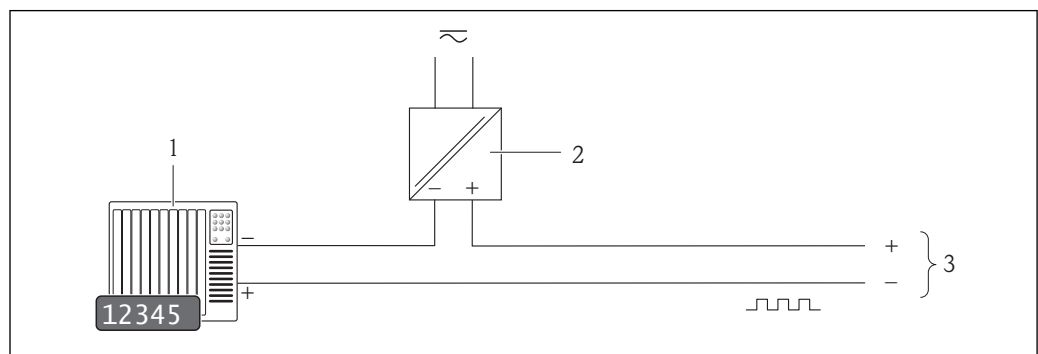


A0016801

9 Connection example for pulse/frequency output (passive)

- 1 Automation system with pulse/frequency input (e.g. PLC)
- 2 Power supply
- 3 Transmitter: observe input values (→ 9)

Switch output

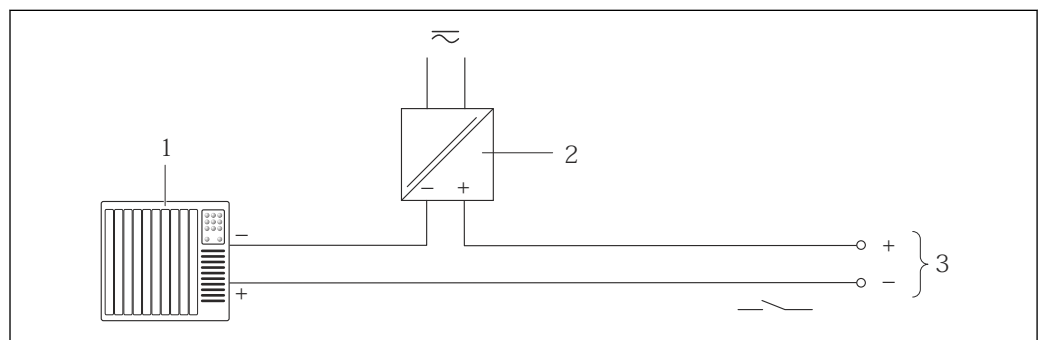


A0016802

10 Connection example for switch output (passive)

- 1 Automation system with switch input (e.g. PLC)
- 2 Power supply
- 3 Transmitter: observe input values (→ 9)

PROFIBUS DP

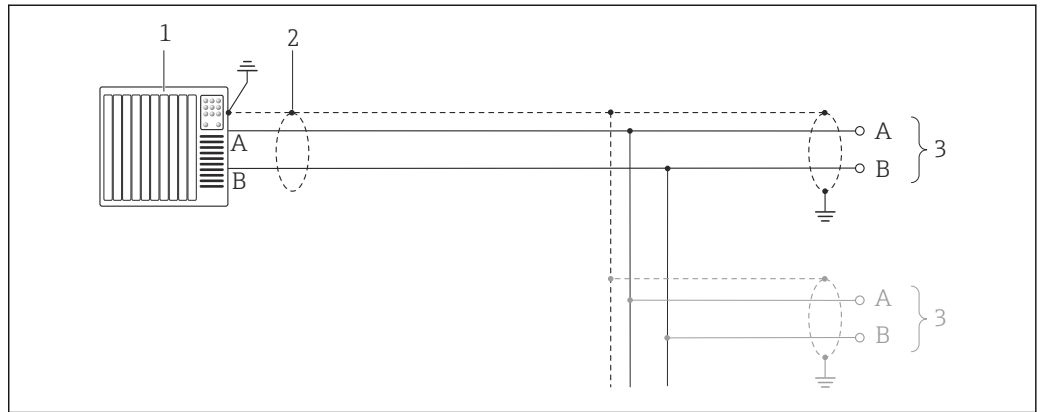


A0021429

11 Connection example for PROFIBUS DP, non-hazardous area and Zone 2/Div. 2

- 1 Control system (e.g. PLC)
- 2 Cable shield: the cable shield must be grounded at both ends to comply with EMC requirements; observe cable specifications (→ 32)
- 3 Distribution box
- 4 Transmitter

i If baud rates > 1.5 MBaud an EMC cable entry must be used and the cable shield must continue as far as the terminal wherever possible.


Modbus RS485

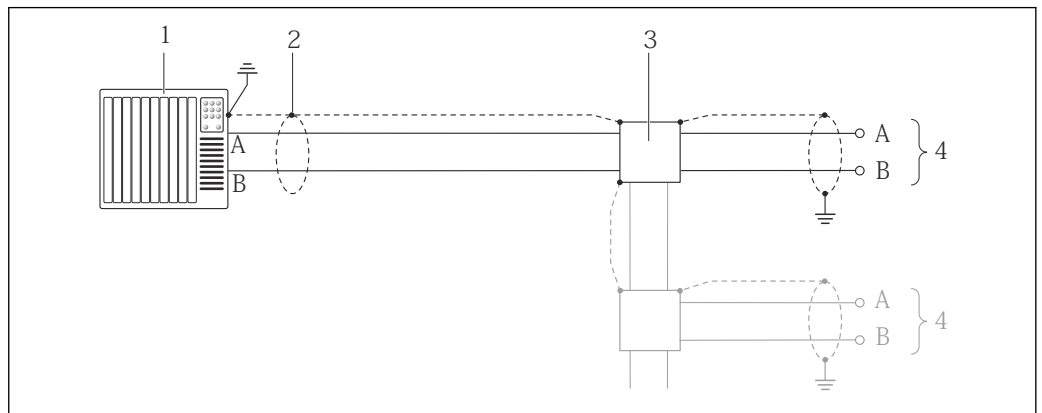

A0016803

12 Connection example for Modbus RS485, non-hazardous area and Zone 2/Div. 2

- 1 Control system (e.g. PLC)
- 2 Cable shield: the cable shield must be grounded at both ends to comply with EMC requirements; observe cable specifications (→ 32)
- 3 Distribution box
- 4 Transmitter

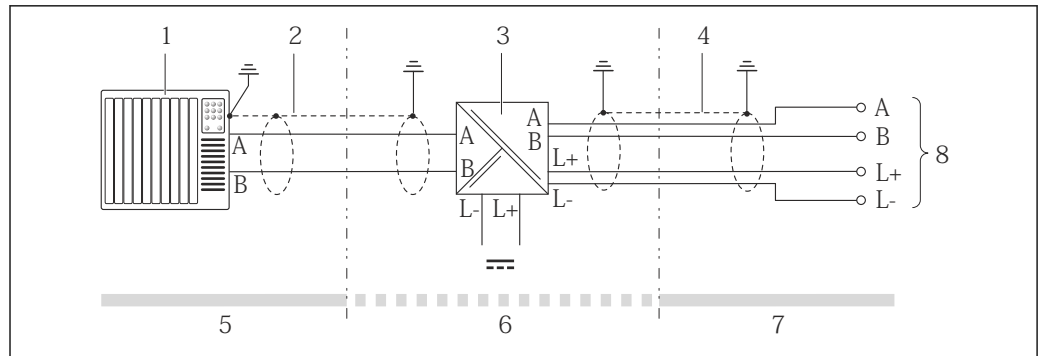


A0016804

13 Connection example for Modbus RS485 intrinsically safe

- 1 Control system (e.g. PLC)
- 2 Cable shield, observe cable specifications (→ 32)
- 3 Safety Barrier Promass 100
- 4 Observe cable specifications (→ 32)
- 5 Non-hazardous area
- 6 Non-hazardous area and Zone 2/Div. 2
- 7 Intrinsically safe area
- 8 Transmitter

EtherNet/IP

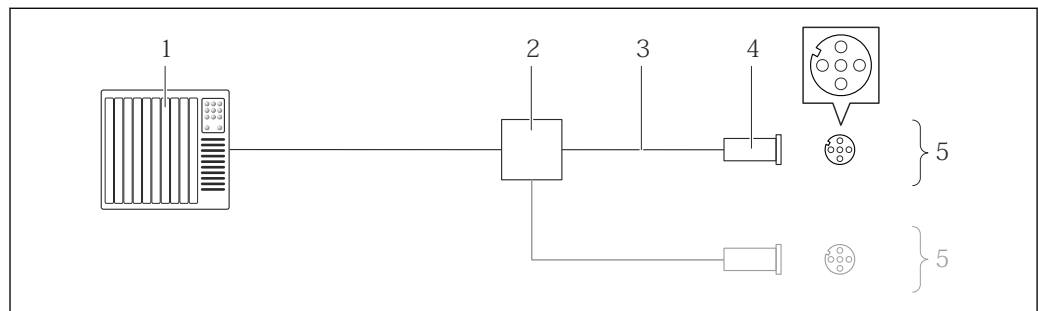


A0016805

14 Connection example for EtherNet/IP


- 1 Control system (e.g. PLC)
- 2 Ethernet switch
- 3 Observe cable specifications (→ 32)
- 4 Connector
- 5 Transmitter

HART input

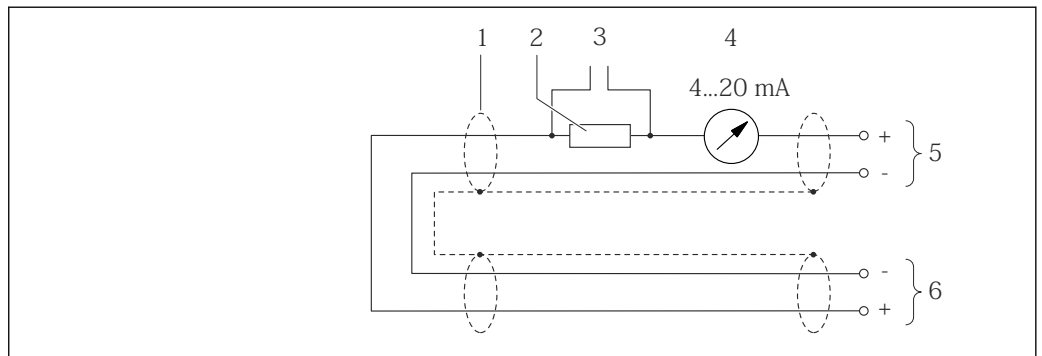

15 Connection example for HART input (burst mode) via current output (active)

- 1 Cable shield, observe cable specifications (→ 32)
- 2 Resistor for HART communication ($\geq 250 \Omega$): observe maximum load (→ 9)
- 3 Connection for HART operating devices (→ 83)
- 4 Analog display unit
- 5 Transmitter
- 6 Sensor for external measured variable

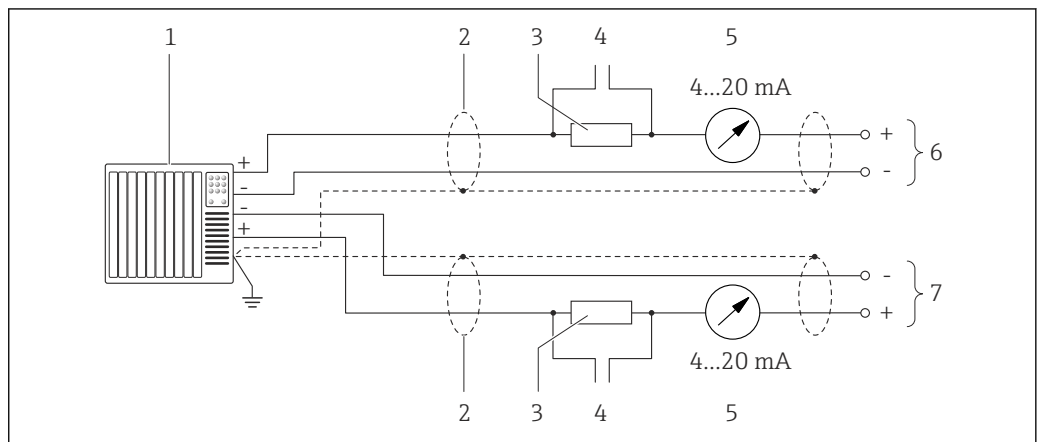

16 Connection example for HART input (master mode) via current output (active)

- 1 Automation system with current input (e.g. PLC).
Prerequisite: automation system with HART version 6, HART commands 113 and 114 can be processed.
- 2 Cable shield, observe cable specifications (→ 32)
- 3 Resistor for HART communication ($\geq 250 \Omega$): observe maximum load (→ 9)
- 4 Connection for HART operating devices (→ 83)
- 5 Analog display unit
- 6 Transmitter
- 7 Sensor for external measured variable

Potential equalization

No special measures for potential equalization are required.



For devices intended for use in hazardous locations, please observe the guidelines in the Ex documentation (XA).

Terminals

Transmitter

Spring terminals for wire cross-sections 0.5 to 2.5 mm^2 (20 to 14 AWG)

Safety Barrier Promass 100

Plug-in screw terminals for wire cross-sections 0.5 to 2.5 mm^2 (20 to 14 AWG)

Cable entries

- Cable gland: M20 × 1.5 with cable ϕ 6 to 12 mm (0.24 to 0.47 in)
- Thread for cable entry:
 - NPT ½"
 - G ½"
 - M20

Cable specification**Permitted temperature range**

- -40 °C (-40 °F) to +80 °C (+176 °F)
- Minimum requirement: cable temperature range \geq ambient temperature +20 K

Power supply cable

Standard installation cable is sufficient.

Signal cable*Current output*

For 4-20 mA HART: Shielded cable recommended. Observe grounding concept of the plant.

Pulse/frequency/switch output

Standard installation cable is sufficient.

PROFIBUS DP

The IEC 61158 standard specifies two types of cable (A and B) for the bus line which can be used for every transmission rate. Cable type A is recommended.

Cable type	A
Characteristic impedance	135 to 165 Ω at a measuring frequency of 3 to 20 MHz
Cable capacitance	<30 pF/m
Wire cross-section	>0.34 mm ² (22 AWG)
Cable type	Twisted pairs
Loop resistance	\leq 110 Ω /km
Signal damping	Max. 9 dB over the entire length of the cable cross-section
Shielding	Copper braided shielding or braided shielding with foil shield. When grounding the cable shield, observe the grounding concept of the plant.

Modbus RS485

The EIA/TIA-485 standard specifies two types of cable (A and B) for the bus line which can be used for every transmission rate. Cable type A is recommended.

Cable type	A
Characteristic impedance	135 to 165 Ω at a measuring frequency of 3 to 20 MHz
Cable capacitance	<30 pF/m
Wire cross-section	>0.34 mm ² (22 AWG)
Cable type	Twisted pairs
Loop resistance	\leq 110 Ω /km
Signal damping	Max. 9 dB over the entire length of the cable cross-section
Shielding	Copper braided shielding or braided shielding with foil shield. When grounding the cable shield, observe the grounding concept of the plant.

EtherNet/IP


The standard ANSI/TIA/EIA-568-B.2 Annex specifies CAT 5 as the minimum category for a cable used for EtherNet/IP. CAT 5e and CAT 6 are recommended.

 For more information on planning and installing EtherNet/IP networks, please refer to the "Media Planning and Installation Manual. EtherNet/IP" of the ODVA Organization.

Connecting cable between Safety Barrier Promass 100 and measuring device

Cable type	Shielded twisted-pair cable with 2x2 wires. When grounding the cable shield, observe the grounding concept of the plant.
Maximum cable resistance	2,5 Ω, one side

► Comply with the maximum cable resistance specifications to ensure the operational reliability of the measuring device.



The maximum cable length for individual wire cross-sections is specified in the table below. Observe the maximum capacitance and inductance per unit length of the cable and connection values for hazardous areas (→  12).

Wire cross-section		Maximum cable length	
[mm ²]	[AWG]	[m]	[ft]
0.5	20	70	230
0.75	18	100	328
1.0	17	100	328
1.5	16	200	656
2.5	14	300	984

Performance characteristics

Reference operating conditions

- Error limits based on ISO 11631
- Water with +15 to +45 °C (+59 to +113 °F) at 2 to 6 bar (29 to 87 psi)
- Specifications as per calibration protocol
- Accuracy based on accredited calibration rigs that are traced to ISO 17025.

 To obtain measured errors, use the *Applicator* sizing tool (→  90)

Maximum measured error

o.r. = of reading; 1 g/cm³ = 1 kg/l; T = medium temperature


Base accuracy

Mass flow and volume flow (liquids)

±0.10 %

Mass flow (gases)

±0.50 % o.r.

 Design fundamentals (→  36)

Density (liquids)

- Reference conditions: ±0.0005 g/cm³
- Standard density calibration: ±0.01 g/cm³
(valid over the entire temperature range and density range)
- Wide-range density specification (order code for "Application package", option EF "Special density and concentration") : ±0.002 g/cm³ (valid range for special density calibration: 0.0 to 2 g/cm³, +5 to +80 °C (+41 to +176 °F))

Temperature

±0.5 °C ± 0.005 · T °C (±0.9 °F ± 0.003 · (T - 32) °F)

Zero point stability

DN		Zero point stability	
[mm]	[in]	[kg/h]	[lb/min]
8	$\frac{3}{8}$	0.20	0.007
15	$\frac{1}{2}$	0.65	0.024
25	1	1.80	0.066
40	$1\frac{1}{2}$	4.50	0.165
50	2	7.0	0.257

Flow values

Flow values as turndown parameter depending on nominal diameter.

SI units


DN	1:1	1:10	1:20	1:50	1:100	1:500
[mm]	[kg/h]	[kg/h]	[kg/h]	[kg/h]	[kg/h]	[kg/h]
8	2 000	200	100	40	20	4
15	6 500	650	325	130	65	13
25	18 000	1 800	900	360	180	36
40	45 000	4 500	2 250	900	450	90
50	70 000	7 000	3 500	1 400	700	140

US units

DN	1:1	1:10	1:20	1:50	1:100	1:500
[inch]	[lb/min]	[lb/min]	[lb/min]	[lb/min]	[lb/min]	[lb/min]
$\frac{3}{8}$	73.50	7.350	3.675	1.470	0.735	0.147
$\frac{1}{2}$	238.9	23.89	11.95	4.778	2.389	0.478
1	661.5	66.15	33.08	13.23	6.615	1.323
$1\frac{1}{2}$	1 654	165.4	82.70	33.08	16.54	3.308
2	2 573	257.3	128.7	51.46	25.73	5.146

Accuracy of outputs

o.r. = of reading; o.f.s. = of full scale value

 The output accuracy must be factored into the measured error if analog outputs are used, but can be ignored for fieldbus outputs (e.g. Modbus RS485, EtherNet/IP).

Current output

Accuracy	Max. ± 0.05 % o.f.s. or ± 5 μ A
-----------------	---

Pulse/frequency output

Accuracy	Max. ± 50 ppm o.r.
-----------------	------------------------

Repeatability

o.r. = of reading; $1 \text{ g/cm}^3 = 1 \text{ kg/l}$; T = medium temperature

Base repeatability

Mass flow and volume flow (liquids)

±0.05 % o.r.

Mass flow (gases)

±0.25 % o.r.



Design fundamentals (→ 36)

Density (liquids)

±0.00025 g/cm³

Temperature

±0.25 °C ± 0.0025 · T °C (±0.45 °F±0.0015 · (T-32) °F)

Response time

The response time depends on the configuration (damping).

Influence of ambient temperature

o.r. = of reading; o.f.s. = of full scale value

Current output

Temperature coefficient	Max. ±50 ppm/°C o.f.s. or ±1 µA/°C
--------------------------------	------------------------------------

Pulse/frequency output

Temperature coefficient	Max. ±50 ppm o.r. /100 °C
--------------------------------	---------------------------

Influence of medium temperature

Mass flow and volume flow


When there is a difference between the temperature for zero point adjustment and the process temperature, the typical measured error of the sensor is ±0.0002 % of the full scale value/°C (±0.0001 % of the full scale value/°F).

Density

When there is a difference between the density calibration temperature and the process temperature, the typical measured error of the sensor is ±0.0001 g/cm³ /°C (±0.00005 g/cm³ /°F). Field density calibration is possible.

Wide-range density specification (special density calibration)

If the process temperature is outside the valid range (→ 33) the measured error is ±0.0001 g/cm³ /°C (±0.00005 g/cm³ /°F)

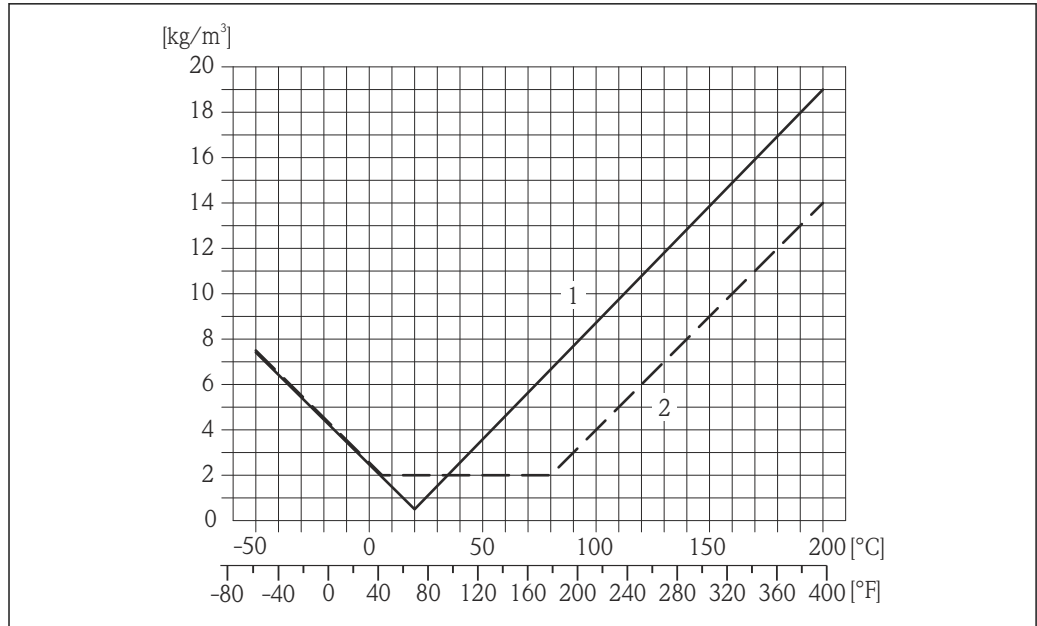

- 1 Field density calibration, for example at +20 °C (+68 °F)
- 2 Special density calibration

Temperature

$$\pm 0.005 \cdot T \text{ } ^\circ\text{C} (\pm 0.005 \cdot (T - 32) \text{ } ^\circ\text{F})$$

Influence of medium pressure

The table below shows the effect on accuracy of mass flow due to a difference between calibration pressure and process pressure.

o.r. = of reading

DN		[% o.r./bar]	[% o.r./psi]
[mm]	[in]		
8	3/8	-0.002	-0.0001
15	1/2	-0.006	-0.0004
25	1	-0.005	-0.0003
40	1 1/2	-0.005	-0.0003
50	2	-0.005	-0.0003

Design fundamentals

o.r. = of reading, o.f.s. = of full scale value

BaseAccu = base accuracy in % o.r., BaseRepeat = base repeatability in % o.r.

MeasValue = measured value; ZeroPoint = zero point stability


Calculation of the maximum measured error as a function of the flow rate

Flow rate	Maximum measured error in % o.r.
$\geq \frac{\text{ZeroPoint}}{\text{BaseAccu}} \cdot 100$ <small>A0021332</small>	$\pm \text{BaseAccu}$ <small>A0021339</small>
$< \frac{\text{ZeroPoint}}{\text{BaseAccu}} \cdot 100$ <small>A0021333</small>	$\pm \frac{\text{ZeroPoint}}{\text{MeasValue}} \cdot 100$ <small>A0021334</small>

Calculation of the maximum repeatability as a function of the flow rate

Flow rate	Maximum repeatability in % o.r.
$\geq \frac{1/2 \cdot \text{ZeroPoint}}{\text{BaseRepeat}} \cdot 100$ <small>A0021335</small>	$\pm \text{BaseRepeat}$ <small>A0021340</small>
$< \frac{1/2 \cdot \text{ZeroPoint}}{\text{BaseRepeat}} \cdot 100$ <small>A0021336</small>	$\pm 1/2 \cdot \frac{\text{ZeroPoint}}{\text{MeasValue}} \cdot 100$ <small>A0021337</small>

Example for max. measured error

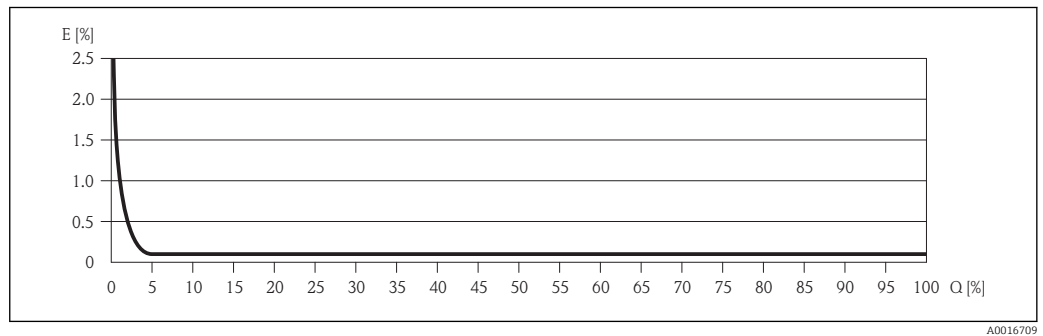

E Error: Maximum measured error as % o.r. (example)
 Q Flow rate as %

Design fundamentals (→ 36)


Installation

No special measures such as supports are necessary. External forces are absorbed by the construction of the device.

Mounting location


To prevent measuring errors arising from accumulation of gas bubbles in the measuring tube, avoid the following mounting locations in the pipe:

- Highest point of a pipeline.
- Directly upstream of a free pipe outlet in a down pipe.

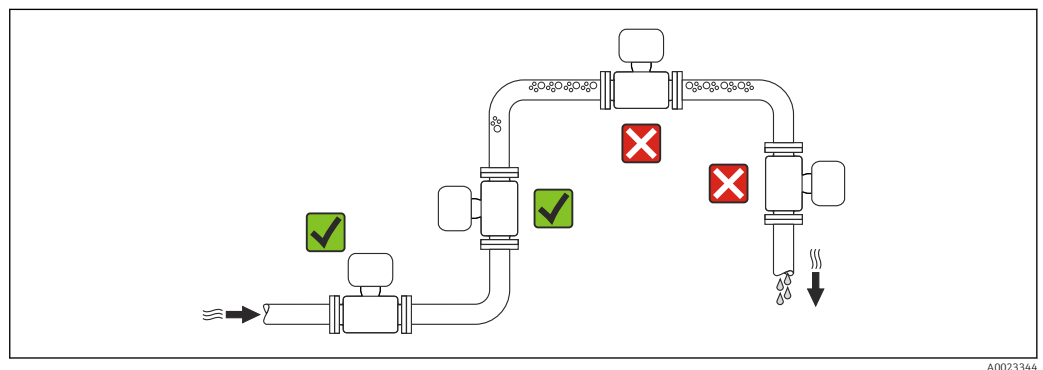

Installation in down pipes

However, the following installation suggestion allows for installation in an open vertical pipeline. Pipe restrictions or the use of an orifice with a smaller cross-section than the nominal diameter prevent the sensor running empty while measurement is in progress.

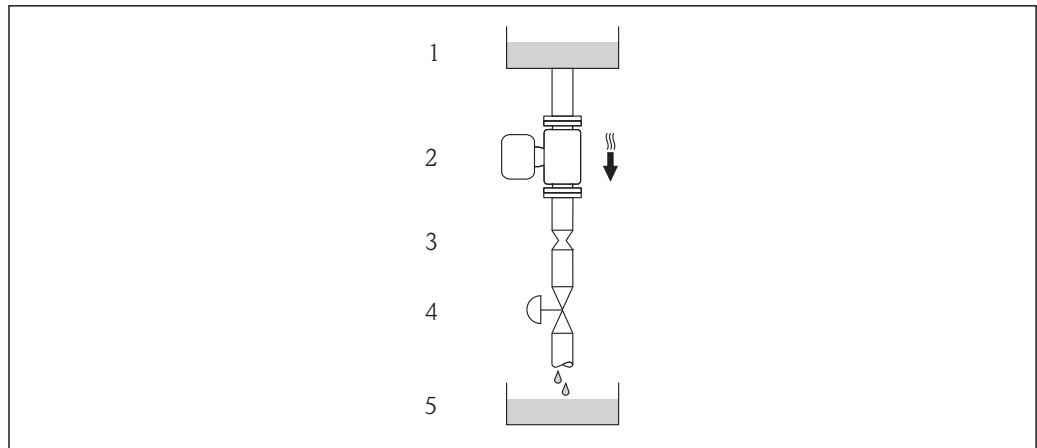

A0015596

17 Installation in a down pipe (e.g. for batching applications)

- 1 Supply tank
- 2 Sensor
- 3 Orifice plate, pipe restriction
- 4 Valve
- 5 Batching tank

DN		Ø orifice plate, pipe restriction	
[mm]	[in]	[mm]	[in]
8	3/8	6	0.24
15	1/2	10	0.40
25	1	14	0.55
40	1 1/2	22	0.87
50	2	28	1.10


Orientation

The direction of the arrow on the sensor nameplate helps you to install the sensor according to the flow direction (direction of medium flow through the piping).

Orientation		Recommendation
A	Vertical orientation	 A0015591
B	Horizontal orientation, transmitter head up	 A0015589
C	Horizontal orientation, transmitter head down	 A0015590
D	Horizontal orientation, transmitter head at side	 A0015592

- 1) Applications with low process temperatures may reduce the ambient temperature. To maintain the minimum ambient temperature for the transmitter, this orientation is recommended.
- 2) Applications with high process temperatures may increase the ambient temperature. To maintain the maximum ambient temperature for the transmitter, this orientation is recommended.

If a sensor is installed horizontally with a curved measuring tube, match the position of the sensor to the fluid properties.

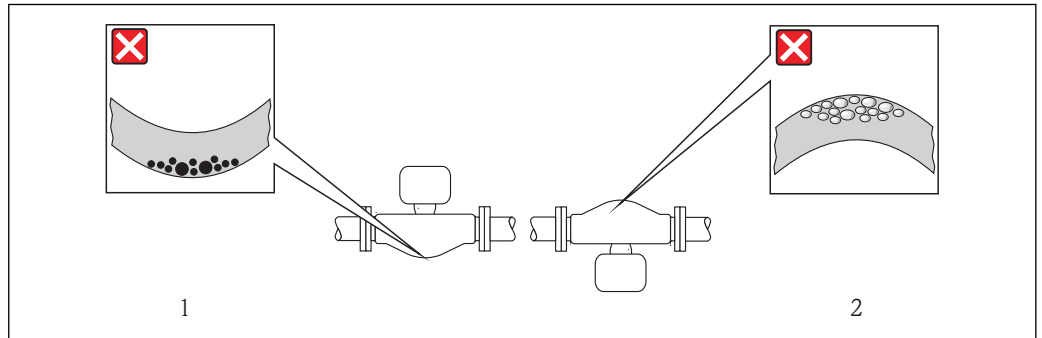

18 Orientation of sensor with curved measuring tube

- 1 Avoid this orientation for fluids with entrained solids: Risk of solids accumulating.
- 2 Avoid this orientation for outgassing fluids: Risk of gas accumulating.

A0014057


Inlet and outlet runs

No special precautions need to be taken for fittings which create turbulence, such as valves, elbows or T-pieces, as long as no cavitation occurs (→ 48).

Special mounting instructions

Guarantees complete drainability

When the sensor is installed in a horizontal line, eccentric clamps can be used to ensure complete drainability. When the system is pitched in a specific direction and at a specific slope, gravity can be used to achieve complete drainability. The sensor must be mounted in the correct position to ensure full drainability in the horizontal position. Markings on the sensor show the correct mounting position to optimize drainability.

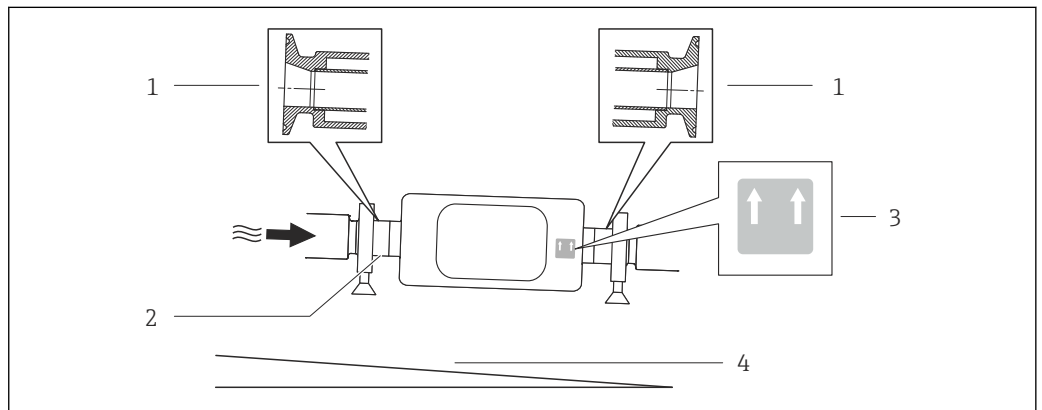


- 1 Eccentric clamp connection
- 2 Line on the underside indicates the lowest point of the eccentric process connection.
- 3 "This side up" label indicates which side is up
- 4 Slope the device in accordance with the hygiene guidelines. Slope: approx. 2 % or 21mm/m (0.24 in/feet)

A0016583

Securing with mounting clamp in the case of hygiene connections

It is not necessary to provide additional support for the sensor for operational performance purposes. If, however, additional support is required for installation purposes, the following dimensions must be observed.

Use mounting clamp with lining between clamp and measuring instrument.

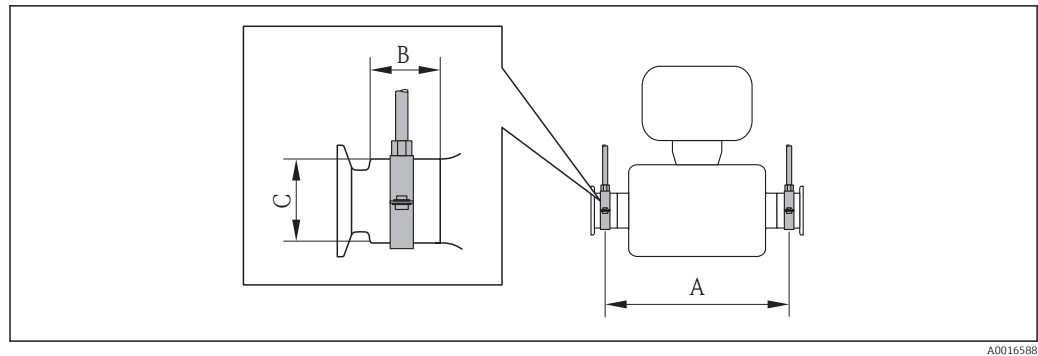

A0016588

SI units

DN [mm]	8	15	25	40	50
A [mm]	298	402	542	750	1019
B [mm]	33	33	33	36.5	44.1
C [mm]	28	28	38	56	75

US units

DN [in]	$\frac{3}{8}$	$\frac{1}{2}$	1	1 $\frac{1}{2}$	2
A [in]	11.73	15.83	21.34	29.53	40.12
B [in]	1.3	1.3	1.3	1.44	1.74
C [in]	1.1	1.1	1.5	2.2	2.95


Zero point adjustment

All measuring devices are calibrated in accordance with state-of-the-art technology. Calibration takes place under reference conditions (\rightarrow 33). Therefore, a zero point adjustment in the field is generally not required.

Experience shows that zero point adjustment is advisable only in special cases:

- To achieve maximum measuring accuracy even with low flow rates
- Under extreme process or operating conditions (e.g. very high process temperatures or very high-viscosity fluids).

**Mounting Safety Barrier
Promass 100**

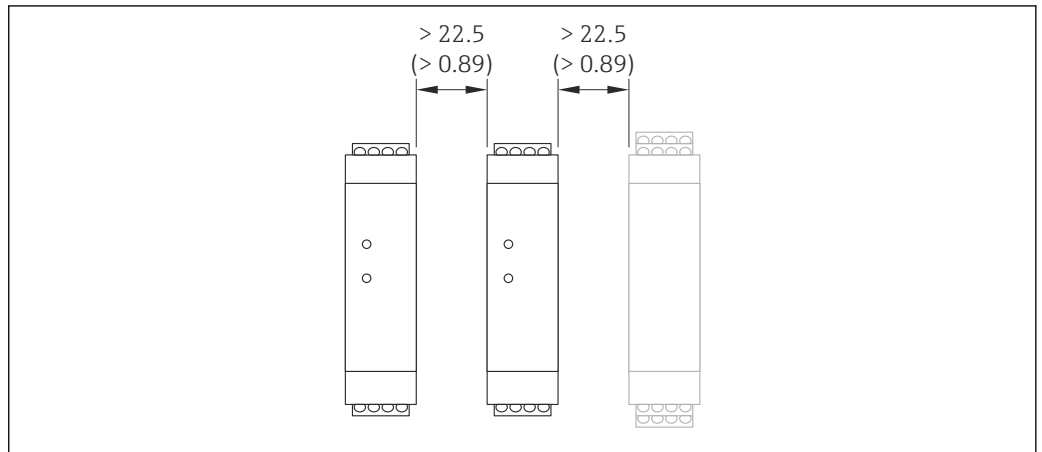

A0016894

19 Minimum distance between additional Safety Barrier Promass 100 or other modules. Engineering unit mm (in)

Environment

Ambient temperature range

Measuring device	Non-Ex	-40 to +60 °C (-40 to +140 °F)
	Ex na, NI version	-40 to +60 °C (-40 to +140 °F)
	Ex ia, IS version	<ul style="list-style-type: none"> ■ -40 to +60 °C (-40 to +140 °F) ■ -50 to +60 °C (-58 to +140 °F) (Order code for "Test, certificate", option JM)
Local display		-20 to +60 °C (-4 to +140 °F) The readability of the display may be impaired at temperatures outside the temperature range.
Safety Barrier Promass 100		-40 to +60 °C (-40 to +140 °F)

- ▶ If operating outdoors:
Avoid direct sunlight, particularly in warm climatic regions.

i Weather protection covers can be ordered from Endress+Hauser: see "Accessories" section

Temperature tables

In the following tables, the following interdependencies between the maximum medium temperature for T1-T6 and the maximum ambient temperature T_a apply when operating the device in hazardous areas.

Ex ia, cCSA_{US} IS

SI units

Order code for "Housing"	T_a [°C]	T6 [85 °C]	T5 [100 °C]	T4 [135 °C]	T3 [200 °C]	T2 [300 °C]	T1 [450 °C]
Option A "Compact coated alu"	35	50	85	120	150 ¹⁾	150 ²⁾	150 ²⁾
Option B "Compact hygienic, stainless"	50	-	85	120	150 ¹⁾	150 ²⁾	150 ²⁾
	60	-	-	120	150 ¹⁾	150 ²⁾	150 ²⁾
Option C "Ultra compact hygienic, stainless, M12 device plug"	35	50	85	120	150 ¹⁾	150 ²⁾	150 ²⁾

Order code for "Housing"	T _a [°C]	T6 [85 °C]	T5 [100 °C]	T4 [135 °C]	T3 [200 °C]	T2 [300 °C]	T1 [450 °C]
	45	–	85	120	150 ¹⁾	150 ²⁾	150 ²⁾
	50	–	–	120	150 ¹⁾	150 ²⁾	150 ²⁾

- 1) The following applies for specified sensors with a maximum medium temperature $T_m = 200\text{ °C}$: $T_m = 170\text{ °C}$
- 2) The following applies for specified sensors with a maximum medium temperature $T_m = 200\text{ °C}$: $T_m = 200\text{ °C}$

US units

Order code for "Housing"	T _a [°F]	T6 [185 °F]	T5 [212 °F]	T4 [275 °F]	T3 [392 °F]	T2 [572 °F]	T1 [842 °F]
Option A "Compact coated alu"	95	122	185	248	302 ¹⁾	302 ²⁾	302 ²⁾
Option B "Compact hygienic, stainless"	122	–	185	248	302 ¹⁾	302 ²⁾	302 ²⁾
	140	–	–	248	302 ¹⁾	302 ²⁾	302 ²⁾
Option C "Ultra compact hygienic, stainless, M12 device plug"	95	122	185	248	302 ¹⁾	302 ²⁾	302 ²⁾
	113	–	185	248	302 ¹⁾	302 ²⁾	302 ²⁾
	122	–	–	248	302 ¹⁾	302 ²⁾	302 ²⁾

- 1) The following applies for specified sensors with a maximum medium temperature $T_m = 392\text{ °F}$: $T_m = 338\text{ °F}$
- 2) The following applies for specified sensors with a maximum medium temperature $T_m = 392\text{ °F}$: $T_m = 392\text{ °F}$

Ex nA, cCSA_{US} NI*SI units*

Order code for "Housing"	T _a [°C]	T6 [85 °C]	T5 [100 °C]	T4 [135 °C]	T3 [200 °C]	T2 [300 °C]	T1 [450 °C]
Option A "Compact coated alu"	35	50	85	120	150 ¹⁾	150 ²⁾	150
Option B "Compact hygienic, stainless"	50	–	85	120	150	150	150
	60	–	–	120	150	150	150
Option C "Ultra compact hygienic, stainless, M12 device plug"	50	–	85	120	150	150	150
	60	–	–	120	150	150	150

- 1) The following applies for specified sensors with a maximum medium temperature $T_m = 200\text{ °C}$: $T_m = 170\text{ °C}$
- 2) The following applies for specified sensors with a maximum medium temperature $T_m = 200\text{ °C}$: $T_m = 200\text{ °C}$

US units

Order code for "Housing"	T _a [°F]	T6 [185 °F]	T5 [212 °F]	T4 [275 °F]	T3 [392 °F]	T2 [572 °F]	T1 [842 °F]
Option A "Compact coated alu"	95	122	185	248	302 ¹⁾	302 ²⁾	302
Option B "Compact hygienic, stainless"	122	–	185	248	302	302	302
	140	–	–	248	302	302	302
Option C "Ultra compact hygienic, stainless, M12 device plug"	122	–	185	248	302	302	302
	140	–	–	248	302	302	302

- 1) The following applies for specified sensors with a maximum medium temperature $T_m = 392\text{ °F}$: $T_m = 338\text{ °F}$
- 2) The following applies for specified sensors with a maximum medium temperature $T_m = 392\text{ °F}$: $T_m = 392\text{ °F}$

Explosion hazards arising from gas and dust

Determining the temperature class and surface temperature with the temperature table

- In the case of gas: Determine the temperature class as a function of the ambient temperature T_a and the medium temperature T_m .
- In the case of dust: Determine the maximum surface temperature as a function of the maximum ambient temperature T_a and the maximum medium temperature T_m .

Example

- Measured maximum ambient temperature: $T_{ma} = 47\text{ °C}$
- Measured maximum medium temperature: $T_{mm} = 108\text{ °C}$

	Ta [°C]	T6 [85°C]	T5 [100°C]	T4 [135°C]	T3 [200°C]	T2 [300°C]	T1 [450°C]
	35	50	85	120	140	140	140
	50	-	85	120	140	140	140
	60	-	-	120	140	140	140
	35	50	85	120	140	140	140
	45	-	85	120	140	140	140
	50	-	-	120	140	140	140

20 Procedure for determining the maximum surface temperature

1. Select device.
2. In the column for the maximum ambient temperature T_a select the temperature that is immediately greater than or equal to the measured maximum ambient temperature T_{ma} that is present.
 - ↳ $T_a = 50\text{ °C}$.
 - The row showing the maximum medium temperature is determined.
3. Select the maximum medium temperature T_m of this row, which is larger or equal to the measured maximum medium temperature T_{mm} .
 - ↳ The column with the temperature class for gas is determined: $108\text{ °C} \leq 120\text{ °C} \rightarrow T4$.
4. The maximum temperature of the temperature class determined corresponds to the maximum surface temperature for dust: $T4 = 135\text{ °C}$

Storage temperature	-50 to +60 °C (-58 to +140 °F) (Order code for "Test, certificate", option JM)
Climate class	DIN EN 60068-2-38 (test Z/AD)
Degree of protection	<p>Transmitter and sensor</p> <ul style="list-style-type: none"> ■ As standard: IP66/67, type 4X enclosure ■ With the order code for "Sensor options", option CM: IP69K can also be ordered ■ When housing is open: IP20, type 1 enclosure ■ Display module: IP20, type 1 enclosure <p>Safety Barrier Promass 100 IP20</p>
Shock resistance	As per IEC/EN 60068-2-31
Vibration resistance	Acceleration up to 1 g, 10 to 150 Hz, based on IEC/EN 60068-2-6
Interior cleaning	<ul style="list-style-type: none"> ■ Sterilization in place (SIP) ■ Cleaning in place (CIP) ■ Cleaning with pigs

Electromagnetic compatibility (EMC)

- As per IEC/EN 61326 and NAMUR Recommendation 21 (NE 21)
- Complies with emission limits for industry as per EN 55011 (Class A)
- Device version with PROFIBUS DP: Complies with emission limits for industry as per EN 50170 Volume 2, IEC 61784



The following applies for PROFIBUS DP: If baud rates > 1.5 Mbaud an EMC cable entry must be used and the cable shield must continue as far as the terminal wherever possible.



For details refer to the Declaration of Conformity.

Process**Medium temperature range****Sensor**

- -50 to +150 °C (-58 to +302 °F)
- -50 to +200 °C (-58 to +392 °F) with extended temperature (order code for "Measuring tube mat.", option TD, TG)

Seals


No internal seals

Density

0 to 5 000 kg/m³ (0 to 312 lb/cf)

Pressure-temperature ratings

The following pressure-temperature ratings refer to the entire device and not just the process connection.

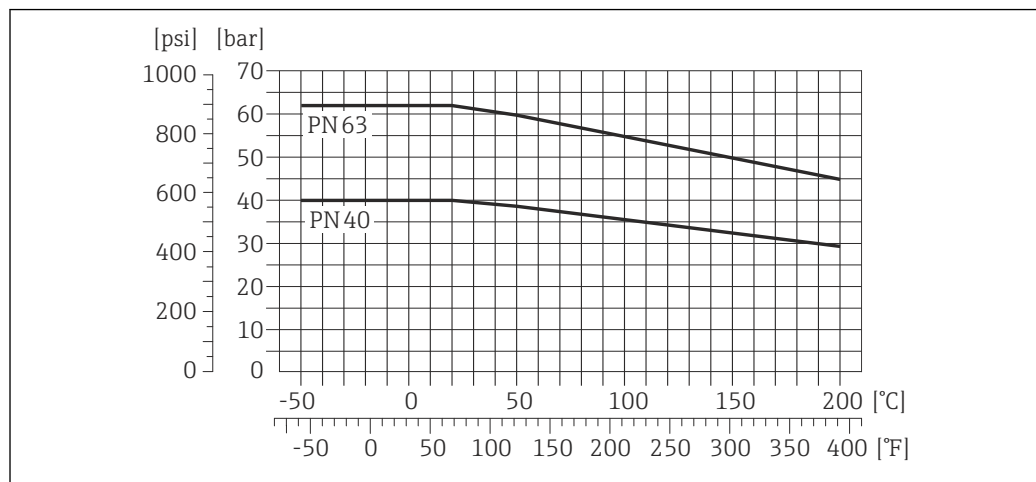
Flange connection according to EN 1092-1 (DIN 2501)

A0020994-EN

21 With flange material 1.4404 (316/316L).

The material load curves for the temperature range +150 to +200 °C (+302 to +392 °F) apply only to the order code for "Measuring tube material", option TD, TG

Flange connection according to ASME B16.5

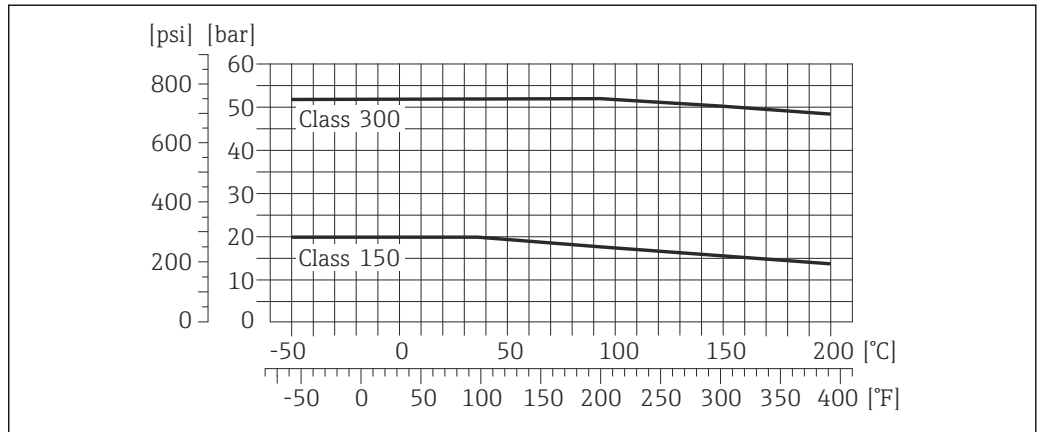


A0020997-EN

22 With flange material 1.4404 (316/316L)

The material load curves for the temperature range +150 to +200 °C (+302 to +392 °F) apply only to the order code for "Measuring tube material", option TD, TG

Flange connection according to JIS B2220

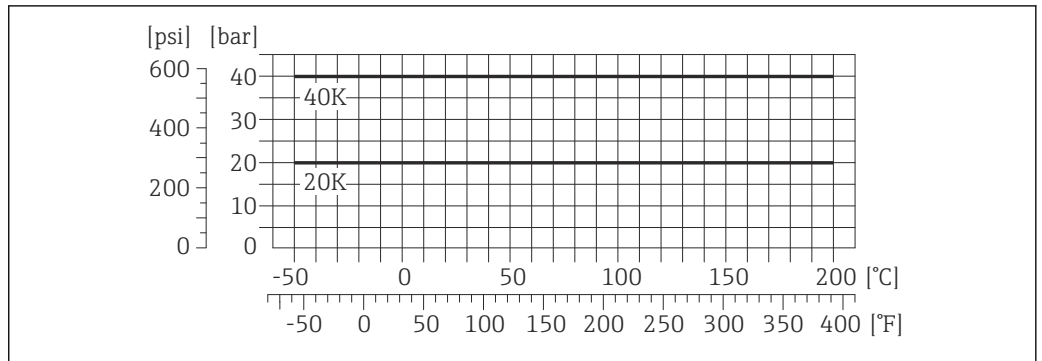


A0020998-EN

23 With flange material 1.4404 (316/316L)

The material load curves for the temperature range +150 to +200 °C (+302 to +392 °F) apply only to the order code for "Measuring tube material", option TD, TG

Process connection to DIN 11851

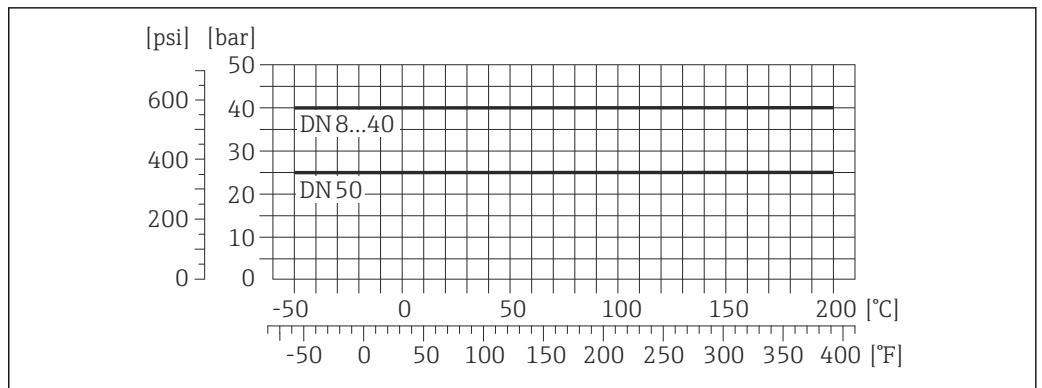


A0020995-EN

24 With connection material 1.4404 (316L)

The material load curves for the temperature range +150 to +200 °C (+302 to +392 °F) apply only to the order code for "Measuring tube material", option TD, TG

DIN 11851 allows for applications up to +140 °C (+284 °F) if suitable sealing materials are used. Please take this into account when selecting seals and counterparts, as these components can limit the pressure and temperature range.


Aseptic threaded hygienic connection according to DIN 11864-1 Form A; aseptic clamp connection according to DIN 11864-3 Form A

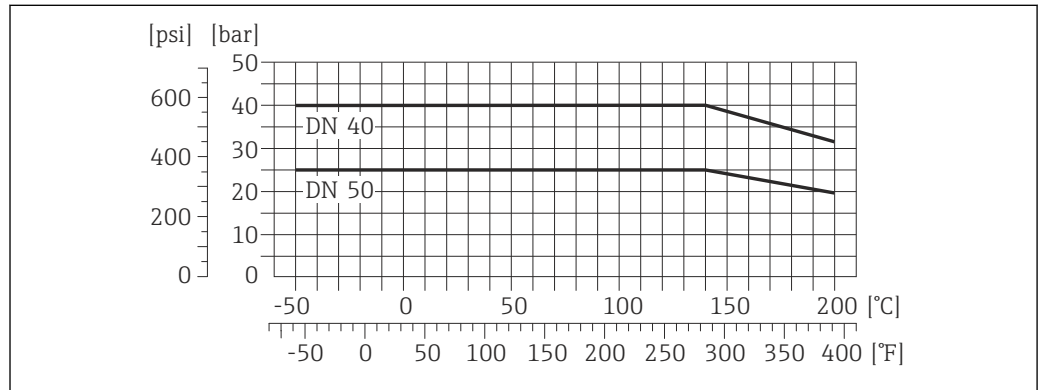

25 With connection material 1.4404 (316L)

The material load curves for the temperature range +150 to +200 °C (+302 to +392 °F) apply only to the order code for "Measuring tube material", option TD, TG


Aseptic flange connection according to DIN 11864-2 Form A

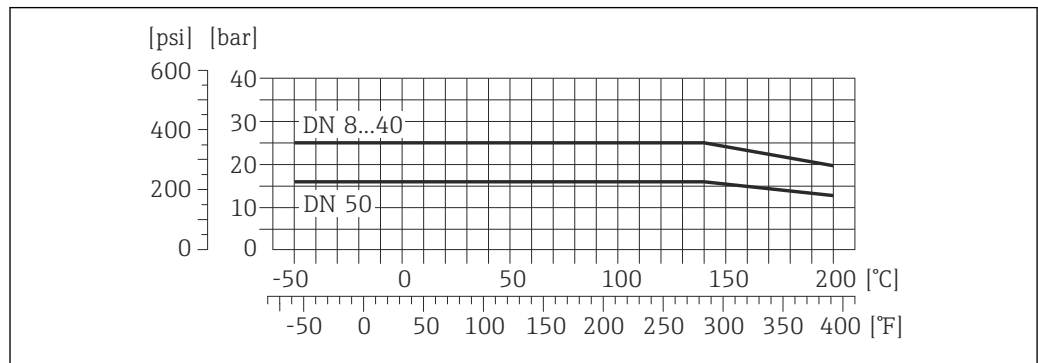

26 With flange material 1.4404 (316L)

The material load curves for the temperature range +150 to +200 °C (+302 to +392 °F) apply only to the order code for "Measuring tube material", option TD, TG

Process connection to SMS 1145

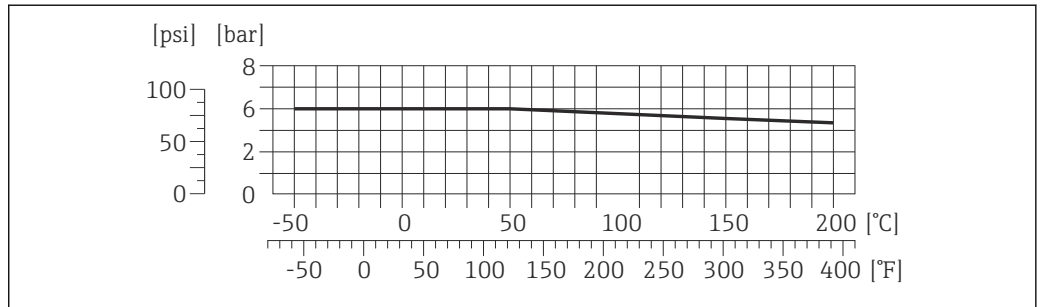


A0020985-EN

27 With connection material 1.4404 (316L)

The material load curves for the temperature range +150 to +200 °C (+302 to +392 °F) apply only to the order code for "Measuring tube material", option TD, TG

SMS 1145 allows for applications up to 6 bar (87 psi) if suitable sealing materials are used. Please take this into account when selecting seals and counterparts, as these components can limit the pressure and temperature range.

Threaded hygienic connection to ISO 2853

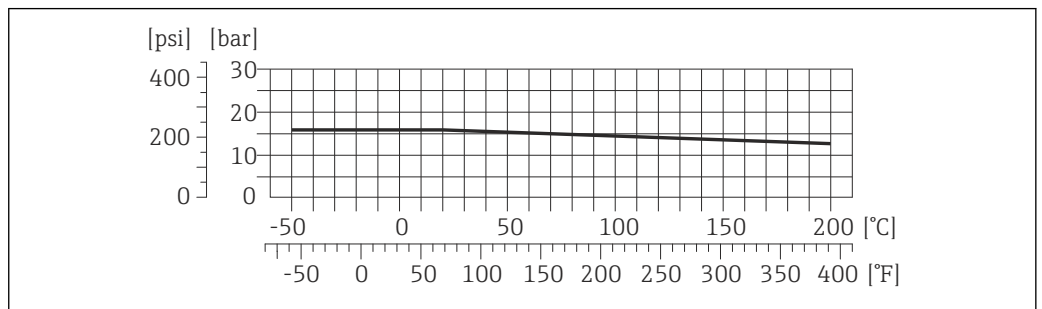


A0020964-EN

28 With connection material 1.4404 (316L)

The material load curves for the temperature range +150 to +200 °C (+302 to +392 °F) apply only to the order code for "Measuring tube material", option TD, TG

Neumo BioConnect (flange with groove), BBS (flange with groove), BBS (aseptic threaded hygienic connection)

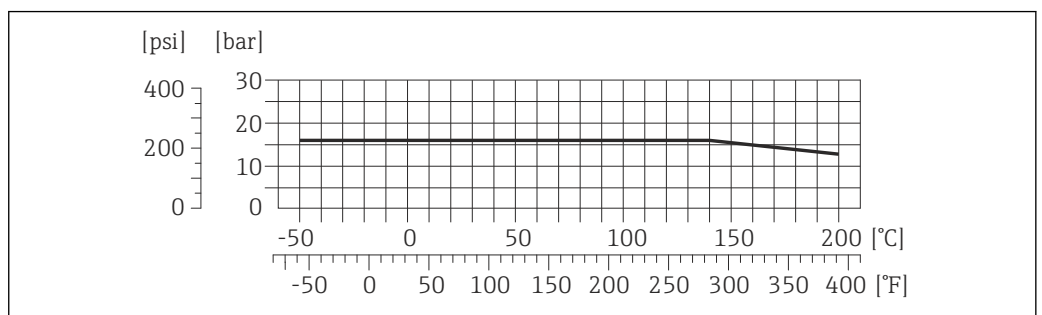

A0021004-EN

29 With flange material 1.4404 (316L)

The material load curves for the temperature range +150 to +200 °C (+302 to +392 °F) apply only to the order code for "Measuring tube material", option TD, TG

Tri-Clamp, ISO 2852 (clamp), DIN 32676 (clamp connection), BBS (clamp with notch), Neumo BioConnect (clamp with notch)

The clamp connections are suitable up to a maximum pressure of 16 bar (232 psi). Please observe the operating limits of the clamp and seal used as they could be under 16 bar (232 psi). The clamp and seal are not included in the scope of supply.

Secondary containment pressure rating

The sensor housing is filled with dry nitrogen and protects the electronics and mechanics inside.


DN		Secondary containment nominal pressure (designed with a safety factor ≥ 4)		Secondary containment burst pressure	
[mm]	[in]	[bar]	[psi]	[bar]	[psi]
8	$\frac{3}{8}$	25	362	190	2755
15	$\frac{1}{2}$	25	362	175	2535
25	1	25	362	165	2930
40	$1\frac{1}{2}$	16	232	64	925
50	2	10	145	54	780

i If there is a risk of measuring tube failure due to process characteristics, e.g. with corrosive process fluids, we recommend the use of sensors whose secondary containment is equipped with special pressure monitoring connections (order code for "Purge connection", option CH).

With the help of these connections, the fluid collected in the secondary containment can be bled off in the event of tube failure. This is especially important in high-pressure gas applications. These connections can also be used for gas purging (gas detection).


Do not open the purge connections unless the containment can be filled immediately with a dry, inert gas. Use only low gauge pressure to purge. Maximum pressure: 5 bar (72.5 psi).


If a device fitted with purge connections is connected to the purge system, the maximum nominal pressure is determined by the purge system itself or by the device, depending on which component has the lower nominal pressure.

Dimensions: (→  79)


Flow limit

Select the nominal diameter by optimizing between the required flow range and permissible pressure loss.

i For an overview of the measuring range full scale values, see the "Measuring range" section (→  8)

- The minimum recommended full scale value is approx. 1/20 of the maximum full scale value
- In most applications, 20 to 50 % of the maximum full scale value can be considered ideal
- Select a lower full scale value for abrasive substances (such as liquids with entrained solids): flow velocity < 1 m/s (< 3 ft/s).
- For gas measurement the following rules apply:
 - The flow velocity in the measuring tubes should not exceed half the sonic velocity (0.5 Mach).
 - The maximum mass flow depends on the density of the gas: formula (→  8)

Pressure loss


i To calculate the pressure loss, use the *Applicator* sizing tool (→  90)

System pressure

It is important that cavitation does not occur, or that gases entrained in the liquids do not outgas. This is prevented by means of a sufficiently high system pressure.

For this reason, the following mounting locations are recommended:


- At the lowest point in a vertical pipe
- Downstream from pumps (no danger of vacuum)

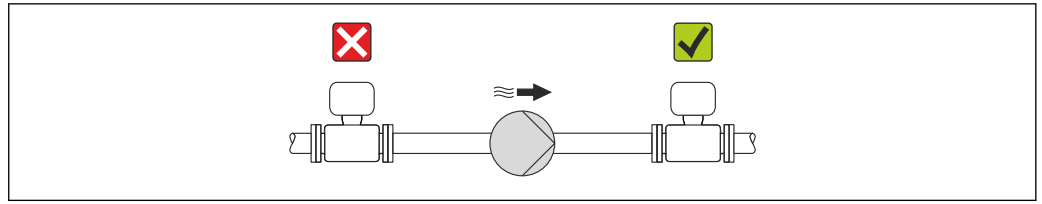

A0015594

Thermal insulation

In the case of some fluids, it is important that the heat radiated from the sensor to the transmitter is kept to a minimum. A wide range of materials can be used for the required insulation.

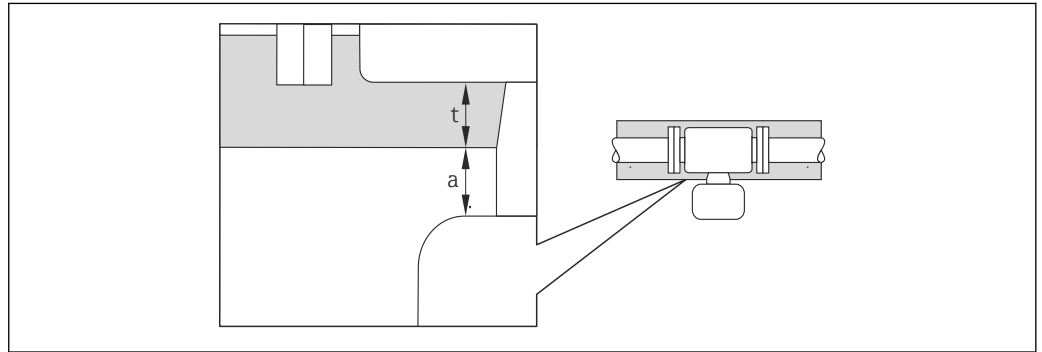


A0019919

- a* Minimum distance to insulation
- t* Insulation thickness

The minimum distance between the transmitter housing and the insulation is 10 mm (0.39 in) so that the transmitter head remains completely exposed.

Insulation thickness for version without neck extension

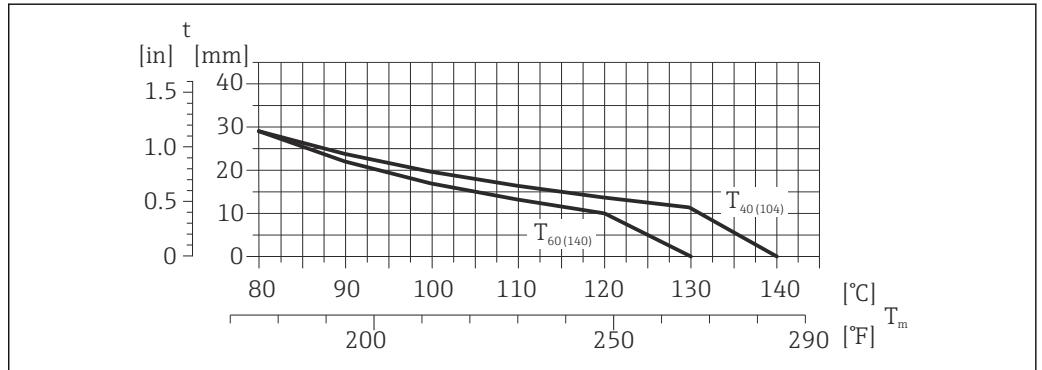


A0023173

30 Recommended insulation thicknesses depending on the medium and ambient temperature

- t* Insulation thickness
- T_m Medium temperature
- $T_{40(104)}$ Insulation thickness with ambient temperature of $T_a = 40^\circ\text{C}$ (104 °F)
- $T_{60(140)}$ Insulation thickness with ambient temperature of $T_a = 60^\circ\text{C}$ (140 °F)

Insulation thickness for version with extension neck

Extension neck for the extended temperature range, order code for "Sensor option", option CG or order code for "Measuring tube material", option TD, TG:

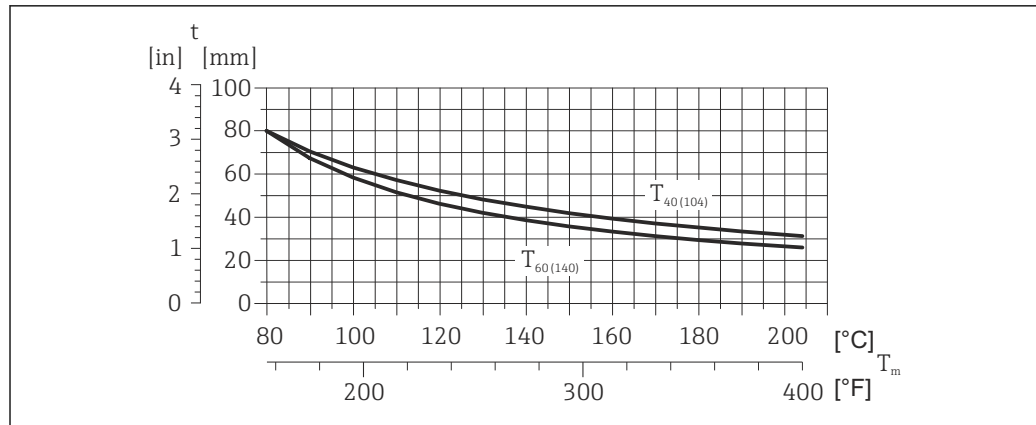

A0023177

31 Recommended insulation thicknesses depending on the medium and ambient temperature

t	Insulation thickness
T_m	Medium temperature
$T_{40(104)}$	Insulation thickness with ambient temperature of $T_a = 40\text{ °C}$ (104 °F)
$T_{60(140)}$	Insulation thickness with ambient temperature of $T_a = 60\text{ °C}$ (140 °F)

NOTICE

The insulation can also be thicker than the recommended insulation thickness.

Prerequisite:


- ▶ The temperature at the lower end of the transmitter housing does not exceed 80 °C (176 °F)
- ▶ Ensure that convection takes place on a sufficiently large scale at the transmitter neck.
- ▶ Ensure that a sufficiently large area of the housing support remains exposed. The uncovered part serves as a radiator and protects the electronics from overheating and excessive cooling.

Heating

Some fluids require suitable measures to avoid loss of heat at the sensor.

Heating options

- Electrical heating, e.g. with electric band heaters
- Via pipes carrying hot water or steam
- Via heating jackets

 Heating jackets for the sensor can be ordered as accessories from Endress+Hauser (→  89).

Vibrations


The high oscillation frequency of the measuring tubes ensures that the correct operation of the measuring system is not influenced by plant vibrations.

Mechanical construction

Design, dimensions

Compact version

Order code for "Housing", option A "Compact coated alu"

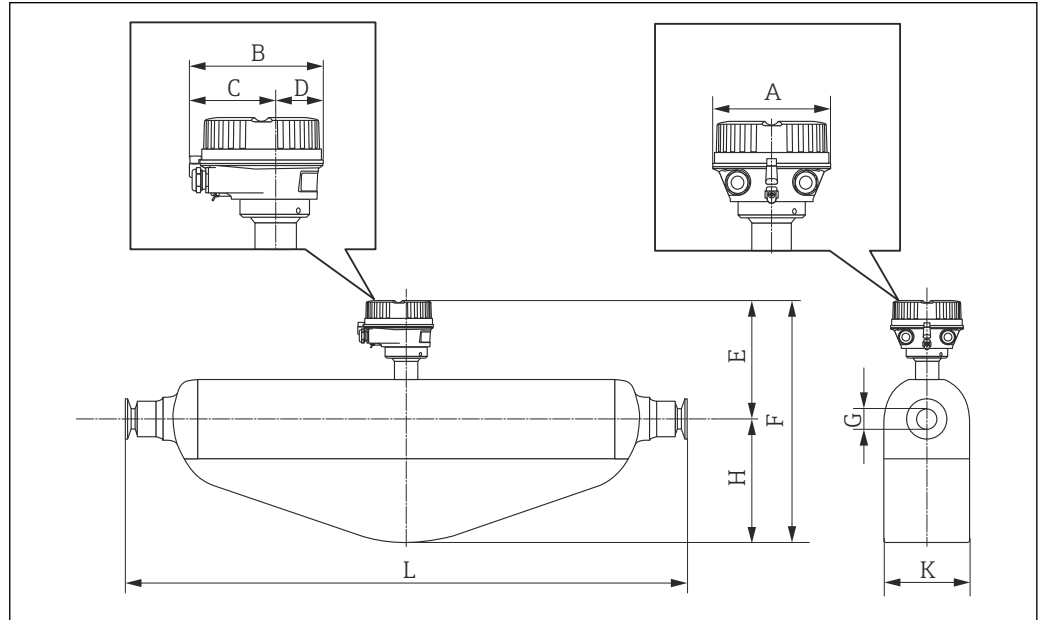

A0016525

Dimensions SI units

DN [mm]	A [mm]	B [mm]	C [mm]	D [mm]	E ^{1) 2)} [mm]	F ^{1) 2)} [mm]	G [mm]	H [mm]	K [mm]	L [mm]
8	136	147.5	93.5	54	191	299	8.31	108	92	³⁾
15	136	147.5	93.5	54	191	299	12.00	108	92	³⁾
25	136	147.5	93.5	54	191	312	17.60	121	92	³⁾
40	136	147.5	93.5	54	215	388	26.00	173	132	³⁾
50	136	147.5	93.5	54	226	467	40.50	241	167	³⁾


- 1) If using an extension neck for the extended temperature range, order code for "Sensor option", option CG and order code for "Measuring tube material", option TD, TG: values +70 mm
- 2) If using a display, order code for "Display; Operation", option B: values +28 mm
- 3) dependent on respective process connection

Dimensions US units

DN [in]	A [in]	B [in]	C [in]	D [in]	E ²⁾ [in]	F ^{1) 2)} [in]	G [in]	H [in]	K [in]	L [in]
³ / ₈	5.35	5.81	3.68	2.13	7.52	11.77	0.33	4.25	3.62	³⁾
¹ / ₂	5.35	5.81	3.68	2.13	7.52	11.77	0.47	4.25	3.62	³⁾
1	5.35	5.81	3.68	2.13	7.52	12.28	0.69	4.76	3.62	³⁾
1½	5.35	5.81	3.68	2.13	8.46	15.28	1.02	6.81	5.2	³⁾
2	5.35	5.81	3.68	2.13	8.9	18.39	1.59	9.49	6.57	³⁾

- 1) If using an extension neck for the extended temperature range, order code for "Sensor option", option CG and order code for "Measuring tube material", option TD, TG: values +2.76 in
- 2) If using a display, order code for "Display; Operation", option B: values +1.1 in
- 3) dependent on respective process connection

Order code for "Housing", option B "Compact hygienic, stainless"

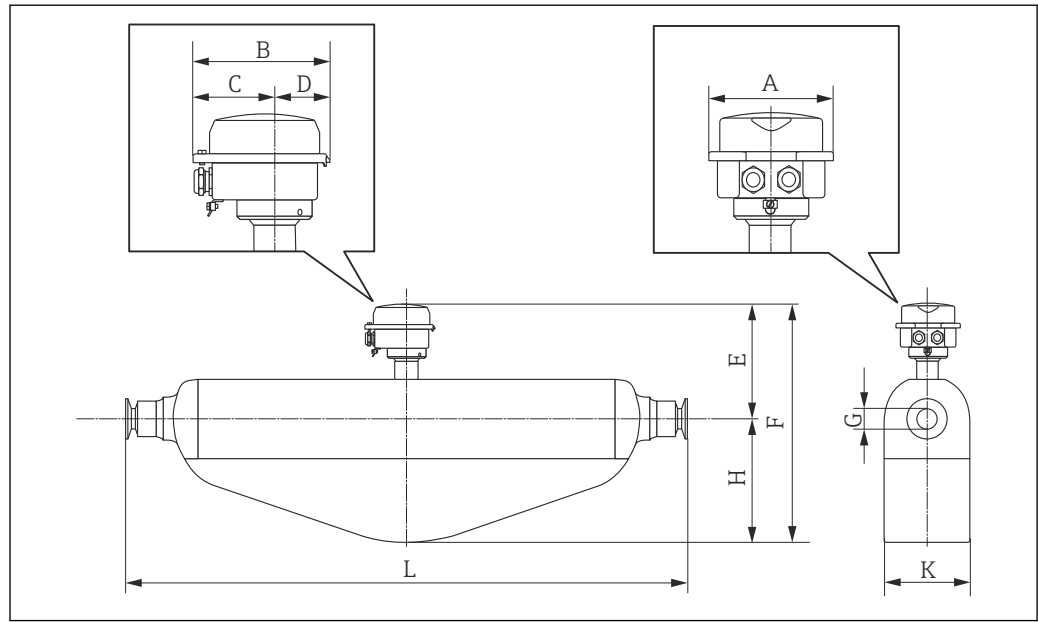

Dimensions SI units

DN [mm]	A [mm]	B [mm]	C [mm]	D [mm]	E ²⁾ [mm]	F ^{1) 2)} [mm]	G [mm]	H [mm]	K [mm]	L [mm]
8	133.5	136.8	78	58.8	186	294	8.31	108	92	³⁾
15	133.5	136.8	78	58.8	186	294	12.00	108	92	³⁾
25	133.5	136.8	78	58.8	186	307	17.60	121	92	³⁾
40	133.5	136.8	78	58.8	210	383	26.00	173	132	³⁾
50	133.5	136.8	78	58.8	221	462	40.50	241	167	³⁾


- 1) If using an extension neck for the extended temperature range, order code for "Sensor option", option CG and order code for "Measuring tube material", option TD, TG: values +70 mm
- 2) If using a display, order code for "Display; Operation", option B: values +14 mm
- 3) dependent on respective process connection

Dimensions US units

DN [in]	A [in]	B [in]	C [in]	D [in]	E ²⁾ [in]	F ^{1) 2)} [in]	G [in]	H [in]	K [in]	L [in]
³ / ₈	5.26	5.39	3.07	2.31	7.32	11.57	0.33	4.25	3.62	³⁾
¹ / ₂	5.26	5.39	3.07	2.31	7.32	11.57	0.47	4.25	3.62	³⁾
1	5.26	5.39	3.07	2.31	7.32	12.09	0.69	4.76	3.62	³⁾
1 ¹ / ₂	5.26	5.39	3.07	2.31	8.27	15.08	1.02	6.81	5.2	³⁾
2	5.26	5.39	3.07	2.31	8.7	18.19	1.59	9.49	6.57	³⁾

- 1) If using an extension neck for the extended temperature range, order code for "Sensor option", option CG and order code for "Measuring tube material", option TD, TG: values +2.76 in
- 2) If using a display, order code for "Display; Operation", option B: values +0.55 in
- 3) dependent on respective process connection

Order code for "Housing", option C "Ultra compact hygienic, stainless, M12 device plug"

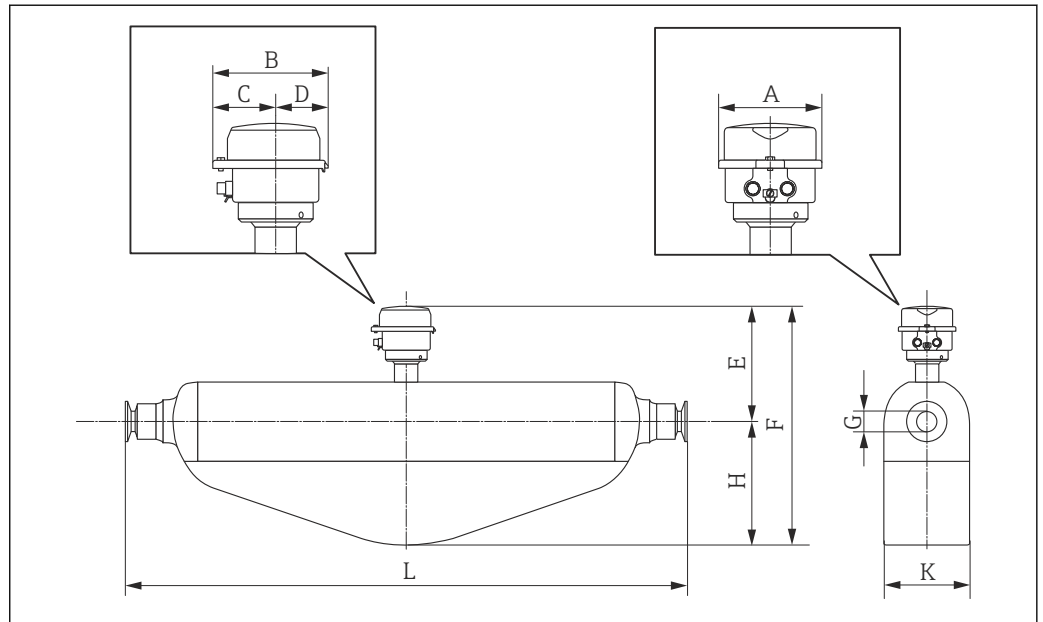

A0016527

Dimensions SI units

DN [mm]	A [mm]	B [mm]	C [mm]	D [mm]	E ²⁾ [mm]	F ^{1) 2)} [mm]	G [mm]	H [mm]	K [mm]	L [mm]
8	111.4	123.6	67.7	55.9	186	294	8.31	108	92	³⁾
15	111.4	123.6	67.7	55.9	186	294	12.00	108	92	³⁾
25	111.4	123.6	67.7	55.9	186	307	17.60	121	92	³⁾
40	111.4	123.6	67.7	55.9	210	383	26.00	173	132	³⁾
50	111.4	123.6	67.7	55.9	221	462	40.50	241	167	³⁾

- 1) If using an extension neck for the extended temperature range, order code for "Sensor option", option CG and order code for "Measuring tube material", option TD, TG: values +70 mm
- 2) If using a display, order code for "Display; Operation", option B: values +14 mm
- 3) dependent on respective process connection


Dimensions US units

DN [in]	A [in]	B [in]	C [in]	D [in]	E ²⁾ [in]	F ^{1) 2)} [in]	G [in]	H [in]	K [in]	L [in]
³ / ₈	4.39	4.87	2.67	2.2	7.32	11.57	0.33	4.25	3.62	³⁾
¹ / ₂	4.39	4.87	2.67	2.2	7.32	11.57	0.47	4.25	3.62	³⁾
1	4.39	4.87	2.67	2.2	7.32	12.09	0.69	4.76	3.62	³⁾
1½	4.39	4.87	2.67	2.2	8.27	15.08	1.02	6.81	5.2	³⁾
2	4.39	4.87	2.67	2.2	8.7	18.19	1.59	9.49	6.57	³⁾

- 1) If using an extension neck for the extended temperature range, order code for "Sensor option", option CG and order code for "Measuring tube material", option TD, TG: values +2.76 in
- 2) If using a display, order code for "Display; Operation", option B: values +0.55 in
- 3) dependent on respective process connection

Process connections in SI units

Flange connections EN (DIN)

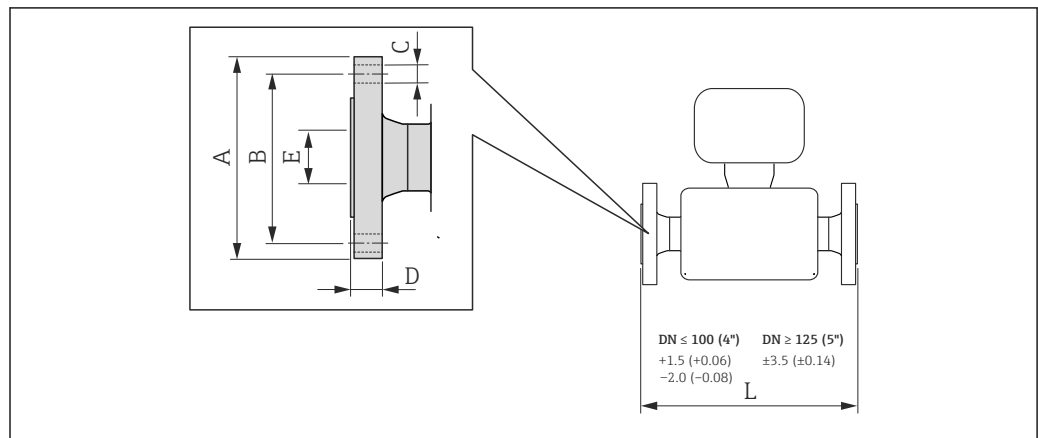

A0015621


32 Engineering unit mm (in)

Flange according to EN 1092-1 (DIN 2501), PN 40: 1.4404 (316/316L) (order code for "Process connection", option D2W)						
Surface roughness (flange): EN 1092-1 Form B1 (DIN 2526 Form C), Ra 3.2 to 12.5 µm						
DN [mm]	A [mm]	B [mm]	C [mm]	D [mm]	E [mm]	L [mm]
8 ¹⁾	95	65	4 × Ø14	17.0	17.3	336
15	95	65	4 × Ø14	20.0	17.3	440
25	115	85	4 × Ø14	19.0	28.5	580
40	150	110	4 × Ø18	21.0	43.1	794
50	165	125	4 × Ø18	25.0	54.5	1071

1) DN 8 with DN 15 flanges as standard

Flange according to EN 1092-1 (DIN 2501), PN 63: 1.4404 (316/316L) (order code for "Process connection", option D3W)						
Surface roughness (flange): EN 1092-1 Form B2 (DIN 2526 Form E), Ra 0.8 to 3.2 µm						
DN [mm]	A [mm]	B [mm]	C [mm]	D [mm]	E [mm]	L [mm]
50	180	135	4 × Ø22	26	54.5	1083

Flange connections ASME B16.5

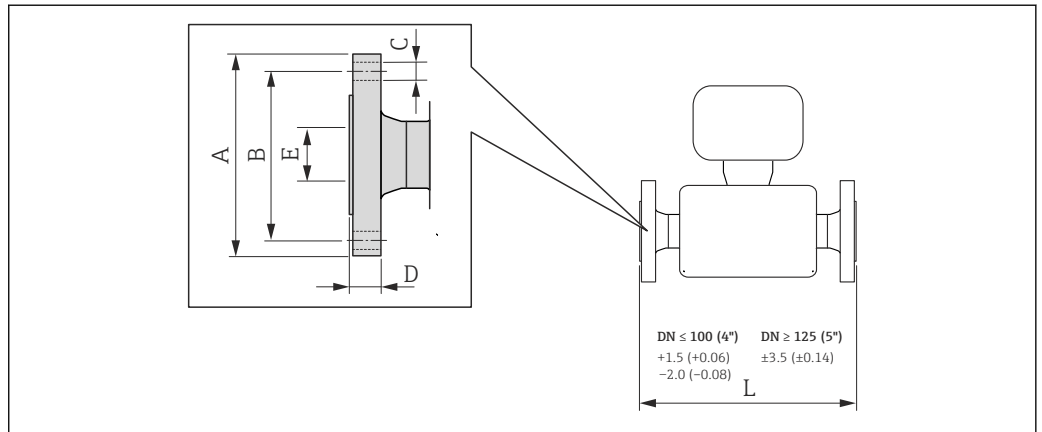

A0015621

33 Engineering unit mm (in)

Flange according to ASME B16.5, Cl 150: 1.4404 (316/316L) (order code for "Process connection", option AAW)
 Surface roughness (flange): Ra 3.2 to 6.3 μm

DN [mm]	A [mm]	B [mm]	C [mm]	D [mm]	E [mm]	L [mm]
8 ¹⁾	90	60.3	4 × Ø15.7	11.2	15.7	336
15	90	60.3	4 × Ø15.7	11.2	15.7	440
25	110	79.4	4 × Ø15.7	14.2	26.7	580
40	125	98.4	4 × Ø15.7	17.5	40.9	794
50	150	120.7	8 × Ø19.1	19.1	52.6	1071


1) DN 8 with DN 15 flanges as standard

Flange according to ASME B16.5, Cl 300: 1.4404 (316/316L) (order code for "Process connection", option ABW)
 Surface roughness (flange): Ra 3.2 to 6.3 μm

DN [mm]	A [mm]	B [mm]	C [mm]	D [mm]	E [mm]	L [mm]
8 ¹⁾	95	66.7	4 × Ø15.7	16.6	15.7	336
15	95	66.7	4 × Ø15.7	16.6	15.7	440
25	125	88.9	4 × Ø19.1	18.1	26.7	580
40	155	114.3	4 × Ø22.3	24.6	40.9	794
50	165	127.0	8 × Ø19.1	27.6	52.6	1071

1) DN 8 with DN 15 flanges as standard

Flange connections JIS

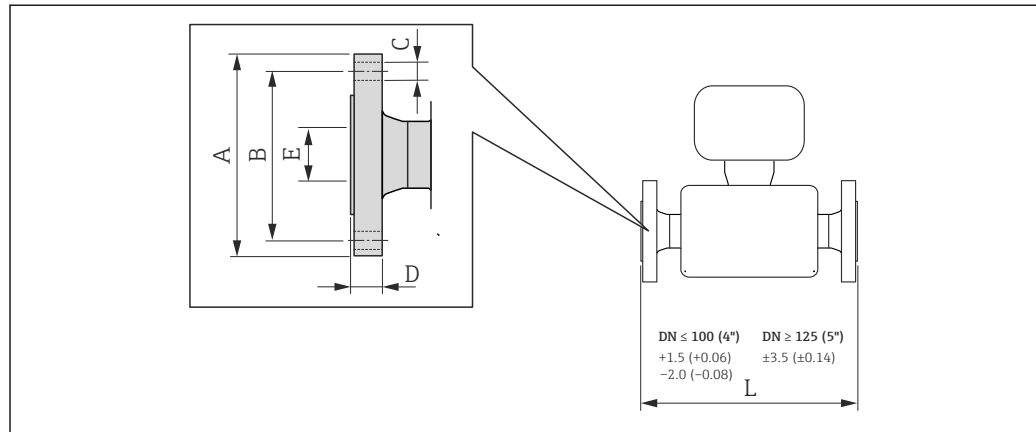

A0015621

34 Engineering unit mm (in)

Flange JIS B2220, 20K: 1.4404 (316/316L) (order code for "Process connection", option NEW)						
Surface roughness (flange): Ra 3.2 to 6.3 µm						
DN [mm]	A [mm]	B [mm]	C [mm]	D [mm]	E [mm]	L [mm]
8 ¹⁾	95	70	4 × Ø15	16.0	15	336
15	95	70	4 × Ø15	16.0	15	440
25	125	90	4 × Ø19	17.5	25	580
40	140	105	4 × Ø19	20.0	40	794
50	155	120	8 × Ø19	27.5	50	1071


1) DN 8 with DN 15 flanges as standard

Flange JIS B2220, 40K: 1.4404 (316/316L) (order code for "Process connection", option NFW)						
Surface roughness (flange): Ra 3.2 to 6.3 µm						
DN [mm]	A [mm]	B [mm]	C [mm]	D [mm]	E [mm]	L [mm]
8 ¹⁾	115	80	4 × Ø19	21.0	15	336
15	115	80	4 × Ø19	21.0	15	440
25	130	95	4 × Ø19	22.0	25	589
40	160	120	4 × Ø23	26.0	38	804
50	165	130	8 × Ø19	26.0	50	1071

1) DN 8 with DN 15 flanges as standard

Tri-Clamp (BS4825, ASME BPE, DIN 11866 line C)

All Tri-Clamp connections (BS4825-3, DN 1", 1½", 2") in accordance with hygienic clamp dimensions as per ASME BPE.

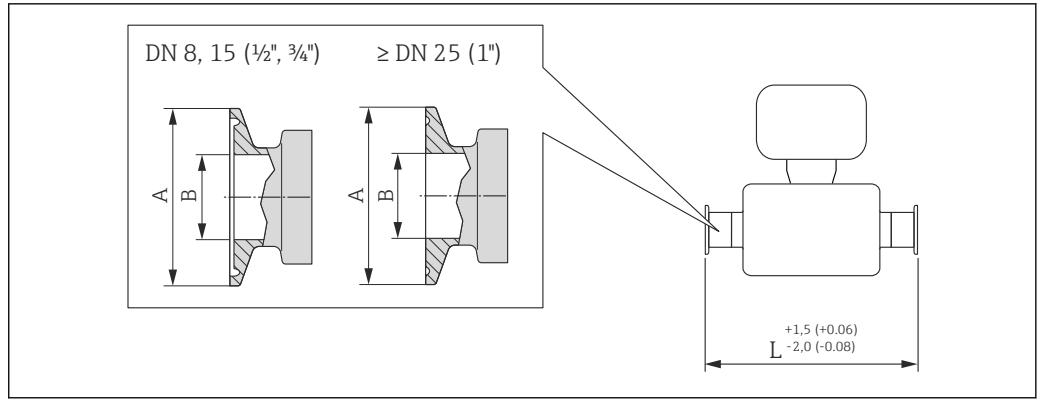

35 Engineering unit mm (in)

Tri-Clamp for pipe size: 1.4435 (316L) (order code for "Process connection", option FHW)

DN [mm]	Clamp [in]	A [mm]	B [mm]	L [mm]
8	1/2	25.0	9.40	362
15	3/4	25.0	15.75	466
25	1	50.4	22.10	606
40	1 1/2	50.4	34.80	818
50	2	63.9	47.50	1096

Available in Ra_{max}0.75 µm mechanically polished or Ra_{max}0.38 µm electropolished (order code for "Measuring tube material", option BB, BC, TD, TG)

1" Tri-Clamp: 1.4435 (316L) (order code for "Process connection", option FPW)

DN [mm]	Clamp [in]	A [mm]	B [mm]	L [mm]
8	1	50.4	22.10	362
15	1	50.4	22.10	466

Available in Ra_{max}0.75 µm mechanically polished or Ra_{max}0.38 µm electropolished (order code for "Measuring tube material", option BB, BC, TD, TG)

3/4" Tri-Clamp: 1.4435 (316L) (order code for "Process connection", option FFW)

DN [mm]	Clamp [in]	A [mm]	B [mm]	L [mm]
8	3/4	25.0	15.75	362

Available in Ra_{max}0.75 µm mechanically polished or Ra_{max}0.38 µm electropolished (order code for "Measuring tube material", option BB, BC, TD, TG)


1/2" Tri-Clamp: 1.4435 (316L) (order code for "Process connection", option FCW)

DN [mm]	Clamp [in]	A [mm]	B [mm]	L [mm]
15	1/2	25.0	9.40	466

Available in Ra_{max}0.75 µm mechanically polished or Ra_{max}0.38 µm electropolished (order code for "Measuring tube material", option BB, BC, TD, TG)

Tri-Clamp, eccentric (BS4825, ASME BPE, DIN 11866 line C)

All Tri-Clamp connections (BS4825-3, DN 1", 1½", 2") in accordance with hygienic clamp dimensions as per ASME BPE.

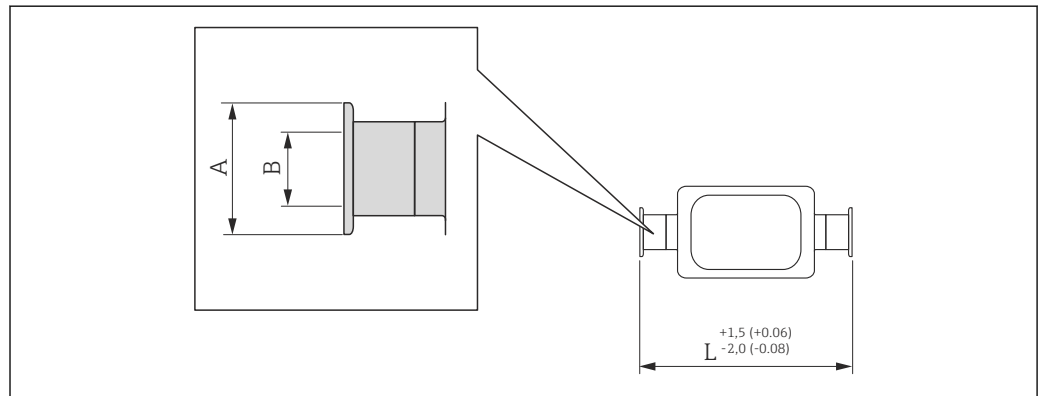


A0016543

36 Engineering unit mm (in)

Tri-Clamp, eccentric: 1.4435 (316L)					
DN [mm]	Order code for "Process connection", option	Clamp [in]	A [mm]	B [mm]	L [mm]
8	FEB	½	25.0	9.40	362
15	FED	¾	25.0	15.75	466
25	FEF	1	50.4	22.10	606
40	FEH	1½	50.4	34.80	825
50	FEK	2	63.9	47.50	1103

Available in Ra_{max}0.75 µm mechanically polished or Ra_{max}0.38 µm electropolished (order code for "Measuring tube material", option BB, BC, TD, TG)
 Additional information on "Eccentric clamps" (→ 39)

DIN 11851 (threaded hygienic connection), DIN 11866 line A

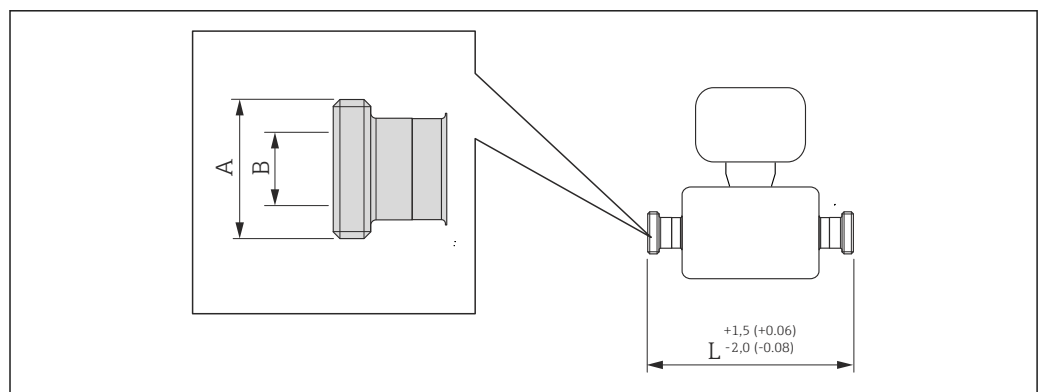

A0015628

37 Engineering unit mm (in)

Threaded hygienic connection DIN 11851: 1.4435 (316L) (order code for "Process connection", option KCW)

DN [mm]	A [in]	B [mm]	L [mm]
8	Rd 34 × 1/8	16	362
15	Rd 34 × 1/8	16	466
25	Rd 52 × 1/6	26	606
40	Rd 65 × 1/6	38	825
50	Rd 78 × 1/6	50	1 107


Available in Ra_{max}0.75 μm mechanically polished or Ra_{max}0.38 μm electropolished (order code for "Measuring tube material", option BB, BC, TD, TG)

Rd 28 × 1/8" - Threaded hygienic connection DIN 11851: 1.4435 (316L) (order code for "Process connection", option KAW)

DN [mm]	A [in]	B [mm]	L [mm]
8	Rd 28 × 1/8	10.00	362
15	Rd 28 × 1/8	10.00	466

Available in Ra_{max}0.75 μm mechanically polished or Ra_{max}0.38 μm electropolished (order code for "Measuring tube material", option BB, BC, TD, TG)

DIN 11864-1 Form A (aseptic threaded hygienic connection), DIN 11866 line A

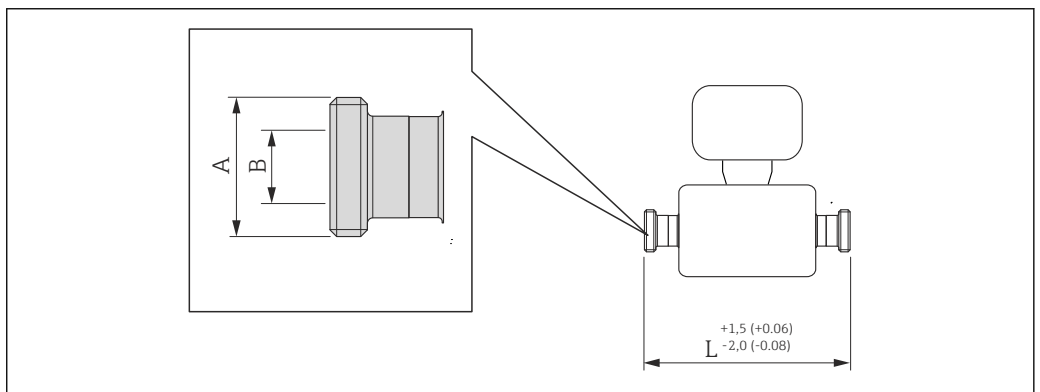


38 Engineering unit mm (in)

Threaded hygienic connection DIN 11864-1 Form A: 1.4435 (316L) (order code for "Process connection", option KGW)

DN [mm]	A [in]	B [mm]	L [mm]
8	Rd 34 × 1/8	16	362
15	Rd 34 × 1/8	16	466
25	Rd 52 × 1/6	26	620
40	Rd 65 × 1/6	38	825
50	Rd 78 × 1/6	50	1 107

Available in Ra_{max}0.75 μm mechanically polished or Ra_{max}0.38 μm electropolished (order code for "Measuring tube material", option BB, BC, TD, TG)

DIN 11864-2 Form A (aseptic flange), DIN 11866 line A

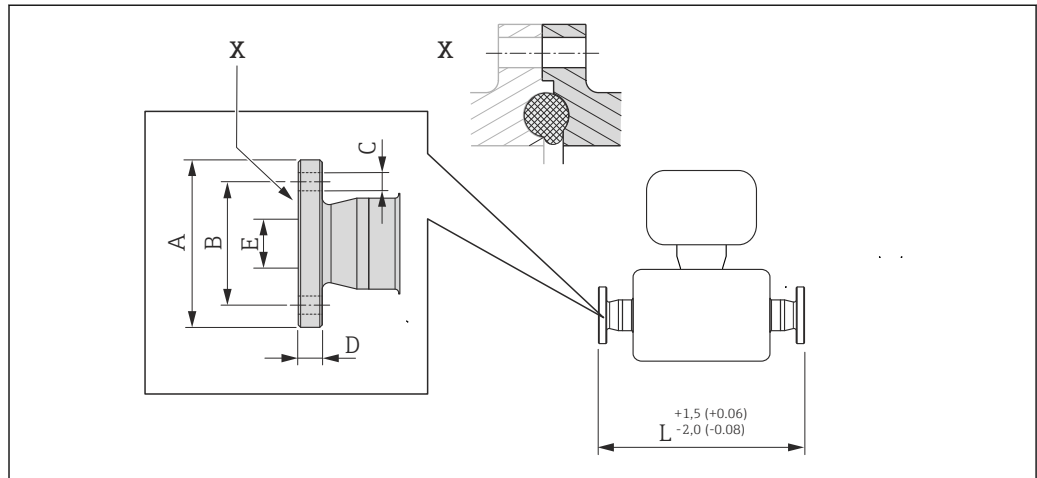

A0015627


39 Detail X: Asymmetrical process connection; the part shown in gray is provided by the supplier. Engineering unit mm (in).

DIN 11864-2 Form A (flange with groove): 1.4435 (316L) (order code for "Process connection", option KJW)

DN [mm]	A [mm]	B [mm]	C [mm]	D [mm]	E [mm]	L [mm]
8	59.0	42	4 × Ø9	10	16.00	384
15	59.0	42	4 × Ø9	10	16.00	488
25	70.0	53	4 × Ø9	10	26.00	626
40	82.0	65	4 × Ø9	10	38.00	840
50	94.0	77	4 × Ø9	10	50.00	1120

Available in Ra_{max}0.76 µm mechanically polished or Ra_{max}0.38 µm electropolished (order code for "Measuring tube material", option BB, BC, TD, TG)

DIN 11864-3 Form A (aseptic clamp with notch), DIN 11866 line A

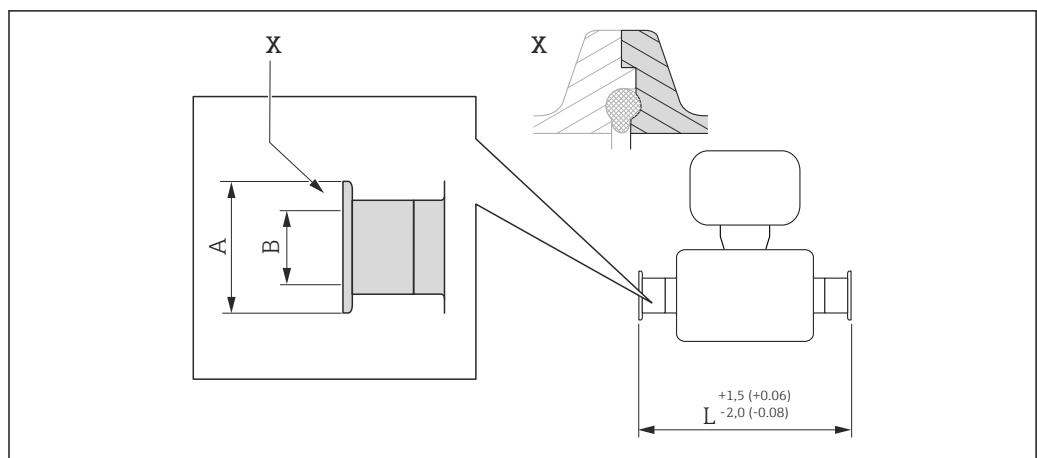


A0016905

40 Detail X: Asymmetrical process connection; the part shown in gray is provided by the supplier. Engineering unit mm (in).

DIN 11864-3 Form A: 1.4435 (316L) (order code for "Process connection", option KLV)			
DN [mm]	A [mm]	B [mm]	L [mm]
8	34.0	16.05	370
15	34.0	16.05	474
25	50.5	26.05	614
40	64.0	38.05	825
50	77.5	50.05	1096

Available in Ra_{max}0.76 µm mechanically polished or Ra_{max}0.38 µm electropolished (order code for "Measuring tube material", option BB, BC, TD, TG)

DIN 11864-3 Form A (aseptic clamp with notch, eccentric), DIN 11866 line A

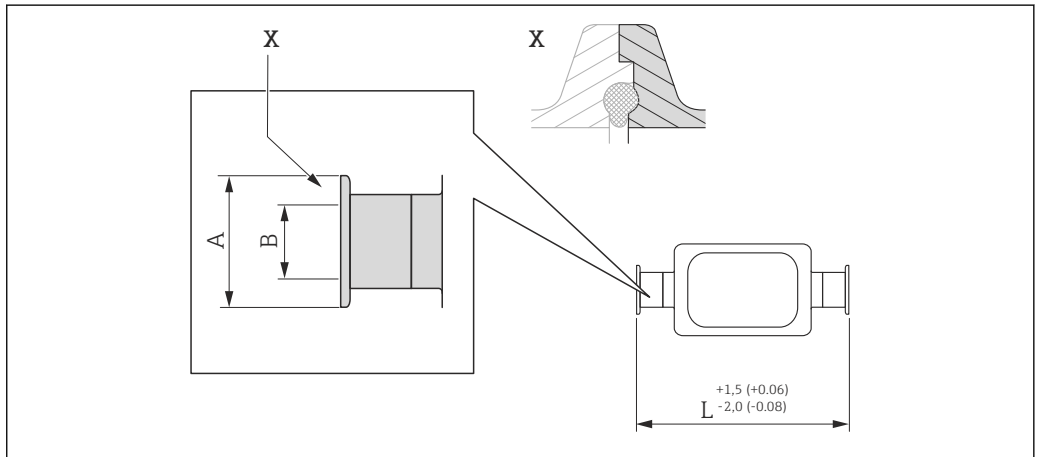


A0016904

41 Detail X: Asymmetrical process connection; the part shown in gray is provided by the supplier. Engineering unit mm (in).

DIN 11864-3 Form A, eccentric: 1.4435 (316L) (order code for "Process connection", option KNW)			
DN [mm]	A [mm]	B [mm]	L [mm]
8	34.0	10.00	370
15	34.0	16.00	474
25	50.5	26.00	624
50	77.5	50.00	1112

Available in Ra_{max}0.76 µm mechanically polished or Ra_{max}0.38 µm electropolished (order code for "Measuring tube material", option BB, BC, TD, TG)
 Additional information on "Eccentric clamps" (→ 39)

DIN 32676 (clamp connection), DIN 11866 line A

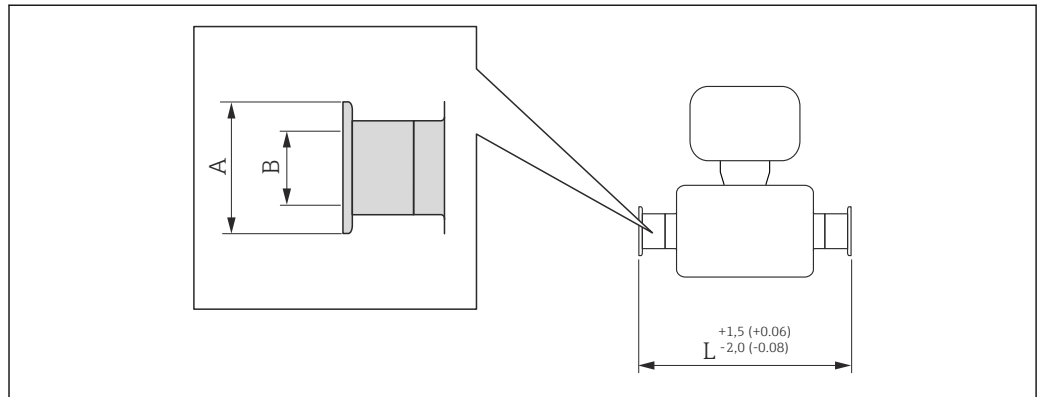


A0015625

42 Engineering unit mm (in)

Clamp connection DIN 32676: 1.4435 (316L) (order code for "Process connection", option KPW)			
DN [mm]	A [mm]	B [mm]	L [mm]
8	34.0	16.00	362
15	34.0	16.00	466
25	50.5	26.00	606
40	50.5	38.00	819
50	64.0	50.00	1097

Available in Ra_{max}0.76 µm mechanically polished or Ra_{max}0.38 µm electropolished (order code for "Measuring tube material", option BB, BC, TD, TG)

DIN 32676 (clamp connection, eccentric), DIN 11866 line A

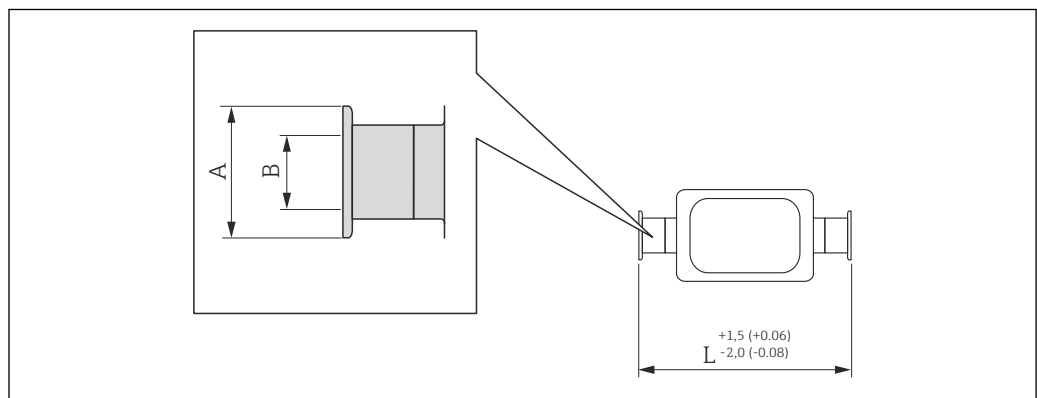


A0016543

43 Engineering unit mm (in)

Eccentric clamp connection DIN 32676: 1.4435 (316L) (order code for "Process connection", option KRW)			
DN [mm]	A [mm]	B [mm]	L [mm]
8	34.0	10.00	362
15	34.0	16.00	466
25	50.5	26.00	606

Eccentric clamp connection DIN 32676: 1.4435 (316L) (order code for "Process connection", option KRW)			
DN [mm]	A [mm]	B [mm]	L [mm]
40	64.0	50.00	1103
50			

ISO 2852 (clamp connection), ISO 2037

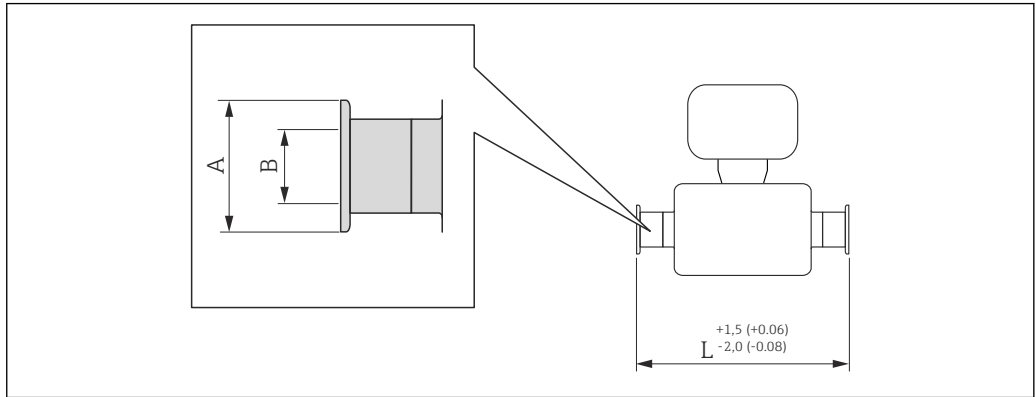


A0015625

44 Engineering unit mm (in)

Clamp connection ISO 2852, ISO 2037: 1.4435 (316L) (order code for "Process connection", option JSA)			
DN [mm]	A [mm]	B [mm]	L [mm]
8	50.5	22.6	362
15	50.5	22.6	466
25	50.5	22.6	606
40	50.5	35.6	818
50	64.0	48.6	1096

Available in Ra_{max}0.76 µm mechanically polished or Ra_{max}0.38 µm electropolished (order code for "Measuring tube material", option BB, BC, TD, TG)

ISO 2852 (clamp connection), DIN11866 line B

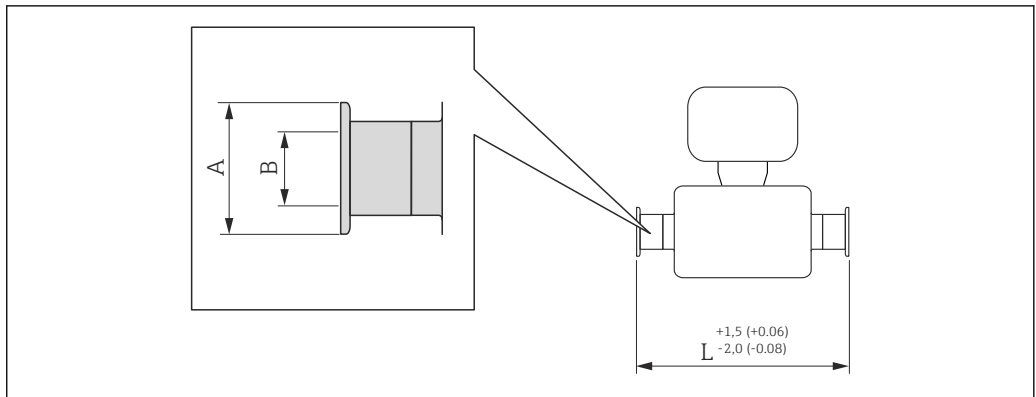


A0015625

45 Engineering unit mm (in)

Clamp connection ISO 2852, DIN11866 line B: 1.4435 (316L) (order code for "Process connection", option JSC)			
DN [mm]	A [mm]	B [mm]	L [mm]
8	34.0	14.00	362
15	34.0	18.10	466
25	50.5	29.70	606
40	64.0	44.30	818
50	77.5	56.30	1096

Available in Ra_{max}0.76 µm mechanically polished or Ra_{max}0.38 µm electropolished (order code for "Measuring tube material", option BB, BC, TD, TG)

ISO 2852 (clamp connection, eccentric), DIN11866 line B

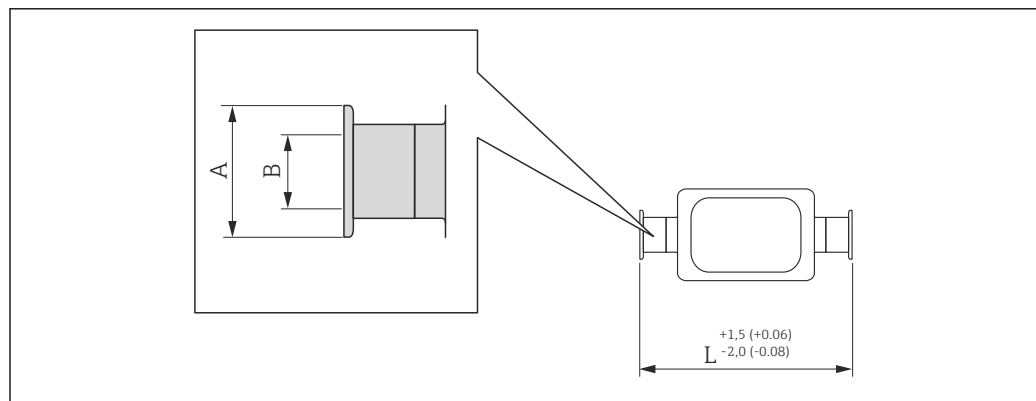


A0016543

46 Engineering unit mm (in)

Eccentric clamp connection ISO 2852, DIN11866 line B: 1.4435 (316L) (order code for "Process connection", option JEC)			
DN [mm]	A [mm]	B [mm]	L [mm]
8	34.0	10.30	362
15	34.0	14.00	466
25	34.0	18.10	606
40	50.5	29.70	825
50	64.0	44.30	1096

Available in Ra_{max}0.76 µm mechanically polished or Ra_{max}0.38 µm electropolished (order code for "Measuring tube material", option BB, BC, TD, TG)
Additional information on "Eccentric clamps" (→ 39)

ISO 2852 (clamp connection, eccentric), DIN11866 line B, for connecting to DN15 pipes

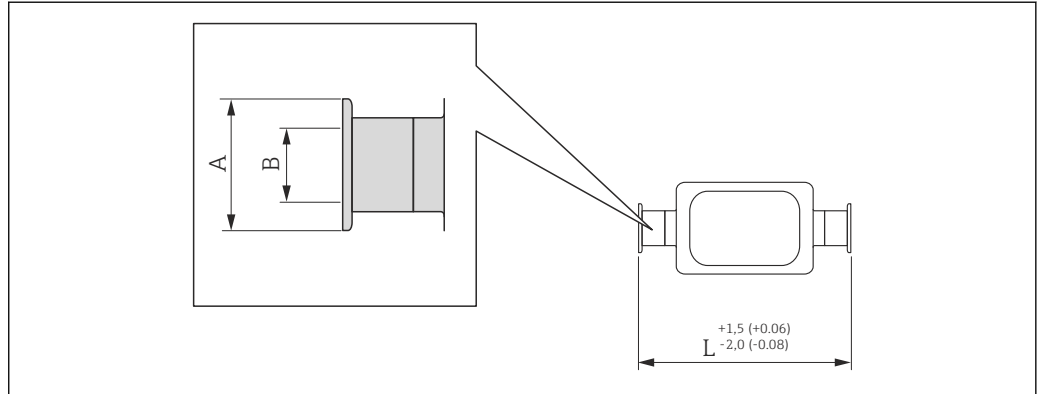

A0016543


47 Engineering unit mm (in)

Eccentric clamp connection ISO 2852, DIN11866 line B, for connecting to DN15 pipes: 1.4435 (316L) (order code for "Process connection", option JED)

DN [mm]	A [mm]	B [mm]	L [mm]
25	50.5	18.10	606

Available in Ra_{max}0.76 µm mechanically polished or Ra_{max}0.38 µm electropolished (order code for "Measuring tube material", option BB, BC, TD, TG)
 Additional information on "Eccentric clamps" (→ 39)

ISO 2853 (threaded hygienic connection), ISO 2037

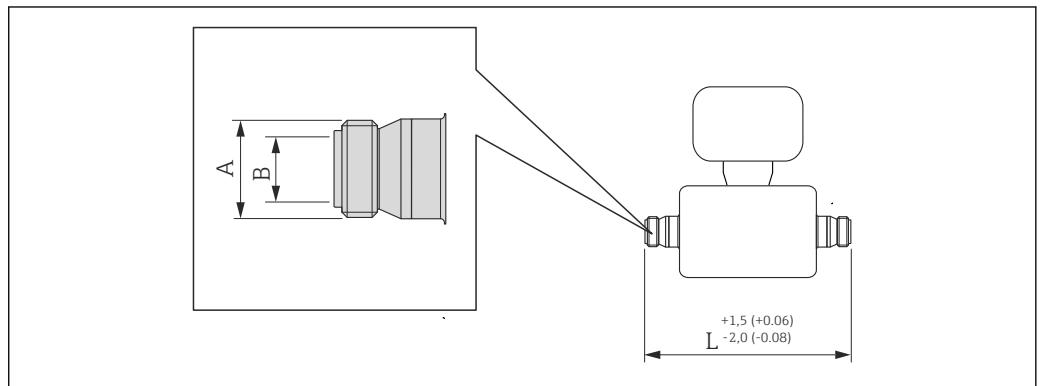

A0015623


48 Engineering unit mm (in)

Threaded hygienic connection ISO 2853: 1.4435 (316L) (order code for "Process connection", option JSD)

DN [mm]	A [mm]	B [mm]	L [mm]
8	37.13	22.60	370
15	37.13	22.60	474
25	37.13	22.60	614
40	50.65	35.60	829
50	64.10	48.60	1107

Available in Ra_{max}0.76 µm mechanically polished or Ra_{max}0.38 µm electropolished (order code for "Measuring tube material", option BB, BC, TD, TG)

Neumo BioConnect (clamp with notch), DIN 11866 line A

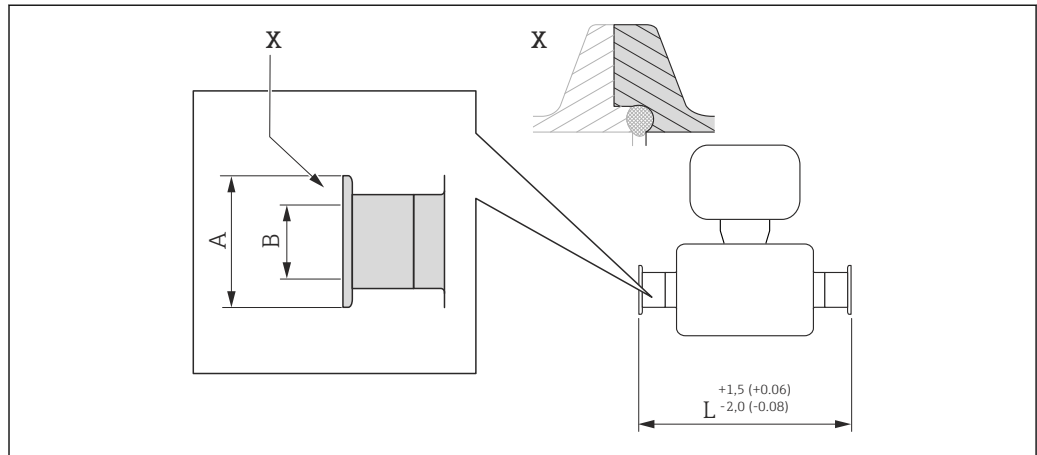


A0016905

49 Detail X: Asymmetrical process connection; the part shown in gray is provided by the supplier. Engineering unit mm (in).

Clamp connection Neumo BioConnect: 1.4435 (316L) (order code for "Process connection", option BSA)			
DN [mm]	A [mm]	B [mm]	L [mm]
8	25.0	10.00	362
15	25.0	16.00	466
25	50.4	26.00	606
40	64.0	38.00	819
50	77.4	50.00	1097

Available in Ra_{max}0.76 µm mechanically polished or Ra_{max}0.38 µm electropolished (order code for "Measuring tube material", option BB, BC, TD, TG)

Neumo BioConnect (clamp with notch, eccentric), DIN 11866 line A

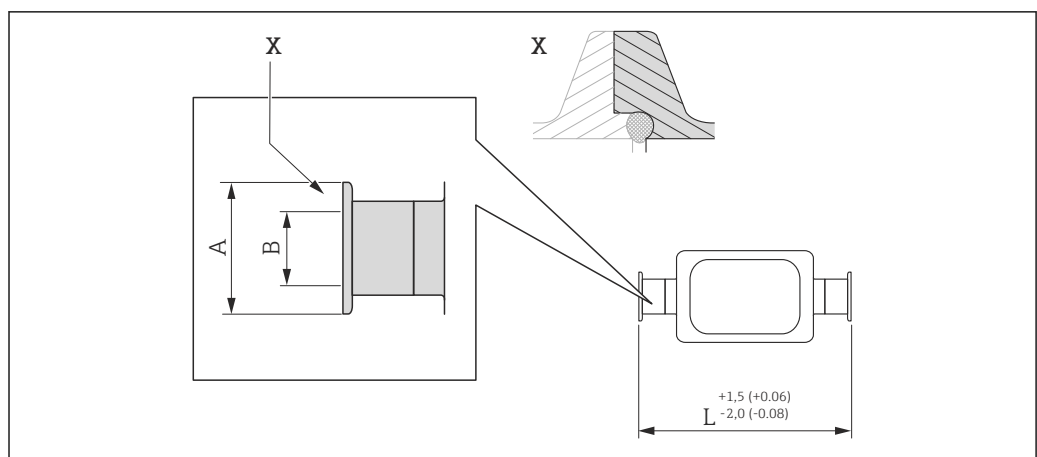


A0016906

50 Detail X: Asymmetrical process connection; the part shown in gray is provided by the supplier. Engineering unit mm (in).

Eccentric clamp connection Neumo BioConnect: 1.4435 (316L) (order code for "Process connection", option BEA)			
DN [mm]	A [mm]	B [mm]	L [mm]
8	25	10.00	362
15	25	16.00	466
25	25	26.00	610
50	25	50.00	1102

Available in Ra_{max}0.76 µm mechanically polished or Ra_{max}0.38 µm electropolished (order code for "Measuring tube material", option BB, BC, TD, TG)
 Additional information on "Eccentric clamps" (→ 39)

Neumo BioConnect (flange), DIN 11866 line A

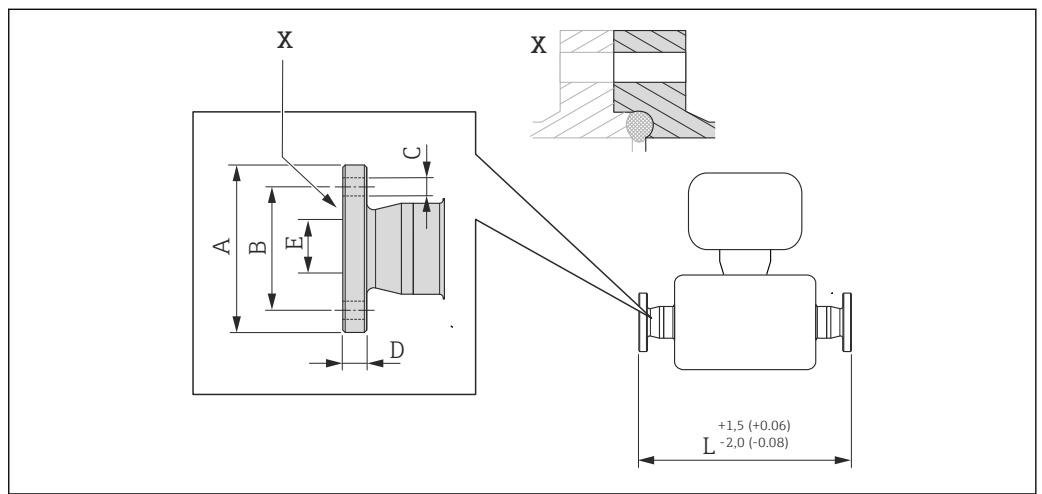


A0016907

51 Detail X: Asymmetrical process connection; the part shown in gray is provided by the supplier. Engineering unit mm (in).

Flange, Neumo BioConnect: 1.4435 (316L) (order code for "Process connection", option BSB)						
DN [mm]	A [mm]	B [mm]	C [mm]	D [mm]	E [mm]	L [mm]
8	65	45	4 × Ø9	10	10.00	384
15	75	55	4 × Ø9	10	16.00	488
25	85	65	4 × Ø9	12	26.00	626
40	100	80	4 × Ø9	12	38.00	840
50	110	90	4 × Ø9	14	50.00	1120

Available in Ra_{max}0.76 µm mechanically polished or Ra_{max}0.38 µm electropolished (order code for "Measuring tube material", option BB, BC, TD, TG)

BBS (clamp with notch), DIN 11866 line A

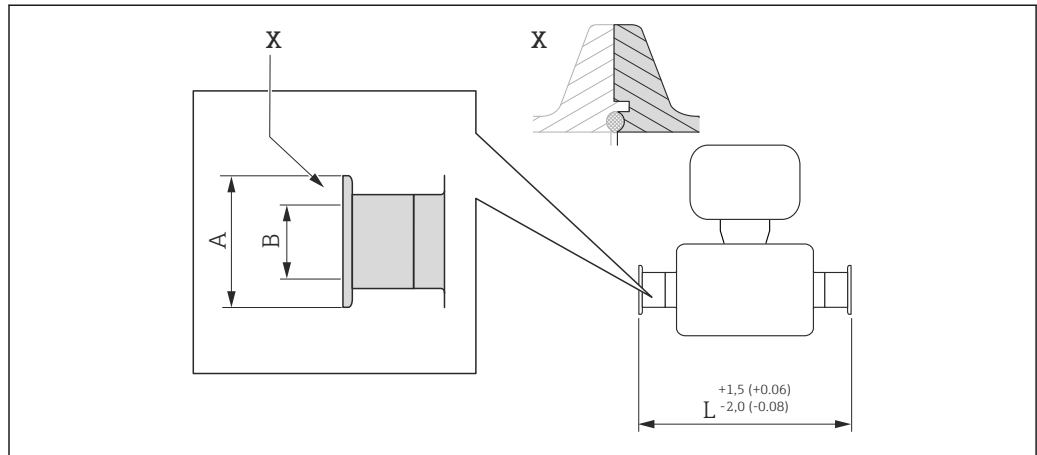


A0016908

52 Detail X: Asymmetrical process connection; the part shown in gray is provided by the supplier. Engineering unit mm (in).

Clamp connection BBS, DIN11866 line A: 1.4435 (316L) (order code for "Process connection", option BSE)			
DN [mm]	A [mm]	B [mm]	L [mm]
8	25.0	10.00	362
15	50.5	16.00	466
25	50.5	26.00	606
40	64.0	38.00	819
50	77.5	50.00	1097

Available in Ra_{max}0.76 µm mechanically polished or Ra_{max}0.38 µm electropolished (order code for "Measuring tube material", option BB, BC, TD, TG)

BBS (clamp with notch, eccentric), DIN 11866 line A

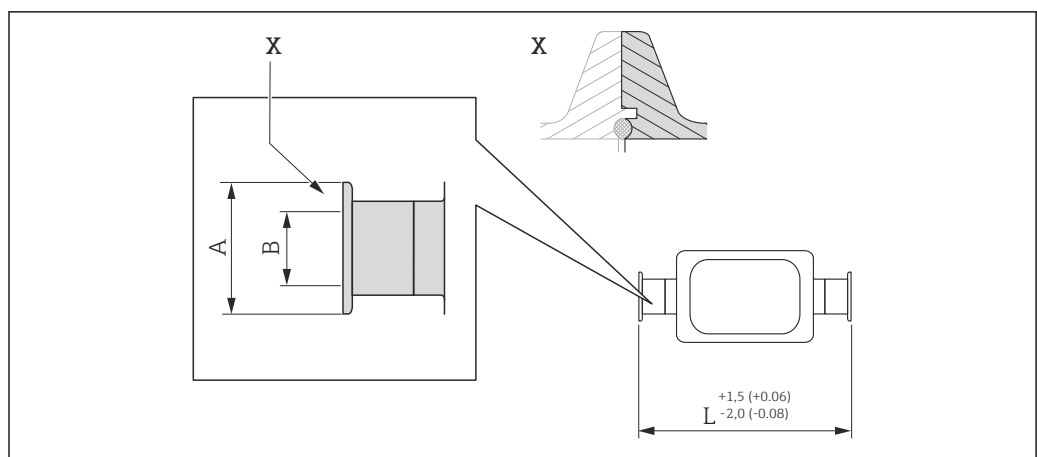


A0016909

53 Detail X: Asymmetrical process connection; the part shown in gray is provided by the supplier. Engineering unit mm (in).

Eccentric clamp connection BBS, DIN11866 line A: 1.4435 (316L) (order code for "Process connection", option BEJ)			
DN [mm]	A [mm]	B [mm]	L [mm]
8	25.0	10.00	362
15	50.5	16.00	466
25	50.5	26.00	606
50	77.5	50.00	1103

Available in Ra_{max}0.76 µm mechanically polished or Ra_{max}0.38 µm electropolished (order code for "Measuring tube material", option BB, BC, TD, TG)
 Additional information on "Eccentric clamps" (→ 39)

BBS (clamp with notch), DIN 11866 line B

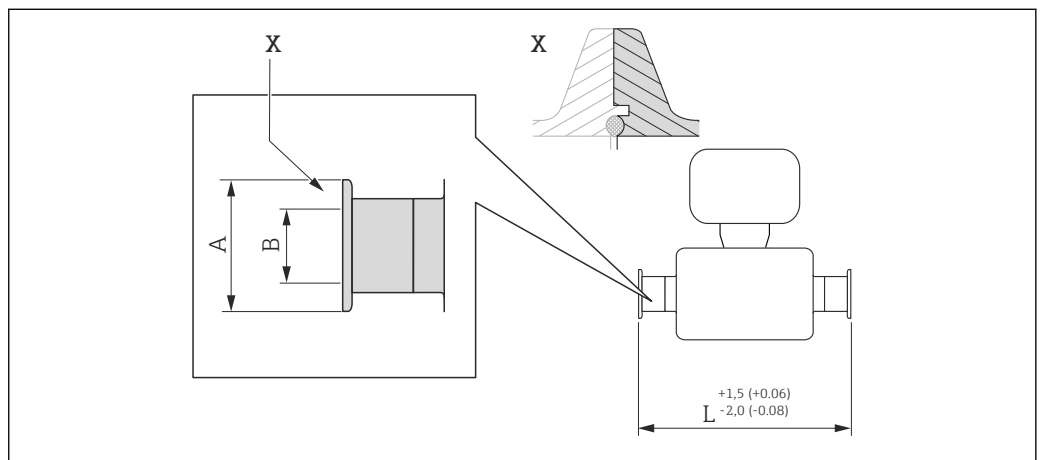


A0016908

54 Detail X: Asymmetrical process connection; the part shown in gray is provided by the supplier. Engineering unit mm (in).

Clamp connection BBS, DIN11866 line B: 1.4435 (316L) (order code for "Process connection", option BSJ)			
DN [mm]	A [mm]	B [mm]	L [mm]
8	25.0	14.00	362
15	50.5	18.10	466
25	50.5	29.70	606
40	64.0	44.30	825
50	77.5	56.30	1103

Available in Ra_{max}0.76 µm mechanically polished or Ra_{max}0.38 µm electropolished (order code for "Measuring tube material", option BB, BC, TD, TG)

BBS (clamp with notch, eccentric), DIN 11866 line B

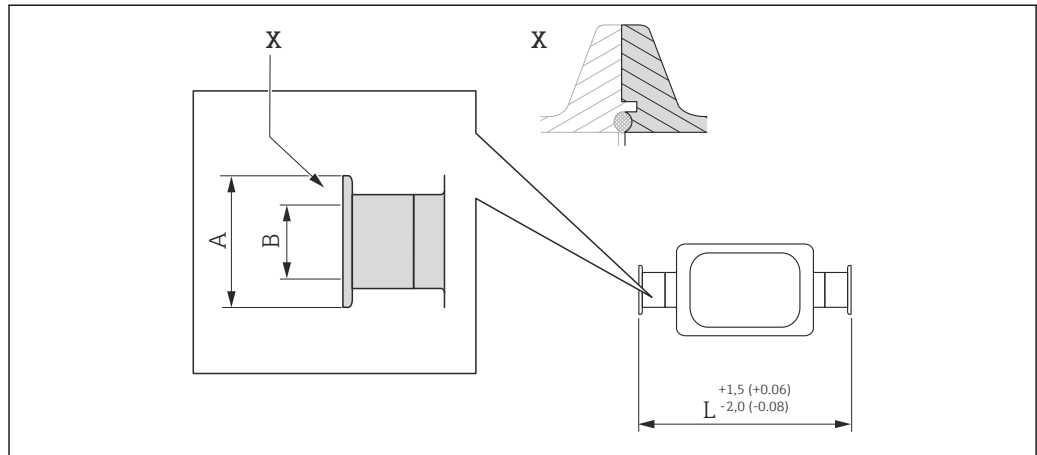


A0016909

55 Detail X: Asymmetrical process connection; the part shown in gray is provided by the supplier. Engineering unit mm (in).

Eccentric clamp connection BBS, DIN11866 line B: 1.4435 (316L) (order code for "Process connection", option BEK)			
DN [mm]	A [mm]	B [mm]	L [mm]
8	25.0	10.30	362
15	50.5	14.00	466
25	50.5	18.10	606
40	50.5	29.70	825
50	64.0	44.30	1103

Available in Ra_{max}0.76 μm mechanically polished or Ra_{max}0.38 μm electropolished (order code for "Measuring tube material", option BB, BC, TD, TG)
 Additional information on "Eccentric clamps" (→ 39)

BBS (flange), DIN 11866 line A

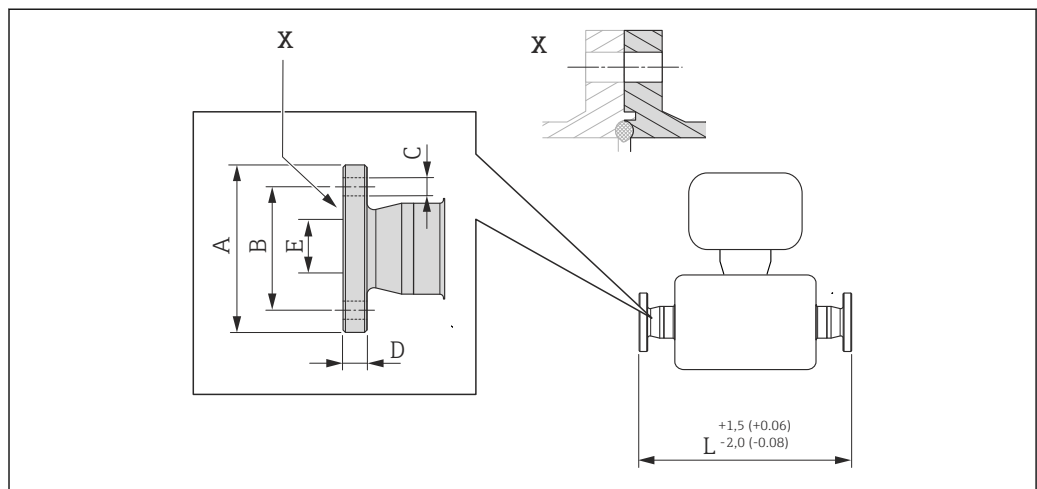


A0016910

56 Detail X: Asymmetrical process connection; the part shown in gray is provided by the supplier. Engineering unit mm (in).

Flange, BBS: 1.4435 (316L) (order code for "Process connection", option BSK)						
DN [mm]	A [mm]	B [mm]	C [mm]	D [mm]	E [mm]	L [mm]
8	59	42	4 × Ø9	10	10.00	384
15	59	42	4 × Ø9	10	16.00	488
25	70	53	4 × Ø9	10	26.00	626
40	82	65	4 × Ø9	10	38.00	840
50	94	77	4 × Ø9	10	50.00	1120

Available in Ra_{max}0.76 µm mechanically polished or Ra_{max}0.38 µm electropolished (order code for "Measuring tube material", option BB, BC, TD, TG)

BBS (flange), DIN 11866 line B

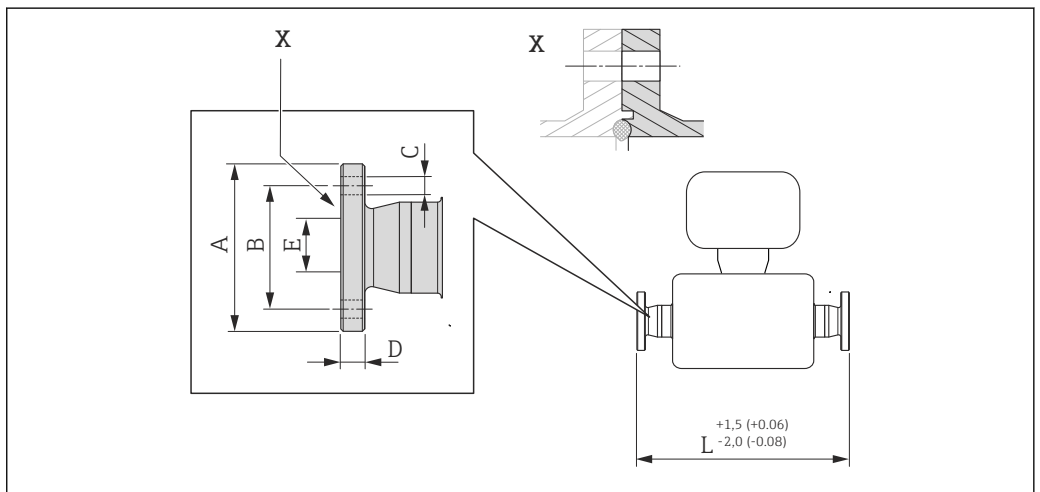


A0016910

57 Detail X: Asymmetrical process connection; the part shown in gray is provided by the supplier. Engineering unit mm (in).

Flange, BBS: 1.4435 (316L) (order code for "Process connection", option BSL)						
DN [mm]	A [mm]	B [mm]	C [mm]	D [mm]	E [mm]	L [mm]
8	59	42	4 × Ø9	10	14.00	384
15	62	45	4 × Ø9	10	18.10	488
25	74	57	4 × Ø9	10	29.70	626
40	88	71	4 × Ø9	10	44.30	840
50	103	85	4 × Ø9	10	56.30	1120

Available in Ra_{max}0.76 µm mechanically polished or Ra_{max}0.38 µm electropolished (order code for "Measuring tube material", option BB, BC, TD, TG)

BBS (aseptic threaded hygienic connection), DIN 11866 line A

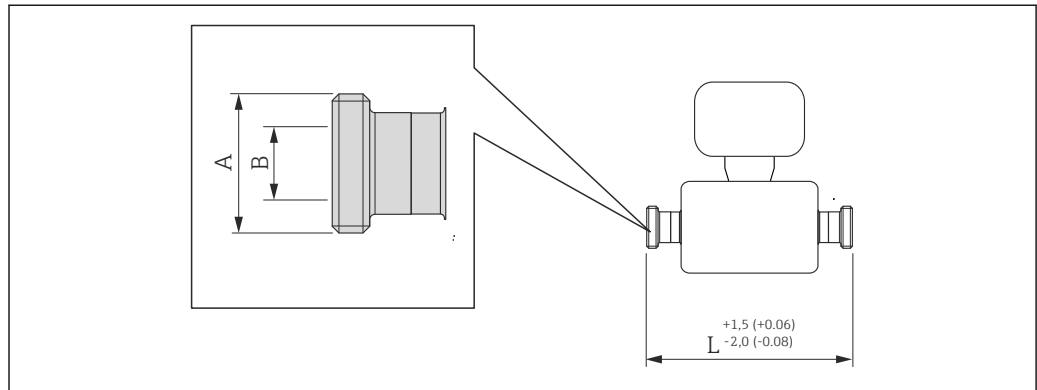


A0015628

58 Engineering unit mm (in)

Aseptic threaded hygienic connection BBS, DIN 11866 line A: 1.4435 (316L) (order code for "Process connection", option BSC)			
DN [mm]	A [in]	B [mm]	L [mm]
8	M22 × 1.5	10.00	362
15	M30 × 2	16.00	466
25	M42 × 2	26.00	606
40	M52 × 2	38.00	819
50	M68 × 2	50.00	1097

Available in Ra_{max}0.76 µm mechanically polished or Ra_{max}0.38 µm electropolished (order code for "Measuring tube material", option BB, BC, TD, TG)

BBS (aseptic threaded hygienic connection), DIN 11866 line B

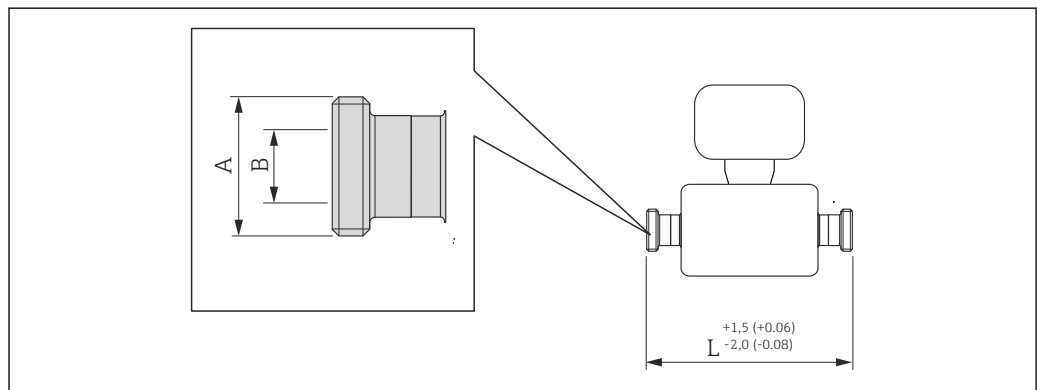

A0015628


59 Engineering unit mm (in)

Aseptic threaded hygienic connection BBS, DIN 11866 line B: 1.4435 (316L) (order code for "Process connection", option BSD)			
DN [mm]	A [in]	B [mm]	L [mm]
8	M26 × 1.5	10.00	362
15	M30 × 2	16.00	466
25	M42 × 2	26.00	606

Aseptic threaded hygienic connection BBS, DIN 11866 line B: 1.4435 (316L) (order code for "Process connection", option BSD)			
DN [mm]	A [in]	B [mm]	L [mm]
40	M56 × 2	38.00	829
50	M68 × 2	50.00	1 103

Available in Ra_{max}0.76 µm mechanically polished or Ra_{max}0.38 µm electropolished (order code for "Measuring tube material", option BB, BC, TD, TG)

SMS 1145 (threaded hygienic connection)

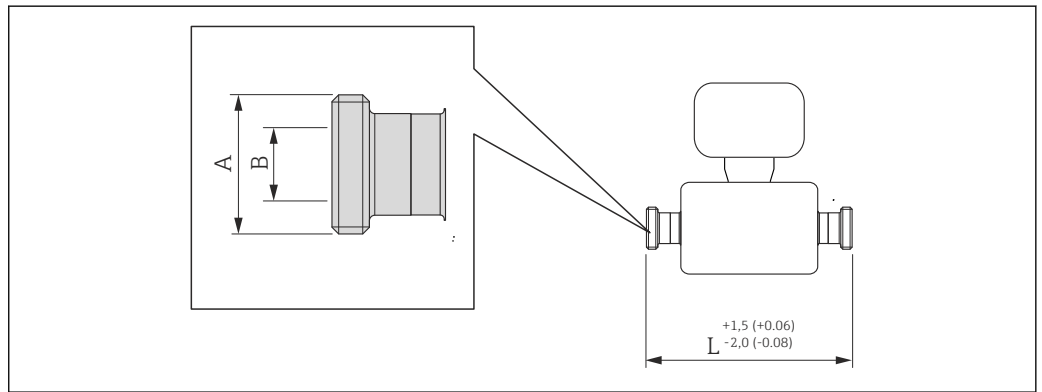

A0015628


60 Engineering unit mm (in)

Threaded hygienic connection SMS 1145: 1.4435 (316L) (order code for "Process connection", option SAW)			
DN [mm]	A [in]	B [mm]	L [mm]
8	Rd 40 × 1/6	22.50	362
15	Rd 40 × 1/6	22.50	466
25	Rd 40 × 1/6	22.50	606
40	Rd 60 × 1/6	35.50	829
50	Rd 70 × 1/6	48.50	1 107

Available in Ra_{max}0.76 µm mechanically polished or Ra_{max}0.38 µm electropolished (order code for "Measuring tube material", option BB, BC, TD, TG)

Process connections in US units

Flange connections ASME B16.5

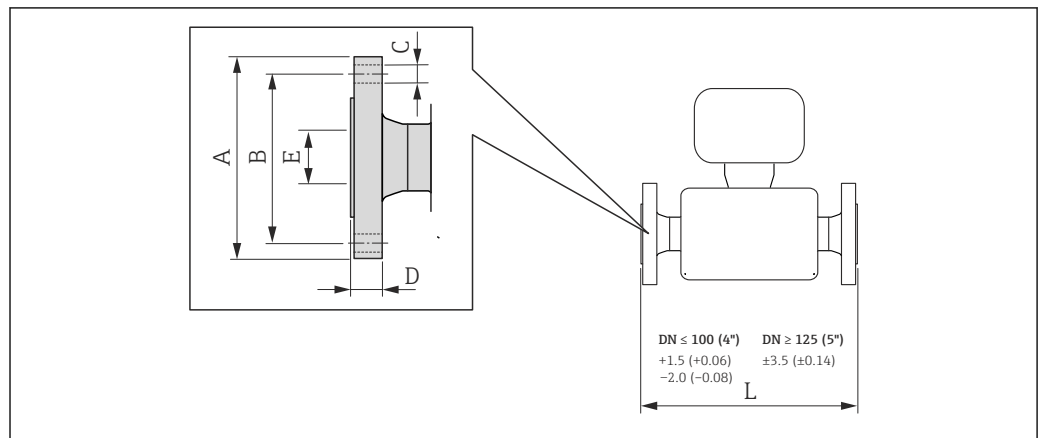

A0015621

61 Engineering unit mm (in)

Flange according to ASME B16.5, Cl 150: 1.4404 (316/316L) (order code for "Process connection", option AAW)

Surface roughness (flange): Ra 3.2 to 6.3 μm

DN [in]	A [in]	B [in]	C [in]	D [in]	E [in]	L [in]
$\frac{3}{8}$ ¹⁾	3.54	2.37	4 × Ø0.62	0.44	0.62	13.23
$\frac{1}{2}$	3.54	2.37	4 × Ø0.62	0.44	0.62	17.32
1	4.33	3.13	4 × Ø0.62	0.56	1.05	22.83
1½	4.92	3.87	4 × Ø0.62	0.69	1.61	31.26
2	5.91	4.75	4 × Ø0.75	0.75	2.07	42.17

1) DN $\frac{3}{8}$ " with DN $\frac{1}{2}$ " flanges as standard

Flange according to ASME B16.5, Cl 300: 1.4404 (316/316L) (order code for "Process connection", option ABW)


Surface roughness (flange): Ra 3.2 to 6.3 μm

DN [in]	A [in]	B [in]	C [in]	D [in]	E [in]	L [in]
$\frac{3}{8}$ ¹⁾	3.74	2.63	4 × Ø0.62	0.65	0.62	13.23
$\frac{1}{2}$	3.74	2.63	4 × Ø0.62	0.65	0.62	17.32
1	4.92	3.50	4 × Ø0.75	0.71	1.05	22.83
1½	6.10	4.50	4 × Ø0.88	0.97	1.61	31.26
2	6.50	5.00	8 × Ø0.75	1.09	2.07	42.17

1) DN $\frac{3}{8}$ " with DN $\frac{1}{2}$ " flanges as standard

Tri-Clamp (BS4825, ASME BPE, DIN 11866 line C)

All Tri-Clamp connections (BS4825-3, DN 1", 1½", 2") in accordance with hygienic clamp dimensions as per ASME BPE.

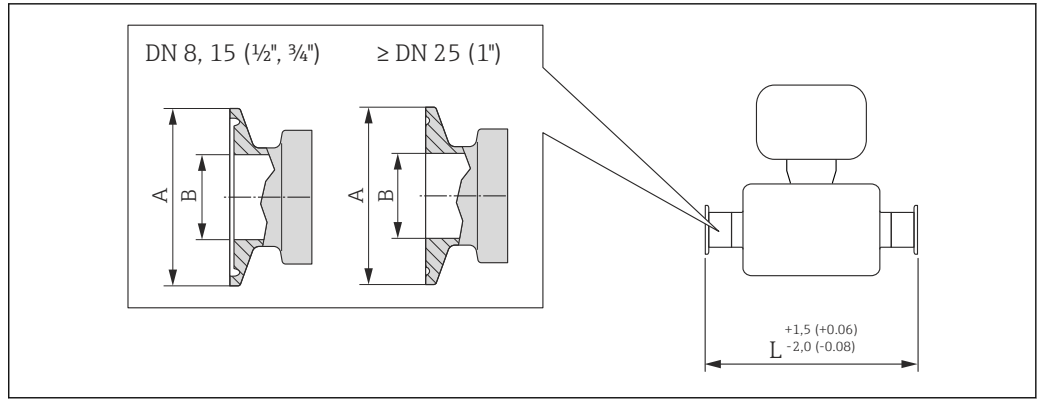

A0023942

62 Engineering unit mm (in)

Tri-Clamp for pipe size: 1.4435 (316L) (order code for "Process connection", option FHW)

DN [in]	Clamp [in]	A [in]	B [in]	L [in]
3/8	1/2	0.98	0.37	14.25
1/2	3/4	0.98	0.62	18.35
1	1	1.98	0.87	23.86
1 1/2	1 1/2	1.98	1.37	32.2
2	2	2.52	1.87	43.15

Available in Ra_{max}30 μm mechanically polished or Ra_{max}15 μm electropolished (order code for "Measuring tube material", option BB, BC, TD, TG)

1" Tri-Clamp: 1.4435 (316L) (order code for "Process connection", option FPW)

DN [in]	Clamp [in]	A [in]	B [in]	L [in]
3/8	1	1.98	0.87	14.25
1/2	1	1.98	0.87	18.35

Available in Ra_{max}30 μm mechanically polished or Ra_{max}15 μm electropolished (order code for "Measuring tube material", option BB, BC, TD, TG)

3/4" Tri-Clamp: 1.4435 (316L) (order code for "Process connection", option FFW)

DN [in]	Clamp [in]	A [in]	B [in]	L [in]
3/8	3/4	0.98	0.62	14.25

Available in Ra_{max}30 μm mechanically polished or Ra_{max}15 μm electropolished (order code for "Measuring tube material", option BB, BC, TD, TG)


1/2" Tri-Clamp: 1.4435 (316L) (order code for "Process connection", option FCW)

DN [in]	Clamp [in]	A [in]	B [in]	L [in]
1/2	1/2	0.98	0.37	18.35

Available in Ra_{max}30 μm mechanically polished or Ra_{max}15 μm electropolished (order code for "Measuring tube material", option BB, BC, TD, TG)

Tri-Clamp, eccentric (BS4825, ASME BPE, DIN 11866 line C)

All Tri-Clamp connections (BS4825-3, DN 1", 1½", 2") in accordance with hygienic clamp dimensions as per ASME BPE.

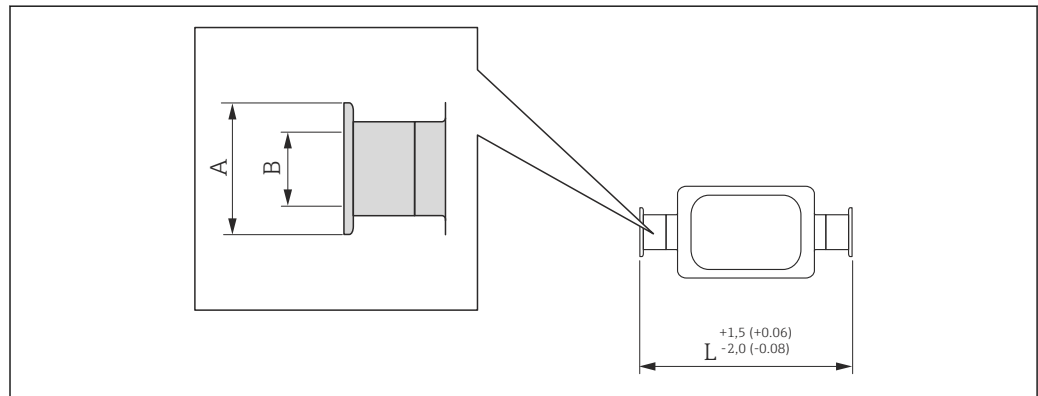

A0016543


63 Engineering unit mm (in)

Tri-Clamp, eccentric: 1.4435 (316L)					
DN [in]	Order code for "Process connection", option	Clamp [in]	A [in]	B [in]	L [in]
3/8	FEB	1/2	0.98	0.37	14.25
1/2	FED	3/4	0.98	0.62	18.35
1	FEF	1	1.98	0.87	23.86
1½	FEH	1½	1.98	1.37	32.48
2	FEK	2	2.52	1.87	43.43

Available in Ra_{max}30 µin mechanically polished or Ra_{max}15 µin electropolished (order code for "Measuring tube material", option BB, BC, TD, TG)

Additional information on "Eccentric clamps" (→ 39)

Neumo BioConnect (clamp with notch), DIN 11866 line A

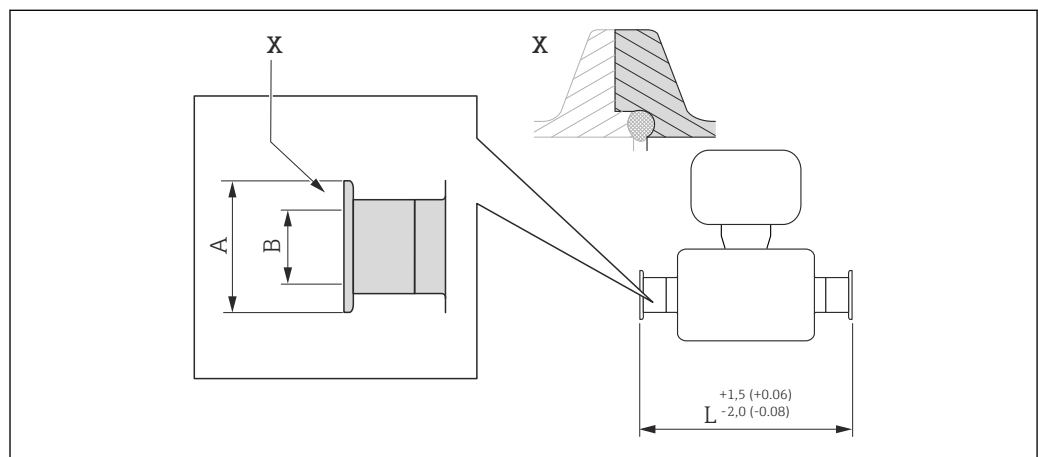


A0016905

64 Detail X: Asymmetrical process connection; the part shown in gray is provided by the supplier. Engineering unit mm (in).

Clamp connection Neumo BioConnect: 1.4435 (316L) (order code for "Process connection", option BSA)			
DN [in]	A [in]	B [in]	L [in]
3/8	0.98	0.39	14.25
1/2	0.98	0.63	18.35
1	1.98	1.02	23.86
1 1/2	2.52	1.5	32.24
2	3.05	1.97	43.19

Available in Ra_{max}30 µin mechanically polished or Ra_{max}15 µin electropolished (order code for "Measuring tube material", option BB, BC, TD, TG)

Neumo BioConnect (clamp with notch, eccentric), DIN 11866 line A

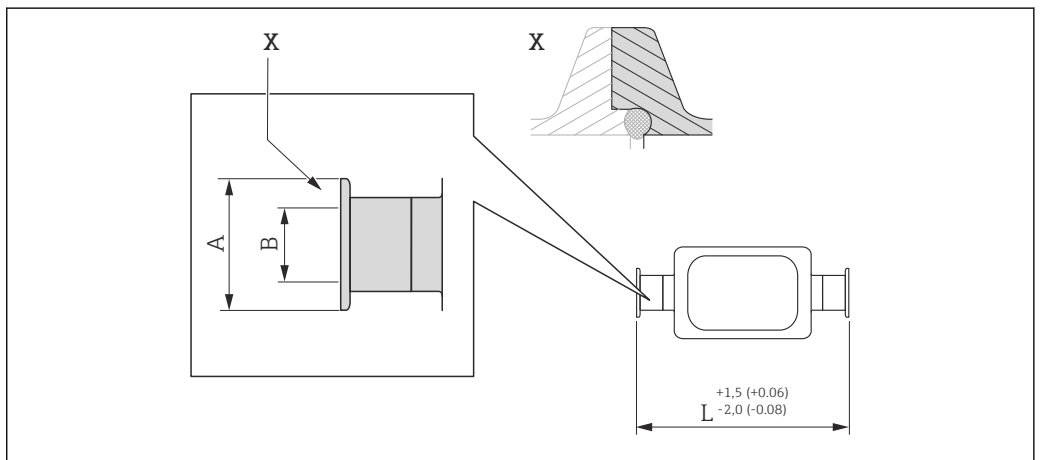


A0016906

65 Detail X: Asymmetrical process connection; the part shown in gray is provided by the supplier. Engineering unit mm (in).

Eccentric clamp connection Neumo BioConnect: 1.4435 (316L) (order code for "Process connection", option BEA)			
DN [in]	A [in]	B [in]	L [in]
3/8	0.98	0.39	14.25
1/2	0.98	0.63	18.35
1	0.98	1.02	24.02
1	0.98	1.97	43.39

Available in Ra_{max}30 µin mechanically polished or Ra_{max}15 µin electropolished (order code for "Measuring tube material", option BB, BC, TD, TG)
 Additional information on "Eccentric clamps" (→ 39)

Neumo BioConnect (flange), DIN 11866 line A

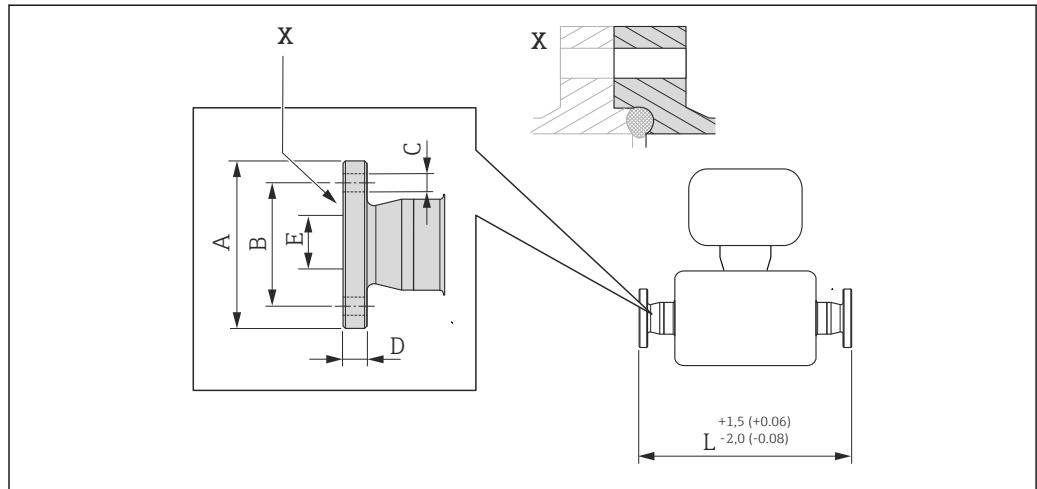


A0016907

66 Detail X: Asymmetrical process connection; the part shown in gray is provided by the supplier. Engineering unit mm (in).

Flange, Neumo BioConnect: 1.4435 (316L) (order code for "Process connection", option BSB)						
DN [in]	A [in]	B [in]	C [in]	D [in]	E [in]	L [in]
3/8	2.56	1.77	4 × Ø0.35	0.39	0.39	15.12
1/2	2.95	2.17	4 × Ø0.35	0.39	0.63	19.21
1	3.35	2.56	4 × Ø0.35	0.47	1.02	24.65
1 1/2	3.94	3.15	4 × Ø0.35	0.47	1.5	33.07
2	4.33	3.54	4 × Ø0.35	0.55	1.97	44.09

Available in Ra_{max}30 µin mechanically polished or Ra_{max}15 µin electropolished (order code for "Measuring tube material", option BB, BC, TD, TG)

SMS 1145 (threaded hygienic connection)

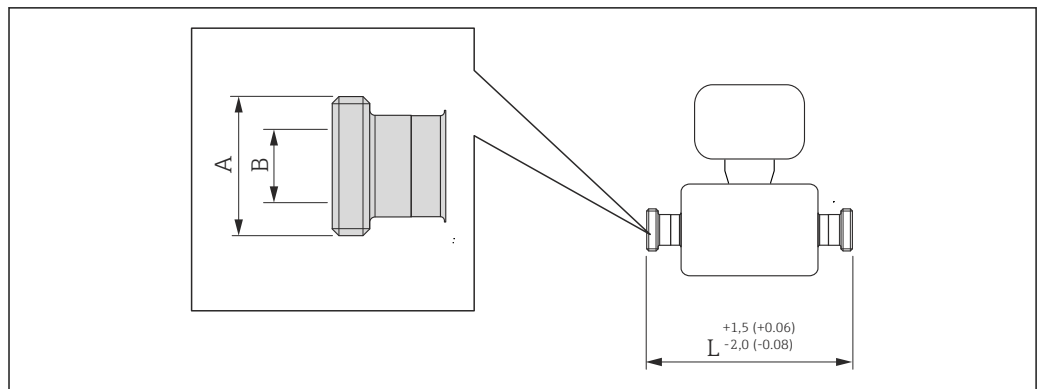

A0015628

67 Engineering unit mm (in)

Threaded hygienic connection SMS 1145: 1.4435 (316L) (order code for "Process connection", option SAW)


DN [in]	A [in]	B [in]	L [in]
3/8	Rd 40 × 1/6	0.89	14.25
1/2	Rd 40 × 1/6	0.89	18.35
1	Rd 40 × 1/6	0.89	23.86
1 1/2	Rd 60 × 1/6	1.4	32.64
2	Rd 70 × 1/6	1.91	43.58

Available in Ra_{max}30 µin mechanically polished or Ra_{max}15 µin electropolished (order code for "Measuring tube material", option BB, BC, TD, TG)

Safety Barrier Promass 100

Top-hat rail EN 60715:

- TH 35 x 7.5
- TH 35 x 15

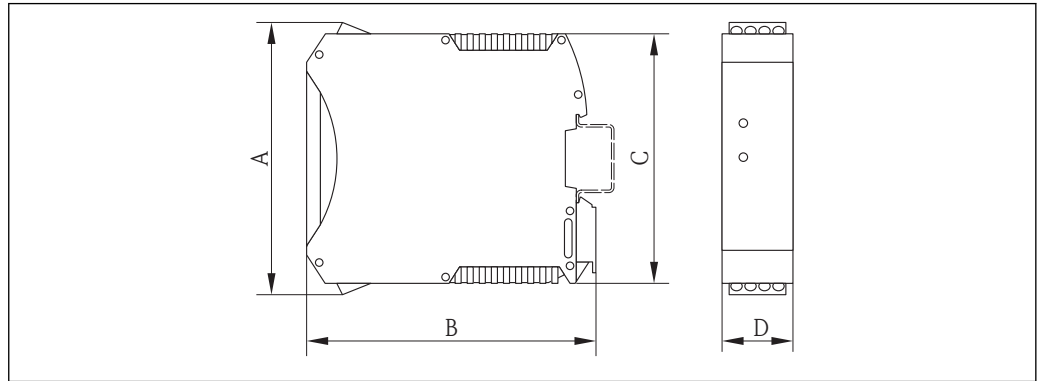


A0016777

A		B		C		D	
[mm]	[in]	[mm]	[in]	[mm]	[in]	[mm]	[in]
108	4.25	114.5	4.51	99	3.9	22.5	0.89

Accessories

Purge connections / secondary containment monitoring

Order code for "Sensor options", option CH

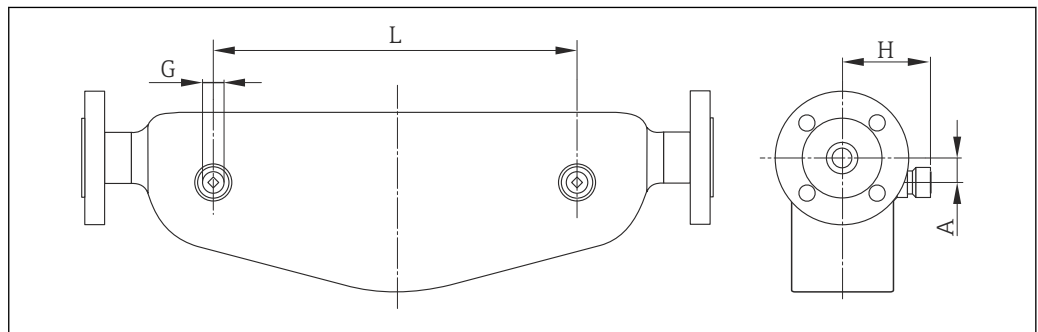

A0003288

DN		G	A		H		L	
[mm]	[in]	[in]	[mm]	[in]	[mm]	[in]	[mm]	[in]
8	3/8	1/2 NPT	25	0.98	82	3.23	110	4.34
15	1/2	1/2 NPT	25	0.98	82	3.23	204	8.04
25	1	1/2 NPT	25	0.98	82	3.23	348	13.54
40	1 1/2	1/2 NPT	45	1.77	102	4.02	526	20.70
50	2	1/2 NPT	58	2.28	119.5	4.70	763	30.04

Weight**Compact version***Weight in SI units*

All values (weight) refer to devices with EN/DIN PN 40 flanges. Weight information in [kg].

DN [mm]	Weight [kg]
8	11
15	13
25	19
40	41
50	78

Weight in US units

All values (weight) refer to devices with EN/DIN PN 40 flanges. Weight information in [lbs].

DN [in]	Weight [lbs]
3/8	24
1/2	29
1	42
1 1/2	90
2	172


Safety Barrier Promass 100

49 g (1.73 ounce)

Materials**Transmitter housing**

- Order code for "Housing", option **A** "Compact, aluminum coated":
Aluminum, AlSi10Mg, coated
- Order code for "Housing", option **B** "Compact, hygienic, stainless":
Hygienic version, stainless steel 1.4301 (304)
- Order code for "Housing", option **C** "Ultra compact, hygienic, stainless":
Hygienic version, stainless steel 1.4301 (304)

Cable entries/cable glands

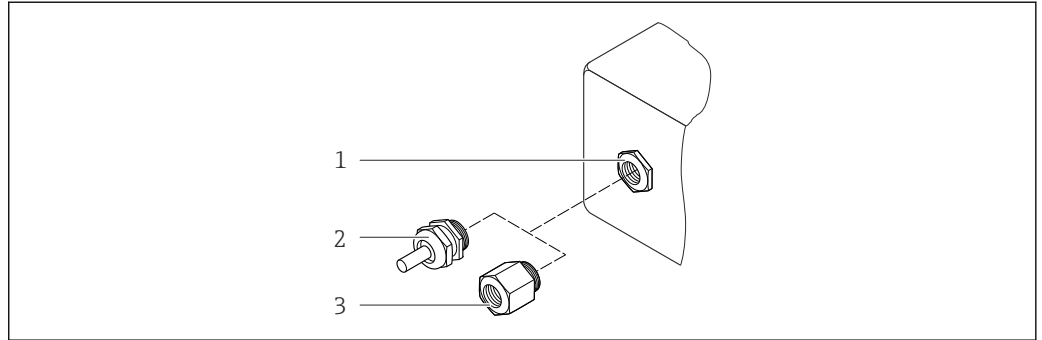

A0020640

▣ 68 Possible cable entries/cable glands

- 1 Cable entry in transmitter housing, wall-mount housing or connection housing with internal thread M20 x 1.5
- 2 Cable gland M20 x 1.5
- 3 Adapter for cable entry with internal thread G 1/2" or NPT 1/2"

Order code for "Housing", option A "Compact, coated aluminum"

The various cable entries are suitable for hazardous and non-hazardous areas.

Cable entry/cable gland	Material
Cable gland M20 x 1.5	Nickel-plated brass
Adapter for cable entry with internal thread G 1/2"	
Adapter for cable entry with internal thread NPT 1/2"	

Order code for "Housing", option B "Compact, hygienic, stainless"

The various cable entries are suitable for hazardous and non-hazardous areas.

Cable entry/cable gland	Material
Cable gland M20 x 1.5	Stainless steel, 1.4404 (316L)
Adapter for cable entry with internal thread G 1/2"	
Adapter for cable entry with internal thread NPT 1/2"	

Device plug

Electrical connection	Material
Plug M12x1	<ul style="list-style-type: none"> ▪ Socket: Stainless steel, 1.4404 (316L) ▪ Contact housing: Polyamide ▪ Contacts: Gold-plated brass

Sensor housing

- Acid and alkali-resistant outer surface
- Stainless steel 1.4301 (304)

Measuring tubes

Stainless steel, 1.4435 (316L)

Surface quality (wetted parts)

- Ra_{max} = 0.76 µm (32 µin) mechanically polished
- Ra_{max} = 0.38 µm (16 µin) electropolished
- Delta ferrite <1%

Process connections

- Flanges according to EN 1092-1 (DIN 2501) / according to ASME B16.5 / according to JIS B2220: Stainless steel, 1.4404 (316/316L)
- All other process connections: Stainless steel, 1.4435 (316L)



List of all available process connections (→ 82)

Seals

Welded process connections without internal seals

Safety Barrier Promass 100

Housing: Polyamide

Process connections

- Flanges:
 - EN 1092-1 (DIN 2501)
 - EN 1092-1 (DIN 2512N)
 - ASME B16.5
 - JIS B2220
- Tri-Clamp (OD tubes)
- Clamp with aseptic connection
DIN 11864-3, Form A
- Clamp with compression fitting
 - DIN 32676
 - ISO 2852
 - BioConnect
- Clamp (eccentric) :
 - Tri-Clamp
 - DIN 11864-3 Form A
 - DIN 32676
 - ISO 2852
 - Neumo BioConnect
 - BBS
- Threaded hygienic connection:
 - DIN 11851
 - SMS 1145
 - ISO 2853
 - DIN 11864-1 Form A
- Flange:
 - DIN 11864-2 Form A



For information on the materials of the process connections (→ 80)

Operability**Operating concept****Operator-oriented menu structure for user-specific tasks**

- Commissioning
- Operation
- Diagnostics
- Expert level

Quick and safe commissioning

- Individual menus for applications
- Menu guidance with brief explanations of the individual parameter functions

Reliable operation

- Operation in the following languages:
 - Via "FieldCare" operating tool: English, German, French, Spanish, Italian, Chinese, Japanese
 - Via Web browser: English, German, French, Spanish, Italian, Dutch, Portuguese, Polish, Russian, Turkish, Chinese, Japanese, Bahasa (Indonesian), Vietnamese, Czech
- Uniform operating philosophy applied to operating tools and Web browser
- If replacing the electronic module, transfer the device configuration via the plug-in memory (HistoROM DAT) which contains the process and measuring device data and the event logbook. No need to reconfigure. For devices with Modbus RS485, the data recovery function is implemented without the plug-in memory (HistoROM DAT).

Efficient diagnostics increase measurement availability

- Troubleshooting measures can be called up via the operating tools and Web browser
- Diverse simulation options
- Status indicated by several light emitting diodes (LEDs) on the electronic module in the housing compartment

Local display

The local display is only available with the following device version:
 Order code for "Display; Operation", option **B**: 4-line; via communication


Display element

- 4-line liquid crystal display with 16 characters per line.
- White background lighting; switches to red in event of device errors.
- Format for displaying measured variables and status variables can be individually configured.
- Permitted ambient temperature for the display: -20 to +60 °C (-4 to +140 °F). The readability of the display may be impaired at temperatures outside the temperature range.

Remote operation

Via HART protocol

This communication interface is present in the following device version:
 Order code for "Output", option **B**: 4-20 mA HART, pulse/frequency/switch output

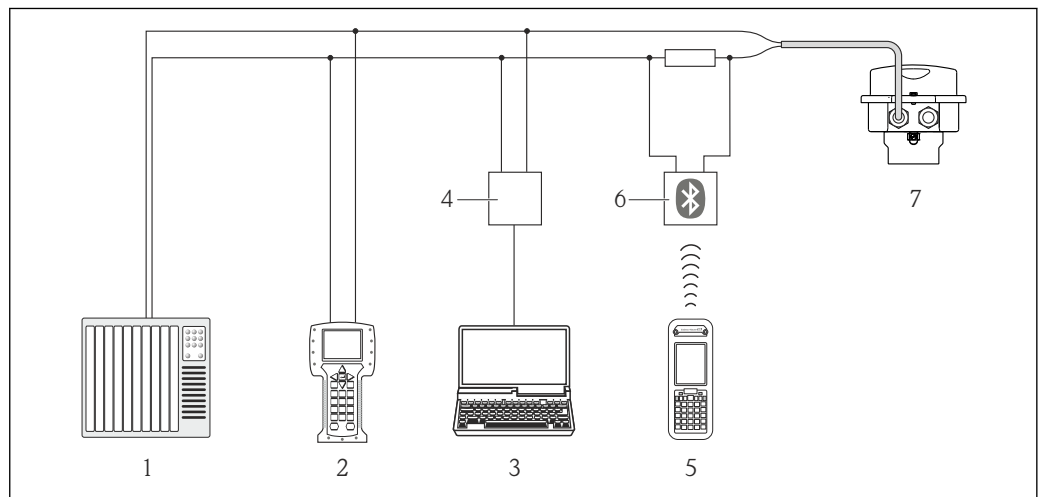


69 Options for remote operation via HART protocol

- 1 Control system (e.g. PLC)
- 2 Field Communicator 475
- 3 Computer with operating tool (e.g. FieldCare, AMS Device Manager, SIMATIC PDM)
- 4 Commubox FXA195 (USB)
- 5 Field Xpert SFX350 or SFX370
- 6 VIATOR Bluetooth modem with connecting cable
- 7 Transmitter

Via Ethernet-based fieldbus

This communication interface is present in the following device version:
 Order code for "Output", option **N**: EtherNet/IP

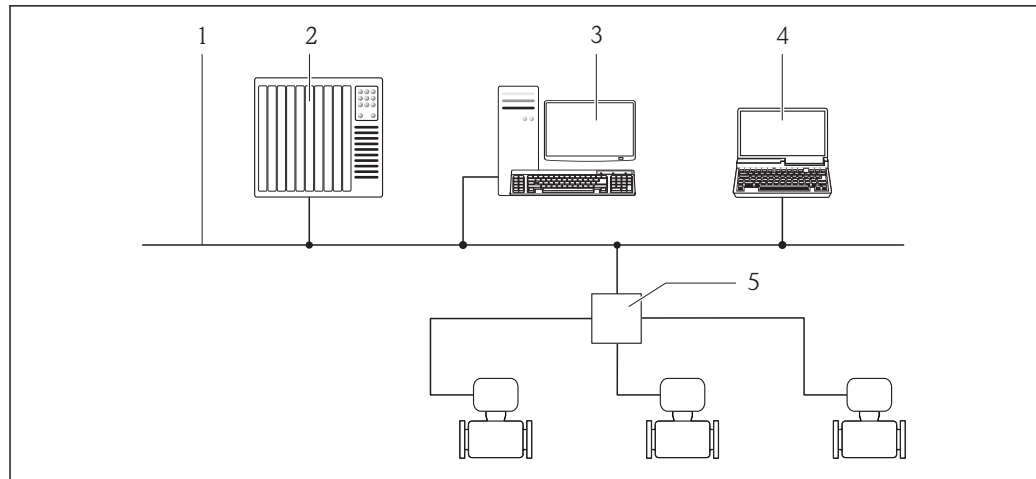


A0016961

- 1 Ethernet network
- 2 Automation system, e.g. "RSLogix" (Rockwell Automation)
- 3 Workstation for measuring device operation: with Add-on Profile Level 3 for "RSLogix 5000" (Rockwell Automation) or with Electronic Data Sheet (EDS)
- 4 Computer with Web browser (e.g. Internet Explorer) for accessing the integrated device Web server or with "FieldCare" operating tool with COM DTM "CDI Communication TCP/IP"
- 5 Ethernet switch

Via PROFIBUS DP network

This communication interface is present in the following device version:
Order code for "Output", option L: PROFIBUS DP

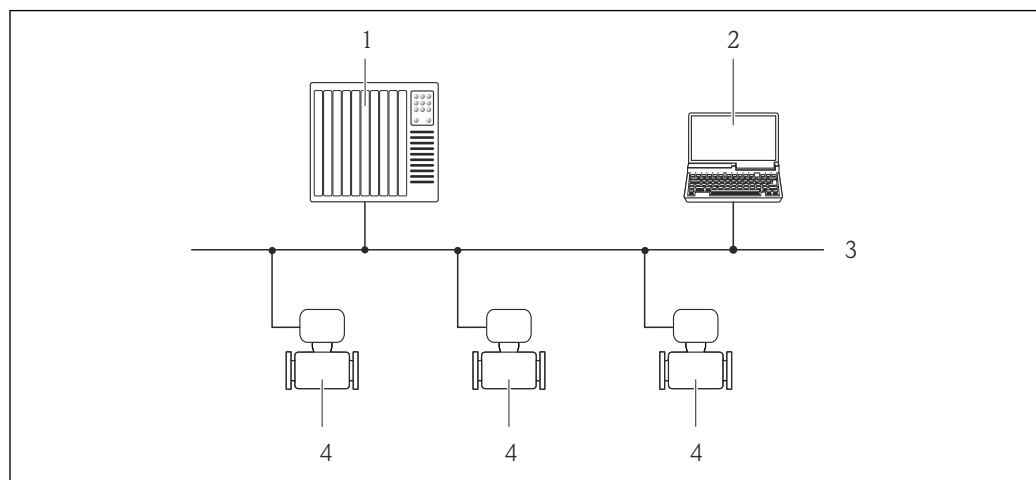

A0020903

- 1 Automation system
- 2 Computer with PROFIBUS network card
- 3 PROFIBUS DP network
- 4 Measuring device


Service interface

Service interface (CDI-RJ45)

This communication interface is present in the following device version:

- Order code for "Output", option B: 4-20 mA HART, pulse/frequency/switch output
- Order code for "Output", option L: PROFIBUS DP
- Order code for "Output", option N: EtherNet/IP

HART

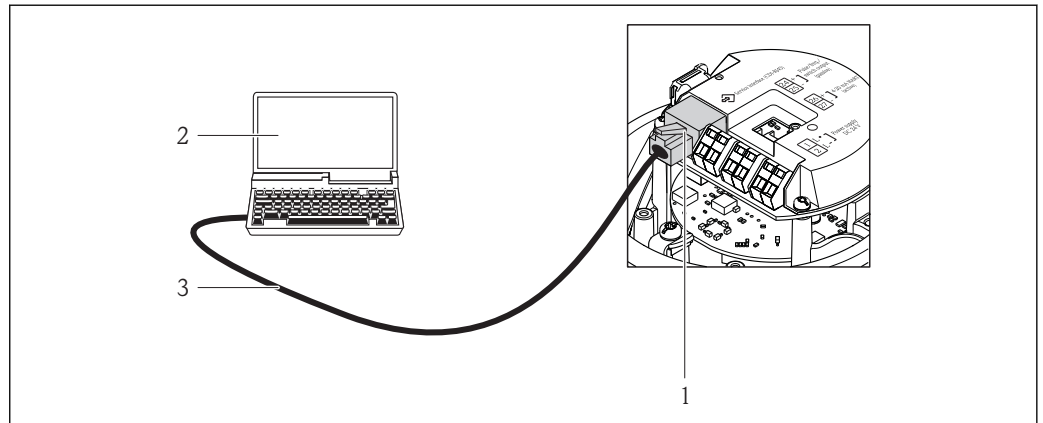


A0016926

70 Connection for the order code for "Output", option B: 4-20 mA HART, pulse/frequency/switch output

- 1 Service interface (CDI -RJ45) of the measuring device with access to the integrated Web server
- 2 Computer with Web browser (e.g. Internet Explorer) for accessing the integrated device Web server or with "FieldCare" operating tool with COM DTM "CDI Communication TCP/IP"
- 3 Standard Ethernet connecting cable with RJ45 plug

PROFIBUS DP

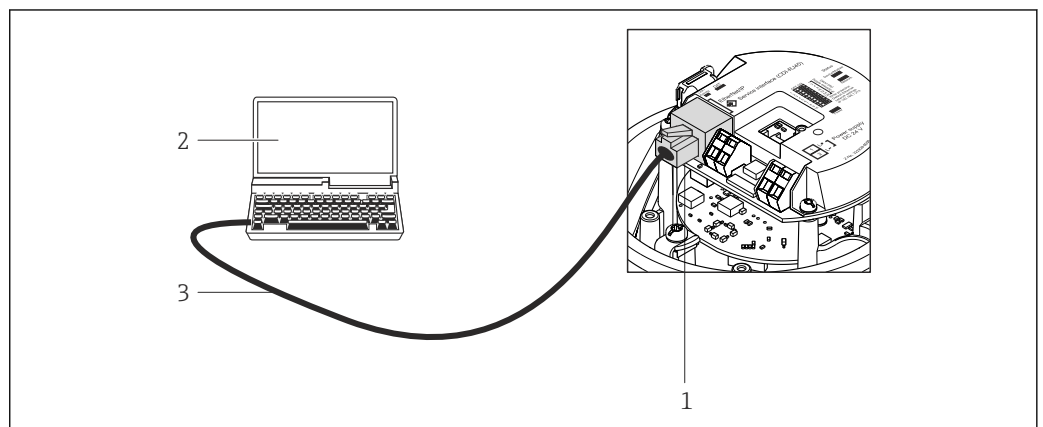


A0021270

71 Connection for order code for "Output", option L: PROFIBUS DP

- 1 Service interface (CDI -RJ45) of the measuring device with access to the integrated Web server
- 2 Computer with Web browser (e.g. Internet Explorer) for accessing the integrated device Web server or with "FieldCare" operating tool with COM DTM "CDI Communication TCP/IP"
- 3 Standard Ethernet connecting cable with RJ45 plug

EtherNet/IP

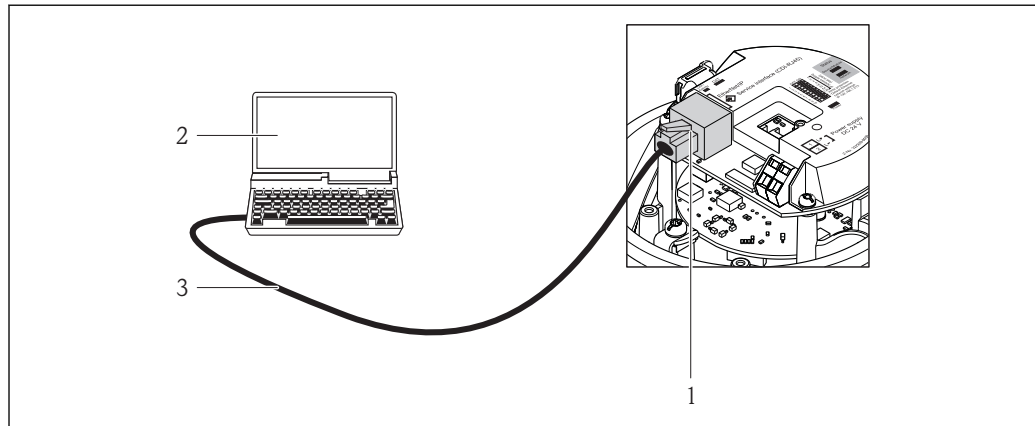


A0016940

72 Connection for order code for "Output", option N: EtherNet/IP

- 1 Service interface (CDI -RJ45) and EtherNet/IP interface of the measuring device with access to the integrated Web server
- 2 Computer with Web browser (e.g. Internet Explorer) for accessing the integrated device Web server or with "FieldCare" operating tool with COM DTM "CDI Communication TCP/IP"
- 3 Standard Ethernet connecting cable with RJ45 plug

Service interface (CDI)

This communication interface is present in the following device version:
Order code for "Output", option **M**: Modbus RS485

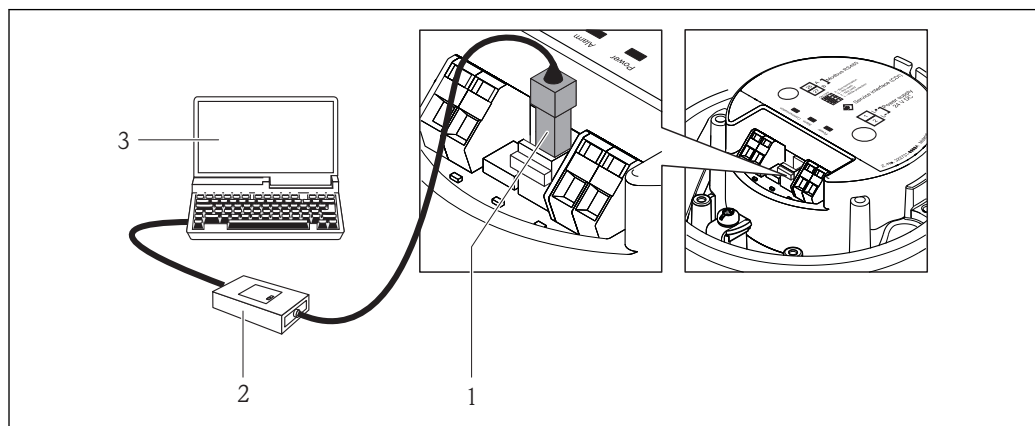

A0016925

- 1 Service interface (CDI) of the measuring device
- 2 Commubox FXA291
- 3 Computer with "FieldCare" operating tool with COM DTM "CDI Communication FXA291"

Certificates and approvals

CE mark

The measuring system is in conformity with the statutory requirements of the applicable EC Directives. These are listed in the corresponding EC Declaration of Conformity along with the standards applied.

Endress+Hauser confirms successful testing of the device by affixing to it the CE mark.

C-Tick symbol

The measuring system meets the EMC requirements of the "Australian Communications and Media Authority (ACMA)".

Ex approval

The measuring device is certified for use in hazardous areas and the relevant safety instructions are provided in the separate "Safety Instructions" (XA) document. Reference is made to this document on the nameplate.



The separate Ex documentation (XA) containing all the relevant explosion protection data is available from your Endress+Hauser sales center.

ATEX/IECEx

Currently, the following versions for use in hazardous areas are available:

Ex ia

Category (ATEX)	Type of protection
II2G	Ex ia IIC T6...T1 Gb
II2G	Ex ia IIC T6...T1 Gb or Ex ia IIB T6...T1 Gb
II1/2G, II2D	Ex ia IIC T6...T1 Ga/Gb or Ex ia IIB T6...T1 Ga/Gb Ex tb IIIC Txxx Db
II2G, II2D	Ex ia IIC T6...T1 Gb or Ex ia IIB T6...T1 Gb Ex tb IIIC Txxx Db

Ex nA

Category (ATEX)	Type of protection
II3G	Ex nA IIC T6...T1 Gc or Ex nA IIC T5-T1 Gc

cCSA_{US}

Currently, the following versions for use in hazardous areas are available:

IS (Ex i)

- Class I Division 1 Groups ABCD
- Class II Division 1 Groups EFG and Class III

NI (Ex nA)

Class I Division 2 Groups ABCD

Hygienic compatibility

- 3A approval
- EHEDG-tested

Certification PROFIBUS

PROFIBUS interface

The measuring device is certified and registered by the PROFIBUS User Organization (PNO). The measuring system meets all the requirements of the following specifications:

- Certified in accordance with PROFIBUS PA Profile 3.02
- The device can also be operated with certified devices of other manufacturers (interoperability)

Modbus RS485 certification

The measuring device meets all the requirements of the MODBUS/TCP conformity test and has the "MODBUS/TCP Conformance Test Policy, Version 2.0". The measuring device has successfully passed all the test procedures carried out and is certified by the "MODBUS/TCP Conformance Test Laboratory" of the University of Michigan.

EtherNet/IP certification

The measuring device is certified and registered by the ODVA (Open Device Vendor Association). The measuring system meets all the requirements of the following specifications:

- Certified in accordance with the ODVA Conformance Test
- EtherNet/IP Performance Test
- EtherNet/IP PlugFest compliance
- The device can also be operated with certified devices of other manufacturers (interoperability)

Pressure Equipment Directive

The devices can be ordered with or without a PED approval. If a device with a PED approval is required, this must be explicitly stated in the order. For devices with nominal diameters less than or equal to DN 25 (1"), this is neither possible nor necessary.

- With the PED/G1/x (x = category) marking on the sensor nameplate, Endress+Hauser confirms compliance with the "Essential Safety Requirements" specified in Annex I of the Pressure Equipment Directive 97/23/EC.
- Devices bearing this marking (PED) are suitable for the following types of medium:
 - Media in Group 1 and 2 with a vapor pressure greater than, or smaller and equal to 0.5 bar (7.3 psi)
 - Unstable gases
- Devices not bearing this marking (PED) are designed and manufactured according to good engineering practice. They meet the requirements of Art.3 Section 3 of the Pressure Equipment Directive 97/23/EC. The range of application is indicated in tables 6 to 9 in Annex II of the Pressure Equipment Directive.

Other standards and guidelines

- EN 60529
Degrees of protection provided by enclosures (IP code)
- IEC/EN 60068-2-6
Environmental influences: Test procedure - Test Fc: vibrate (sinusoidal).
- IEC/EN 60068-2-31
Environmental influences: Test procedure - Test Ec: shocks due to rough handling, primarily for devices.
- EN 61010-1
Safety requirements for electrical equipment for measurement, control and laboratory use
- IEC/EN 61326
Emission in accordance with Class A requirements. Electromagnetic compatibility (EMC requirements).
- NAMUR NE 21
Electromagnetic compatibility (EMC) of industrial process and laboratory control equipment
- NAMUR NE 32
Data retention in the event of a power failure in field and control instruments with microprocessors
- NAMUR NE 43
Standardization of the signal level for the breakdown information of digital transmitters with analog output signal.
- NAMUR NE 53
Software of field devices and signal-processing devices with digital electronics
- NAMUR NE 80
The application of the pressure equipment directive to process control devices
- NAMUR NE 105
Specifications for integrating fieldbus devices in engineering tools for field devices
- NAMUR NE 107
Self-monitoring and diagnosis of field devices
- NAMUR NE 131
Requirements for field devices for standard applications
- NAMUR NE 132
Coriolis mass meter

Ordering information

Detailed ordering information is available from the following sources:

- In the Product Configurator on the Endress+Hauser website: www.endress.com → Select country → Instruments → Select device → Product page function: Configure this product
- From your Endress+Hauser Sales Center: www.endress.com/worldwide

Product Configurator - the tool for individual product configuration

- Up-to-the-minute configuration data
 - Depending on the device: Direct input of measuring point-specific information such as measuring range or operating language
 - Automatic verification of exclusion criteria
 - Automatic creation of the order code and its breakdown in PDF or Excel output format
 - Ability to order directly in the Endress+Hauser Online Shop

Application packages

Many different application packages are available to enhance the functionality of the device. Such packages might be needed to address safety aspects or specific application requirements.


The application packages can be ordered from Endress+Hauser either directly with the device or subsequently. Detailed information on the order code in question is available from your local Endress+Hauser sales center or on the product page of the Endress+Hauser website: www.endress.com.

Heartbeat Technology	
Package	Description
Heartbeat Verification +Monitoring	<p>Heartbeat Monitoring: Continuously supplies monitoring data, which are characteristic of the measuring principle, for an external condition monitoring system. This makes it possible to:</p> <ul style="list-style-type: none"> ▪ Draw conclusions - using these data and other information - about the impact the measuring application has on the measuring performance over time. ▪ Schedule servicing in time. ▪ Monitor the product quality, e.g. gas pockets. <p>Heartbeat Verification: Makes it possible to check the device functionality on demand when the device is installed, without having to interrupt the process.</p> <ul style="list-style-type: none"> ▪ Access via onsite operation or other operating interfaces, such as FieldCare for instance. ▪ Documentation of device functionality within the framework of manufacturer specifications, for proof testing for instance. ▪ End-to-end, traceable documentation of the verification results, including report. ▪ Makes it possible to extend calibration intervals in accordance with operator's risk assessment.








Concentration	
Package	Description
Concentration measurement and special density	<p>Calculation and outputting of fluid concentrations Many applications use density as a key measured value for monitoring quality or controlling processes. The device measures the density of the fluid as standard and makes this value available to the control system. The "Special Density" application package offers high-precision density measurement over a wide density and temperature range particularly for applications subject to varying process conditions.</p> <p>With the help of the "Concentration Measurement" application package, the measured density is used to calculate other process parameters:</p> <ul style="list-style-type: none"> ▪ Temperature-compensated density (reference density). ▪ Percentage mass of the individual substances in a two-phase fluid. (Concentration in %). ▪ Fluid concentration is output with special units (°Brix, °Baumé, °API, etc.) for standard applications. <p>The measured values are output via the digital and analog outputs of the device.</p>

Accessories

Various accessories, which can be ordered with the device or subsequently from Endress+Hauser, are available for the device. Detailed information on the order code in question is available from your local Endress+Hauser sales center or on the product page of the Endress+Hauser website: www.endress.com.



Device-specific accessories	
For the sensor	
Accessories	Description
Heating jacket	<p>Is used to stabilize the temperature of the fluids in the sensor. Water, water vapor and other non-corrosive liquids are permitted for use as fluids. If using oil as a heating medium, please consult with Endress+Hauser.</p> <p> For details, see Operating Instructions BA00099D</p>

Communication-specific accessories



Accessories	Description
Commubox FXA195 HART	For intrinsically safe HART communication with FieldCare via the USB interface.  For details, see "Technical Information" TI00404F
HART Loop Converter HMX50	Is used to evaluate and convert dynamic HART process variables to analog current signals or limit values.  For details, see "Technical Information" TI00429F and Operating Instructions BA00371F
Wireless HART adapter SWA70	Is used for the wireless connection of field devices. The WirelessHART adapter can be easily integrated into field devices and existing infrastructures, offers data protection and transmission safety and can be operated in parallel with other wireless networks with minimum cabling complexity.  For details, see Operating Instructions BA00061S
Fieldgate FXA320	Gateway for the remote monitoring of connected 4-20 mA measuring devices via a Web browser.  For details, see "Technical Information" TI00025S and Operating Instructions BA00053S
Fieldgate FXA520	Gateway for the remote diagnostics and remote configuration of connected HART measuring devices via a Web browser.  For details, see "Technical Information" TI00025S and Operating Instructions BA00051S
Field Xpert SFX350	Field Xpert SFX350 is a mobile computer for commissioning and maintenance. It enables efficient device configuration and diagnostics for HART and FOUNDATION Fieldbus devices in the non-Ex area .  For details, see Operating Instructions BA01202S
Field Xpert SFX370	Field Xpert SFX370 is a mobile computer for commissioning and maintenance. It enables efficient device configuration and diagnostics for HART and FOUNDATION Fieldbus devices in the non-Ex area and the Ex area .  For details, see Operating Instructions BA01202S

Service-specific accessories


Accessories	Description
Applicator	Software for selecting and sizing Endress+Hauser measuring devices: <ul style="list-style-type: none"> ■ Calculation of all the necessary data for identifying the optimum flowmeter: e.g. nominal diameter, pressure loss, accuracy or process connections. ■ Graphic illustration of the calculation results Administration, documentation and access to all project-related data and parameters over the entire life cycle of a project. Applicator is available: <ul style="list-style-type: none"> ■ Via the Internet: https://wapps.endress.com/applicator ■ On CD-ROM for local PC installation.
W@M	Life cycle management for your plant W@M supports you with a wide range of software applications over the entire process: from planning and procurement, to the installation, commissioning and operation of the measuring devices. All the relevant device information, such as the device status, spare parts and device-specific documentation, is available for every device over the entire life cycle. The application already contains the data of your Endress+Hauser device. Endress+Hauser also takes care of maintaining and updating the data records. W@M is available: <ul style="list-style-type: none"> ■ Via the Internet: www.endress.com/lifecyclemanagement ■ On CD-ROM for local PC installation.

FieldCare	FDT-based plant asset management tool from Endress+Hauser. It can configure all smart field units in your system and helps you manage them. By using the status information, it is also a simple but effective way of checking their status and condition.  For details, see Operating Instructions BA00027S and BA00059S
Commubox FXA291	Connects Endress+Hauser field devices with a CDI interface (= Endress+Hauser Common Data Interface) and the USB port of a computer or laptop.  For details, see "Technical Information" TI00405C

System components

Accessories	Description
Memograph M graphic display recorder	The Memograph M graphic display recorder provides information on all relevant measured variables. Measured values are recorded correctly, limit values are monitored and measuring points analyzed. The data are stored in the 256 MB internal memory and also on a SD card or USB stick.  For details, see "Technical Information" TI00133R and Operating Instructions BA00247R
iTEMP	The temperature transmitters can be used in all applications and are suitable for the measurement of gases, steam and liquids. They can be used to read in the fluid temperature.  For details, see "Fields of Activity", FA00006T

Documentation

-  For an overview of the scope of the associated Technical Documentation, refer to the following:
- The CD-ROM provided for the device (depending on the device version, the CD-ROM might not be part of the delivery!)
 - The *W@M Device Viewer* : Enter the serial number from the nameplate (www.endress.com/deviceviewer)
 - The *Endress+Hauser Operations App*: Enter the serial number from the nameplate or scan the 2-D matrix code (QR code) on the nameplate.

Standard documentation

Brief Operating Instructions

Measuring device	Documentation code
Promass P 100	KA01118D

Operating Instructions

Measuring device	Documentation code			
	HART	PROFIBUS DP	Modbus RS485	EtherNet/IP
Promass P 100	BA01192D	BA01253D	BA01059D	BA01067D

Supplementary device-dependent documentation

Safety Instructions


Contents	Documentation code
ATEX/IECEX Ex i	XA00159D
ATEX/IECEX Ex nA	XA01029D
cCSAus IS	XA00160D
INMETRO Ex i	XA01219D
INMETRO Ex nA	XA01220D

Contents	Documentation code
NEPSI Ex i	XA01249D
NEPSI Ex nA	XA01262D

Special Documentation

Contents	Documentation code
Information on the Pressure Equipment Directive	SD00142D
Modbus RS485 Register Information	SD00154D
Concentration Measurement	SD01152D
Heartbeat Technology	SD01153D

Installation Instructions

Contents	Documentation code
Installation Instructions for spare part sets	Specified for each individual accessory (→  89)

Registered trademarks

HART®

Registered trademark of the HART Communication Foundation, Austin, USA

PROFIBUS®

Registered trademark of the PROFIBUS User Organization, Karlsruhe, Germany

Modbus®

Registered trademark of SCHNEIDER AUTOMATION, INC.

EtherNet/IP™

Trademark of ODVA, Inc.

Microsoft®

Registered trademark of the Microsoft Corporation, Redmond, Washington, USA

TRI-CLAMP®

Registered trademark of Ladish & Co., Inc., Kenosha, USA

Applicator®, FieldCare®, Field Xpert™, HistoROM®, TMB®, Heartbeat Technology™

Registered or registration-pending trademarks of the Endress+Hauser Group





www.addresses.endress.com
